



**PASCO** capstone™

# User's Guide

Version 1.1.3

## PASCO Capstone User's Guide

Document Number: 012-14016B

Capstone Version: 1.1.3

### Limited Warranty

For a description of the product warranty, see the PASCO scientific catalog.

### Copyright

This User's Guide is copyrighted with all rights reserved. Permission is granted to non-profit educational institutions for reproduction of any part of this manual, providing the reproductions are used only in their laboratories and classrooms, and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited.

### Trademarks

PASCO, PASCO scientific, PASCO Capstone, DataStudio, PASPORT, SPARK, SPARK Science Learning System, SPARKlab, SPARKbook, SPARKvue, SPARKvue HD, Xplorer, Xplorer GLX are trademarks or registered trademarks of PASCO scientific, in the United States and/or in other countries. All other brands, products, or service names are or may be trademarks or service marks of, and are used to identify products or services of their respective owners. For more information visit [www.pasco.com/legal](http://www.pasco.com/legal).

### Software Licenses

All required licenses for software components of Capstone can be found on the Flash Drive, CD-ROM, or DVD-ROM included with the product or are included in the software download. To obtain source code for GPL/LGPL licensed software components, contact PASCO scientific by phone: 1-800-772-8700 (in the U.S), +1 916 786 3800 (worldwide), or by email: [support@pasco.com](mailto:support@pasco.com)

## Teacher and Technical Support

For questions about PASCO Capstone and other PASCO scientific products that are not answered in the Help or User's Guide, you can contact the PASCO Teacher and Technical Support staff by phone, email, or on the Web.

Phone:	<b>1-800-772-8700 (in the U.S.)</b>
	<b>+1 916 786 3800 (worldwide)</b>
Email:	<a href="mailto:support@pasco.com"><u>support@pasco.com</u></a>
Web:	<a href="http://www.pasco.com/support/"><u>www.pasco.com/support/</u></a>
Tech Notes:	<a href="http://www.pasco.com/support/technical-support/technote/"><u>www.pasco.com/support/technical-support/technote/</u></a>

 See "Where can I find help documentation and curriculum for PASCO Capstone?":  
147

# Contents

PASCO Capstone User's Guide .....	i
Teacher and Technical Support .....	ii
<b>1 Tour the Interface .....</b>	<b>1</b>
Tour the Menu and Toolbar .....	2
Tour the Page Tools .....	5
Tour the Tools Palette .....	6
Tour the Controls Palette .....	8
Tour the Displays Palette .....	10
Tour the Display Area .....	13
Tour Context Help .....	14
Tour the Scope .....	15
<b>2 Getting Started .....</b>	<b>17</b>
Connecting an interface .....	17
Connecting an 850 Universal or 750 USB Interface .....	21
Connecting Bluetooth Interfaces: such as the SPARKlink Air or AirLink2 .....	22
Connecting a 750 SCSI or 500 serial Interface .....	24
Connecting USB PASPORT Interfaces: such as the SPARK SLS, SPARKlink, or Power Link .....	25
Connecting an Xplorer GLX or Xplorer .....	26
Connecting a sensor .....	27
Connecting a sensor with the 850 Interface .....	36
Connecting a sensor with the 750 Interface .....	41
Connecting a sensor with the SPARK SLS, SPARKlink, SPARKlink Air .....	45
Connecting a sensor with the Xplorer GLX .....	50
Connecting a sensor with the Digital Adapter or Photogate Port .....	55
Connecting a sensor with the Analog Adapter .....	56



Connecting a sensor with the AirLink2 .....	56
Adding a display .....	56
Starting an experiment .....	60
Downloading Capstone experiments .....	61
<b>3 How do I...</b> .....	<b>63</b>
How do I... Installation and Update .....	66
How do I automate a network installation? .....	66
How do I force an update to an interface? .....	66
How do I... Experiment Files .....	67
How do I download and open PASCO ready-made experiments? .....	67
How do I save my experiment? .....	68
How do I export all data? .....	68
How do I open DataStudio experiments and data? .....	68
How do I open Xplorer GLX, SPARK SLS, and other non-Capstone exper- iments and data? .....	69
How do I import data from Microsoft Excel*? .....	69
How do I export data to Microsoft Excel*? .....	70
How do I import Xplorer GLX data? .....	71
How do I... Data and Measurements .....	73
How do I record data in PASCO Capstone? .....	73
How do I delete data runs? .....	79
How do I choose one or more runs to display? .....	79
How do I identify a measurement when I have multiple of the same type? .....	80
How do I change the numerical format of data? .....	81
How do I change the numerical format of statistics? .....	83
How do I exclude, delete, and restore selected data? .....	84
How do I edit a measurement? .....	87
How do I create a pure mathematical function? .....	87

How do I... Workbooks .....	89
How do I show, hide, and lock the palettes? .....	89
How do I change what display toolbar buttons are visible? .....	90
How do I... Sensors .....	92
How do I zero a measurement? .....	92
How do I switch the sign on a sensor? .....	93
How do I calibrate a sensor? .....	94
How do I calibrate a motion sensor? .....	95
How do I calibrate a drop counter? .....	96
How do I calibrate a pH sensor? .....	97
How do I... Calculations .....	99
How do I create a calculation? .....	99
How do I use my calculation? .....	102
How do I select a measurement in the calculator? .....	103
How do I see the result of a calculation? .....	103
How do I reorder calculations? .....	105
How do I... Plots .....	106
How do I plot multiple measurements on separate plot areas with shared X-axis? .....	107
How do I zoom to a selection? .....	108
How do I pan and scale? .....	110
How do I select active data? .....	111
How do I plot multiple measurements on same plot with multiple Y-axes? ....	113
How do I plot multiple measurements on same plot with same Y-axes? .....	114
How do I apply a curve fit to a portion of the data? .....	117
How do I change curve fit parameters? .....	118
How do I adjust the weights for the weighted linear fit? .....	119
How do I lock and unlock the axes of a graph? .....	121


How do I change layer order of a display? .....	123
How do I... Tables .....	124
How do I reorder columns in a table? .....	124
How do I use a table to build data in Keep Mode? .....	124
How do I... Text Boxes .....	127
How do I enter subscripts, superscripts, Greek letters, and symbols? .....	127
How do I enter an equation into a text display? .....	127
How do I... Scope .....	130
How do I compare scope traces from the same sensor when monitoring? ....	130
How do I select the scope trigger? .....	131
How do I pre-trigger? .....	134
How do I freeze and record the scope display? .....	135
How do I... Video .....	137
How do I add a video source? .....	137
How do I analyze a video? .....	138
How do I manually synchronize video with data? .....	140
How do I adjust the frame rate of a recorded video? .....	141
How do I adjust the resolution of a video? .....	141
How do I make files with videos smaller? .....	141
How do I troubleshoot video problems? .....	142
<b>4 Frequently Asked Questions (FAQ) .....</b>	<b>145</b>
FAQ... PASCO Capstone .....	147
What is PASCO Capstone? .....	147
Where can I find help documentation and curriculum for PASCO Capstone?	147
What are the new features in the latest PASCO Capstone? .....	148
Can PASCO Capstone open DataStudio, SPARKlab, and GLX files? .....	151
Can Capstone files be opened on computers with different operating systems? .....	151

How does Capstone compare to other PASCO applications? .....	152
FAQ... Installation and Updates .....	160
What licenses are required? .....	160
Is an MSI available for PASCO Capstone? .....	160
Do I need to update the interfaces? .....	161
Do I need to update the sensor datasheets? .....	161
FAQ... Setup .....	162
What do I need to have a complete measurement system? .....	162
Can sensors sample at different rates? .....	164
FAQ... Data files .....	165
Where is my data saved? .....	165
What is the file export format? .....	165
What is the format for text import? .....	166
FAQ... Data and Measurements .....	167
Which data is recorded? .....	167
Can sensors sample at different rates? .....	167
What is Continuous Mode? .....	168
What is Fast Monitor Mode? .....	169
Is there a way to delete data points? .....	170
Which data is active? .....	172
How are the curve fit and smoothing calculations determined? .....	172
FAQ... Calculations and Calibration .....	174
What is the format for calculations? .....	174
What are the equations for calibration? .....	174
1 Point Offset .....	175
1 Point Slope .....	175
2 Point Linear .....	175
Can multiple measurements be calibrated simultaneously? .....	175

---

FAQ... Displays .....	176
What is panning and scaling? .....	176
Why are the display toolbars not active? .....	177
FAQ... Scope .....	178
What is the scope? .....	178
What does the Single Shot button do? .....	178
What is pre-triggering in the scope? .....	179
FAQ... Video and Text .....	180
What cameras and video formats are supported? .....	180
Why does the text color remain black after picking a color? .....	180

# 1 Tour the Interface

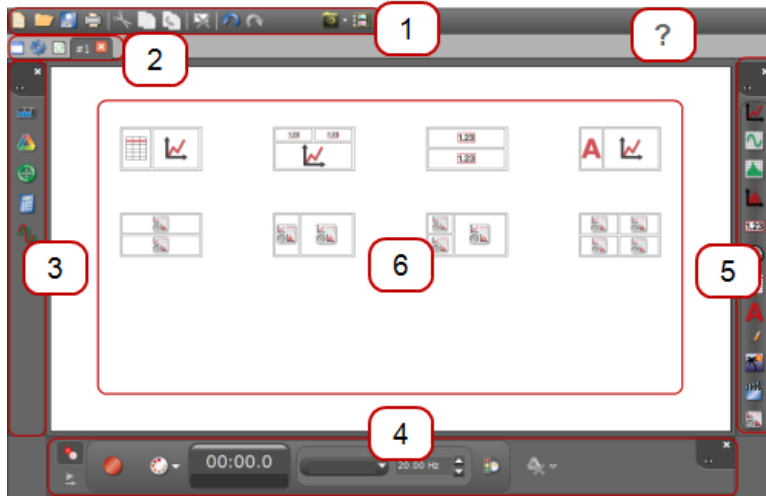

Click to learn more

[1 Menu and Toolbar](#)

[2 Page Tools](#)

[3 Tools Palette](#)

[? Context Help](#)

[4 Controls Palette](#)

[5 Displays Palette](#)


[6 Display Area](#)

 [Online Videos](#)

## Tour the Menu and Toolbar










The Capstone menu includes common application functions and application specific functions. Some functions are available as toolbar buttons.

Descriptions for Capstone Menu and Toolbar functions:  **Tip<sup>1</sup>**

### 1 File

Menu item	Description
Import Data...	Import data into the Capstone data summary.
Export Data...	Export collect data to a text file.
Preferences	Set global preferences including: <ul style="list-style-type: none"> <li>• Preference File</li> <li>• General Application</li> <li>• Language</li> <li>• Workbook</li> <li>• Displays Palette</li> <li>• Tools Palette</li> <li>• Controls Palette</li> </ul>

### 2 Edit





Icon	Menu item	Description
	Undo	Undo the last operation.  <b>Tip<sup>2</sup></b>
	Redo	Redo an Undo operation  <b>Tip<sup>2</sup></b>
	Cut	Cut the selected display element
	Copy	Copy the selected display element
	Paste	Paste a Cut or Copied display element

<sup>1</sup>Capstone includes descriptions within dialog boxes. Click on an item for a brief description.

<sup>2</sup>The operation that will be undone is listed in the menu.

<sup>2</sup>The operation that will be re-done is listed in the menu.

### 3 Workbook










Icon	Menu item	Description
	Add Page	Add Page To Workbook
	Delete Page	Delete this Page
	Rename Page	Rename Current Workbook Page
	Lock All Pages	Lock all current pages for editing. Pages can still be added and will not be locked; to lock added pages, click Lock All Pages.
	Unlock All Pages	Unlock pages to allow editing.
	Edit Master Page	Opens the Master Page Add text and graphics to appear in the background of other pages.
	Workbook Page Properties	Set workbook preferences, including: <ul style="list-style-type: none"> <li>• Page Name</li> <li>• Snap to Grid</li> <li>• Show Tools Palette</li> <li>• Show Displays Palette</li> <li>• Show Controls Palette</li> <li>• Page Shape</li> </ul>

### 4 Display

Icon	Menu item	Description
	Display	Shows tools that apply to the selected display.




5 Journal

Icon	Menu item	Description
	Take Journal snapshot   Select snapshot mode	Takes an image and places it in the journal. Select: Snapshot Entire Workbook Page Snapshot Workbook Page Content Snapshot Workbook Page Display
	Show or hide Journal	Lists the journal snapshots and allows:  Show or hide annotations  Print Journal  Export Journal to HTML  View full-size snapshot  Delete selected snapshot  Move selected snapshot up  Move selected snapshot down


6 Help

Icon	Menu item	Description
	PASCO Capstone Help	Opens the Capstone Help System in a separate browser. A PDF version of the User's Guide is also available from the Help System. <a href="#">See "Where can I find help documentation and curriculum for PASCO Capstone?": 147</a>
	About PASCO Capstone	Opens a dialog with version information and the ability enter a license key.

7 Delete Display

Icon	Menu item	Description
	Delete Selected Display	Deletes the active display on the workbook page. Data are not affected.

## Tour the Page Tools

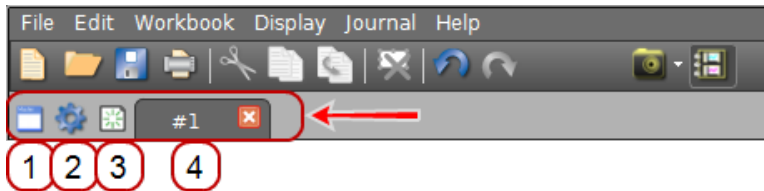

See "How do I... Workbooks": 89.

### 1 Edit Master Page

Icon	Description
	Master Page Add text and graphics to appear in the background of other pages.

### 2 Properties

Icon	Description
	Change properties of current page and Tools Palette

### 3 Add Page

Icon	Description
	Adds a new page to the workbook
	Drag and drop page tabs to rearrange the workbook pages.

See "Adding a display ": 56.


### 4 Rename / Delete

Icon	Description


- 1 Double click to rename the page.
- 2 Deletes the current page from the workbook.

## Tour the Tools Palette

×




**1 Hardware Setup**

Icon	Description
	Configure interface and sensor hardware
	<a href="#">See "Connecting an interface": 17.</a>


---

**2 Timer Setup**

Icon	Description
	Create timers for photogates and other timing devices


---

**3 Data Summary**

Icon	Description
	Manage experiment data and data properties
	<a href="#">See "How do I... Data and Measurements": 73.</a>


---

**4 Calibration**

Icon	Description
	Calibrate sensor measurements
	<a href="#">See "How do I... Sensors": 92.</a>


---

**5 Calculator**

Icon	Description
	Define new data sets as functions of measurements and other data
	<a href="#">See "How do I... Calculations": 99.</a>


---

**6 Signal Generator**


Icon	Description
	Control interface Signal Generator output

---

**7 Digital Output**


Icon	Description
	Specify behavior for expansion port inputs and/or outputs

**8** Curve Fit Editor



Icon	Description
	Edit curve fits created in Graph displays <a href="#">See "How do I change curve fit parameters?": 118.</a>

## Tour the Controls Palette

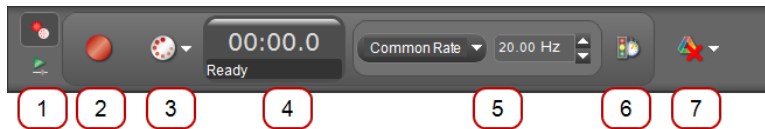

The Controls Palette is used to start/stop data collection, set the data monitoring mode, set the sampling rate, delete data runs, and define Start/Stop conditions.







### 1 Record / Play Selection

Icon	Description
Toggle between:	
	Recording Mode Use to switch to recording Controls Palette
	Playback Mode Use to switch to playback Controls Palette

### 2 Record / Monitor

Icon	Description
Capstone toggles between:	
	Record Begin recording all measurements from each connected sensor
	Monitor Start monitoring measurements from connected sensors

### 3 Record Mode Selection

Icon	Description
Toggle between:	
	Continuous Mode Set to record data at the current sample rate settings
	Keep Mode Set to keep data points when commanded
	Fast Monitor Mode Set to acquire data without recording; required for high speed sampling in Scope and FFT displays
 See " <a href="#">How do I record data in PASCO Capstone?</a> ": 73	

4 Record Timer

Description



- 1 Displays the recording time as reported by the sensors in **minutes:seconds** format.
- 2 Displays the recording status: Ready or Recording

5 Sampling Source and Sample Rate

Description



- 1 Sampling rate source: single sensor or Common Rate

**Note:** Common Rate is applied to all sensors. The Sample Rate range for Common Rate is intersection of all sensor ranges.

- 2 Adjust the Sample Rate



6 Start / Stop Conditions

Icon Description


-  Set automatic Start/Stop conditions

7 Delete Runs

Icon Description



-  Delete most recent run
-  Delete All Runs or a specific run

## Tour the Displays Palette



Drag to page or double-click to create a display ( See "Adding a display ": 56).




### 1 Graph

Icon	Description
	Create a line graph to plot captured data or user defined data. Several tools are available for graphs including statistics, curve fit, slope tool, and smoothing.   See "How do I... Plots": 106.


### 2 Scope

Icon	Description
	Create a display that functions as a digital storage oscilloscope, which is useful for viewing measurements that oscillate rapidly in time.   See "Tour the Scope": 15


### 3 FFT

Icon	Description
	Create a spectral analysis display, which uses a Fast Fourier Transform to measure the relative signal intensity as a function of frequency.


### 4 Histogram

Icon	Description
	Create a histogram bar graph. The height of each of the bars is proportional to the number of data that fall within the range of values that the bar represents. Several tools are available for histograms including statistics, setting bin sizes, and overlaying.


### 5 Digits

Icon	Description
	Create a digits display to show the current data value or a statistical value.

## 6 Meter




Icon	Description
	Create a meter display to indicate the current data value (or statistical value) by the angular displacement of a needle against a dial indicator. Several meter styles are available.

## 7 Table




Icon	Description
	Create a table to list captured data or user defined data in columns.  <b>Note:</b> Tables are used to enter user-defined numerical or text data. Several tools are available for tables including statistics

 See "How do I... Tables": 124.


## 8 Text Box

Icon	Description
	Create a text box for explanatory text and images. The text box can be locked for editing.  <b>Tip</b> <sup>1</sup>   See "How do I... Text Boxes": 127.

## 9 Text Entry Box

Icon	Description
	Create a text entry box for user input. The input can be linked to other text entry boxes.  <b>Tip</b> <sup>1</sup>   See "How do I... Text Boxes": 127.

## 10 Image



Icon	Description
	Create a box to insert an image from a file. Note: Use the Movie icon to capture an image from an attached USB camera.)

<sup>1</sup>Use a Text Entry Box for user student input on a locked experiment.


<sup>1</sup>Use the linkage to present a previously entered student response on a subsequent page of the Workbook. For example, to combine all the answers on a single page.



11 Movie


Icon	Description
	<p>Create a box to insert a movie file, capture an image from a video camera or microscope connected through a USB port, record a stand-alone video using a USB webcam, or use a USB video camera as a sensor to record a movie along with other sensor data.</p>
	<p>A special video analysis mode accessed from the toolbar provides tools for analyzing videos and image snapshots.</p>
	<p> See "<a href="#">How do I... Video</a>": 137.</p>

12 Placeholder

Icon	Description
	<p>Create a spacer for a display to be defined at a later time or create white space.</p>

## Tour the Display Area


The PASCO Capstone Workbook pages facilitate quick and easy layouts and support highly flexible page designs.

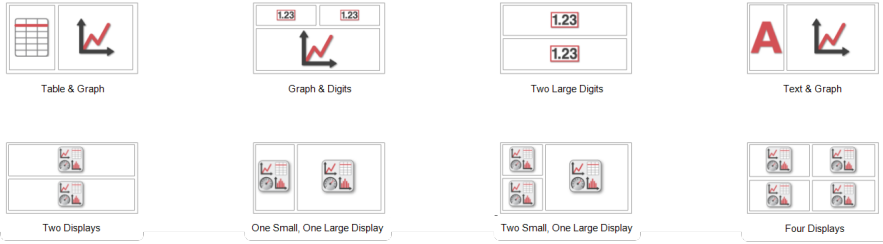Features	See
Drag and drop targets	<a href="#">"Adding a display": 56</a>
Easy resize and alignment tools	<a href="#">"Adding a display": 56</a>
QuickStart templates	<a href="#">"Adding a display": 56</a>
<p data-bbox="224 611 553 646"><b>QuickStart Templates</b></p> 	<a href="#">"Adding a display": 56</a>
Easy table creation, including calculated data	<a href="#">"How do I... Tables": 124</a> <a href="#">"How do I... Calculations": 99</a>
Text boxes with equations copied from word processor	<a href="#">"How do I... Text Boxes": 127</a>

## Tour Context Help

Some items in PASCO Capstone include helpful information at the bottom of dialog windows. To see the help: Click on an item for a brief description.

### Context Help Example: File Properties dialog

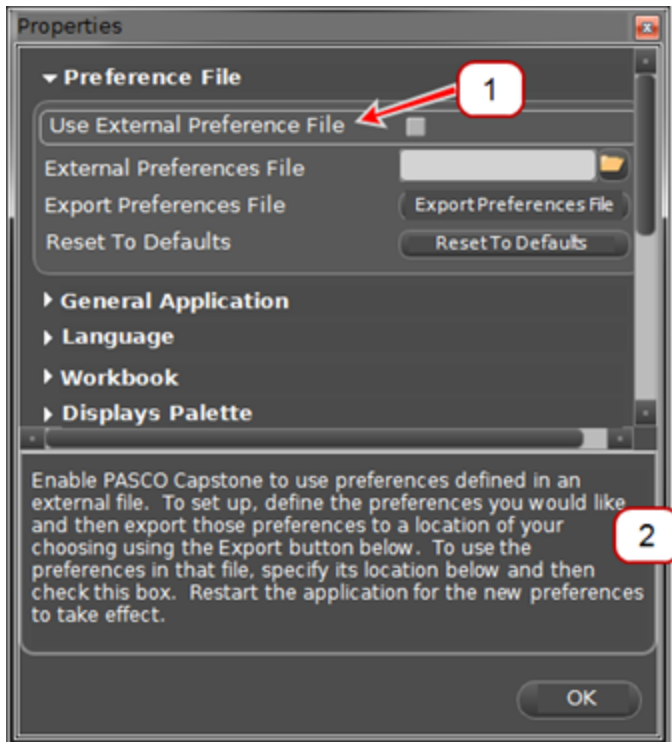


- 1 Click on an item for a brief description.
- 2 Information displays here.

## Tour the Scope

















### Overview

The scope is designed to view high-frequency oscillating data or single high-speed events that the graph display may not be able to capture.

**Maximum Sampling Rate:** The maximum supported sampling rate is a function of the computer-sensor interface that is connected to the computer (for example: 850 Universal Interface → 10 MHz, ScienceWorkshop 750 → 250 kHz).

Features	 See
Continuous Mode and Fast Monitor Mode	"What is Continuous Mode?": 168 "What is Fast Monitor Mode?": 169
Single shot	"What does the Single Shot button do?": 178
Multiple triggers	"How do I select the scope trigger?": 131
Pre-trigger	"How do I pre-trigger?": 134
Image freeze	"How do I freeze and record the scope display?": 135

### Toolbar

Icon	Description
	Adjust y-axis scale to fit data
	Activate and control scope trigger
	Stop collection after one trace
	Automatically adjust sample rate based on time-axis scale
	Activate to view multiple runs; Select visible run(s)
	Increase trace offset
	Set trace offset value to zero
	Decrease trace offset
	Show data coordinates and access Delta Tool
	Creates a data set from active traces
	Increase number of data points in trace
	Decrease number of data points in trace
	Add new y-axis to scope display
	Remove active element or axis
	Allow rearrangement of axes
	Pin toolbar to display




Properties



Show or hide tools

---

 See "Adding a display ": 56

## 2 Getting Started

See the following topics setting up and run a simple experiment:

"Connecting an interface": 17

"Connecting a sensor": 27


"Adding a display ": 56

"Starting an experiment": 60

"Downloading Capstone experiments": 61

 See "How do I...": 63 for detailed instructions.

### Connecting an interface

Capstone software running on your computer receives data from one or more connected sensors through one or more interfaces. (At least one interface and sensor is required to collect data.  See "What do I need to have a complete measurement system? ": 162)

Capstone is compatible with all PASCO USB interfaces and some non-USB interfaces using a serial-to-USB convertor.  **Tip<sup>1</sup>**

#### Automatically detect an interface

1. Connect one or more PASCO USB interfaces to the computer.  **Tip<sup>2</sup>**




"Connecting an 850 Universal or 750 USB Interface": 21

"Connecting a 750 SCSI or 500 serial Interface": 24



"Connecting Bluetooth Interfaces: such as the SPARKlink Air or AirLink2": 22

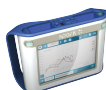

"Connecting USB PASPORT Interfaces: such as the SPARK SLS, SPARKlink, or Power Link": 25



"Connecting an Xplorer GLX or Xplorer": 26


2. Add one or sensors.  See "Connecting a sensor": 27)


<sup>1</sup>You can add interfaces that are not connected to the computer. This is useful to author or prepare experiments in advance.

<sup>2</sup>Interfaces can be connected to your computer with Capstone running. Capstone will automatically add recognized interfaces such as the 850 Universal Interface.

## Manually add an interface


1. Click **Hardware Setup** in the Tools Palette: .

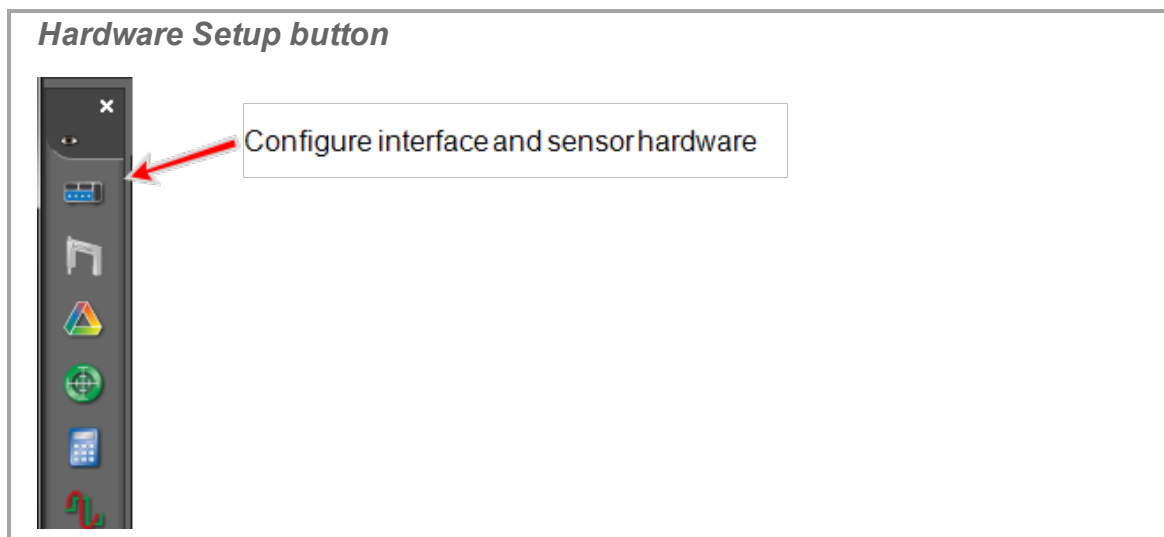

2. Add the interface if not displayed in the **Hardware Setup** list.  **Tip<sup>1</sup>**
3. If the interface is not automatically detected, click the **Choose Interface** button.
4. The Choose Interface dialog opens.


<sup>1</sup>Some PASCO interfaces, such as the PASCO 850 and SPARK, are automatically detected by Capstone when plugged in and turned on.

**Choose Interface**

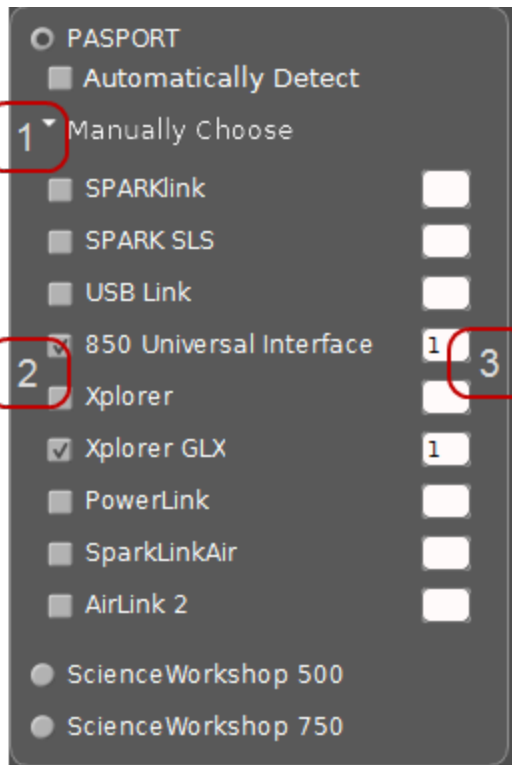

- 1 Click **Manually Choose**.
- 2 Click the interfaces to add or remove.
- 3 Optional: change the default of 1 interface to add multiple interfaces of the same type.



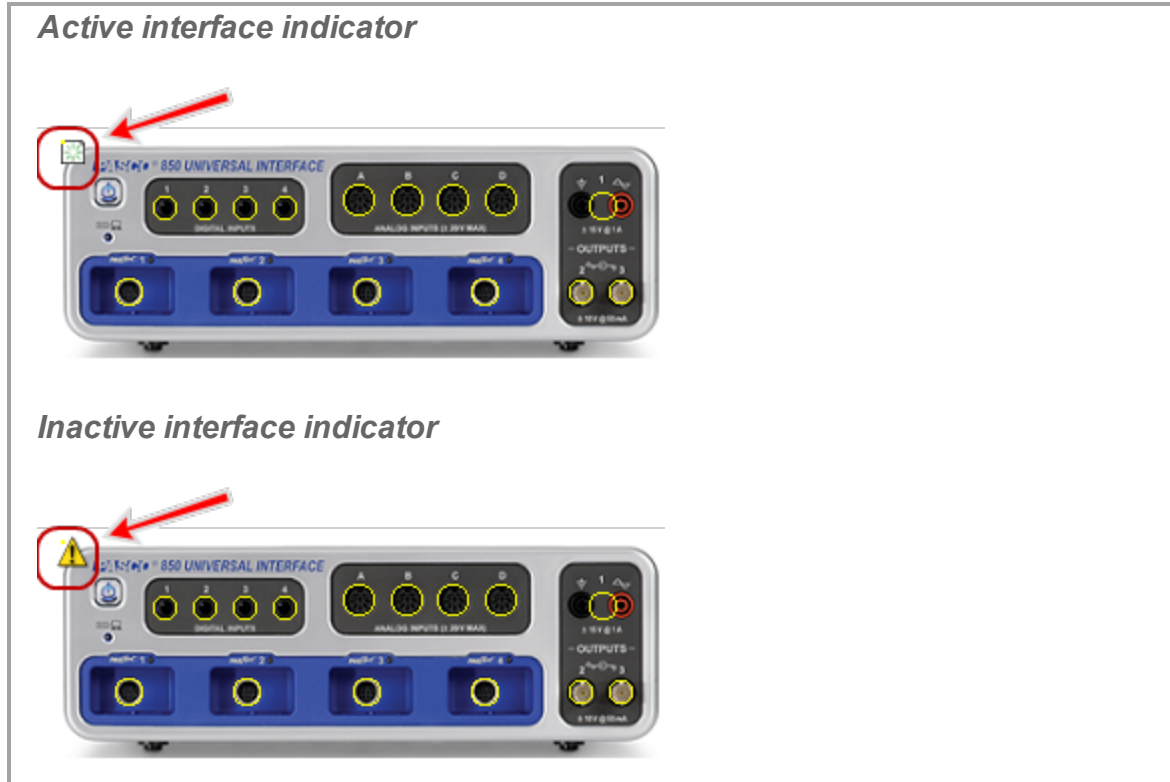


5. Add one or sensors. (See "Connecting a sensor": 27)


Connected and added interfaces are displayed in the **Hardware Setup** dialog.

**Note:** Interfaces that are connected to the computer, powered on, and recognized by Capstone display (☒) and interfaces that are not active display (⚠).

*Active and inactive interfaces and sensors in Hardware Setup*

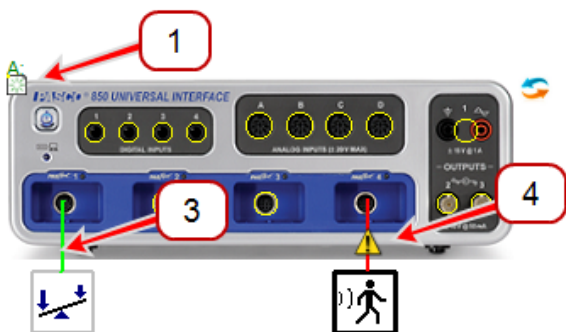

- 1 Active interface
- 2 Inactive interface (disconnected or off)
- 3 Active sensor
- 4 Inactive sensor (disconnected or off)

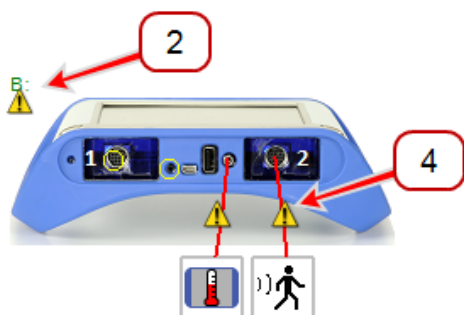

➔ Next step: **"Connecting a sensor": 27**

## Connecting an 850 Universal or 750 USB Interface

The PASCO 850 Universal Interface includes four digital inputs, four analog ports, four PASPORT sensor ports, a 15 W function generator, dual high speed function generators.

### Procedure

1. Use the A-to-mini-B USB cable to connect the smaller USB port of the 850 Interface to a USB port on your computer (or a USB hub connected to the computer).
2. Connect the AC adapter to a wall outlet and to the AC adapter port on the rear of the 850 Universal Interface.

**850 Universal Interface: USB and Power**



3. Press and hold the power button () until the green lights on the front flash.

**850 Universal Interface: On button**



4. Configure the interface for Capstone. ([See "Connecting an interface": 17](#))
5. Add sensors. ([See "Connecting a sensor with the 850 Interface": 36](#))

**Next step: "Connecting a sensor": 27**

## Connecting Bluetooth Interfaces: such as the SPARKlink Air or AirLink2


SPARKlink Air is a sensor interface with two ports to accept PASPORT sensors and ports for the included Voltage Probe and Temperature Probe.

The SPARKlink Air interface connects to a computer or tablet device wirelessly via Bluetooth (radio), or can also connect to a computer's USB port with the included USB cable. The SPARKlink Air operates on battery power or can be connected to the included AC adapter. The SPARKlink Air battery can also be recharged using the USB cable connected to a USB port.

### Procedure

1. Optional: Connect the AC adapter (included with the SPARKlink Air) to a wall outlet and to the AC adapter port on the bottom of the device.

➡ The Battery Status LED may blink red to indicate that the battery is charging through the power cable.

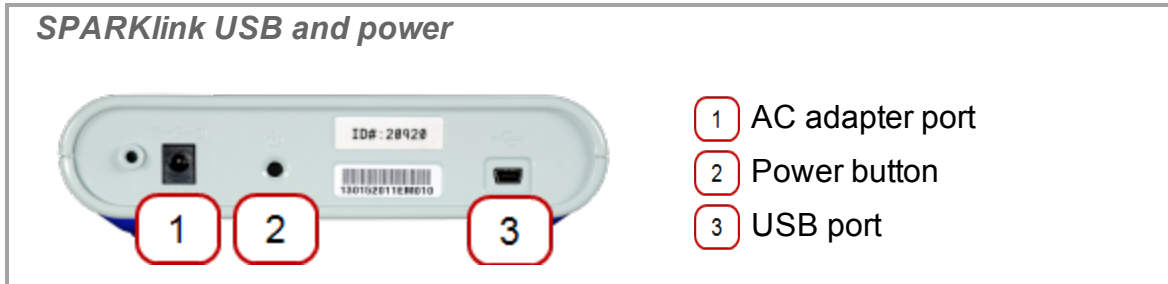

2. Connect the SPARKlink Air to the computer via bluetooth or USB.


#### Pairing bluetooth

- i. Press and hold the ON/OFF button on the SPARKlink Air.
- ii. After the status LEDs blink in sequence, release the ON/OFF button.
  - ➡ The Bluetooth status LED will blink blue.
- iii. Enable the Windows or Macintosh computer for Bluetooth communication.

#### USB connection

- i. Use an A-to-mini-B USB cable (such as PASCO part PS-2528) to connect the smaller USB port of the SPARK to a USB port on your computer (or a USB hub connected to the computer).
- ii. Press and hold the ON/OFF button on the SPARKlink Air.
- iii. After the status LEDs blink in sequence, release the ON/OFF button.
  - ➡ The USB Status LED will shine green. The Battery Status LED may blink red to indicate that the battery is charging through the USB.


**LED status definitions**



1	Bluetooth status	Slow Blinking – Not connected
1	Bluetooth status	Solid – Connected
2	USB Status	Fast Blinking – Sampling data
2	USB Status	Off – Not connected
2	USB Status	On – Connected
3	Battery status	Solid red – Low battery
3	Battery status	Blinking red – Charging
3	Battery status	Solid green – Fully charged

3. Plug in PASPORT sensors, temperature probe, or voltage leads.

**SPARKlink ports**

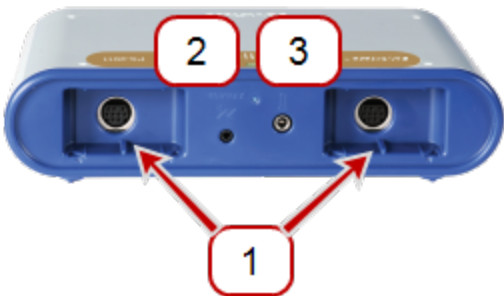

1	PASPORT ports
2	Voltage port
3	Temperature port

4. Configure the interface for Capstone. (See ["Connecting an interface": 17](#))

Next step: **"Connecting a sensor": 27**

### Connecting a 750 SCSI or 500 serial Interface

1. Use the A-to-mini-B USB cable (included with the 750 Interface) to connect the USB port of the 750 Interface to a USB port on your computer (or a USB hub connected to the computer).

2. Connect the AC adapter to a wall outlet and to the AC adapter port on the rear of the 750 Interface.
3. Turn on the interface and hold the power button.
4. Configure the interface for Capstone. ([See "Connecting an interface": 17](#))
5. Add sensors. ([See "Connecting a sensor with the 750 Interface": 41](#))


➔ **Next step: "Connecting a sensor": 27**

### Connecting USB PASPORT Interfaces: such as the SPARK SLS, SPARKlink, or Power Link

1. Use an A-to-mini-B USB cable (such as PASCO part PS-2528) to connect the smaller USB port of the SPARK SLS or SPARKlink to a USB port on your computer (or a USB hub connected to the computer).
2. Optionally connect the AC adapter (included with the SPARK SLS and SPARKlink) to a wall outlet and to the AC adapter port on the bottom of the device.



3. Plug in PASPORT sensors, temperature probe, or voltage leads.

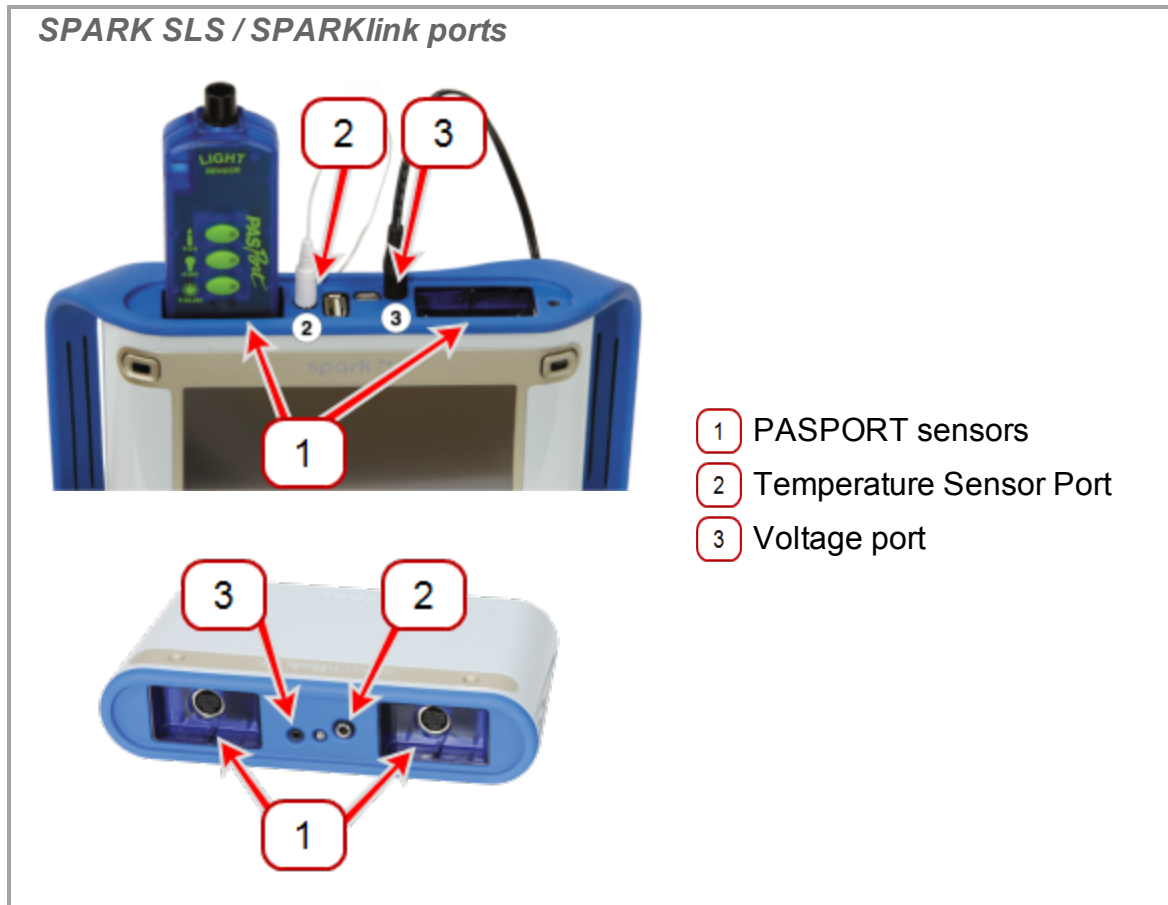

4. SPARK SLS: Press and hold the power button until the SPARK SLS screen turns on.
5. Configure the interface for Capstone. ([🔗 See "Connecting an interface": 17](#))
6. SPARK SLS: The SPARK SLS states it is connected to a computer. Disconnect the SPARK SLS from the computer to use it in standalone mode.

🔗 **Next step: "Connecting a sensor": 27**

## Connecting an Xplorer GLX or Xplorer

1. Use an A-to-mini B USB cable (included) to connect the smaller USB port of the Xplorer GLX to a USB port on your computer (or a USB hub connected to the computer).
2. Optionally, connect the included AC adapter to a wall outlet and to the AC adapter port on the right side of the Xplorer GLX.
3. If the Xplorer GLX is not on, press and hold the power button until the green LED

on the front flashes.

4. Configure the interface for Capstone. (📎 See "Connecting an interface": 17)

➔ Next step: **"Connecting a sensor": 27**

## Connecting a sensor

Capstone software running on your computer receives data from one or more connected sensors through one or more interfaces. (At least one interface and sensor is required to collect data. (📎 See "What do I need to have a complete measurement system? ": 162)

Follow these steps to connect a sensor to your computer.

1. If not already, connect or add an interface to your computer. (📎 See "Connecting an interface": 17)

**Note:** Capstone requires an interface to communicate with the sensor.

2. Plug in one or more sensors into the interface.

### Click for instructions on select interfaces



"Connecting a sensor with the 850 Interface": 36



"Connecting a sensor with the 750 Interface": 41



"Connecting a sensor with the SPARK SLS, SPARKlink, SPARKlink Air": 45



"Connecting a sensor with the Xplorer GLX": 50




"Connecting a sensor with the Digital Adapter or Photogate Port": 55



"Connecting a sensor with the Analog Adapter": 56



"Connecting a sensor with the AirLink2": 56

3. Add the sensor in Capstone  **Tip<sup>1</sup>**


### PASPORT sensor

<sup>1</sup>You can add sensors that are not connected to the interface. This is useful to author or prepare experiments in advance.




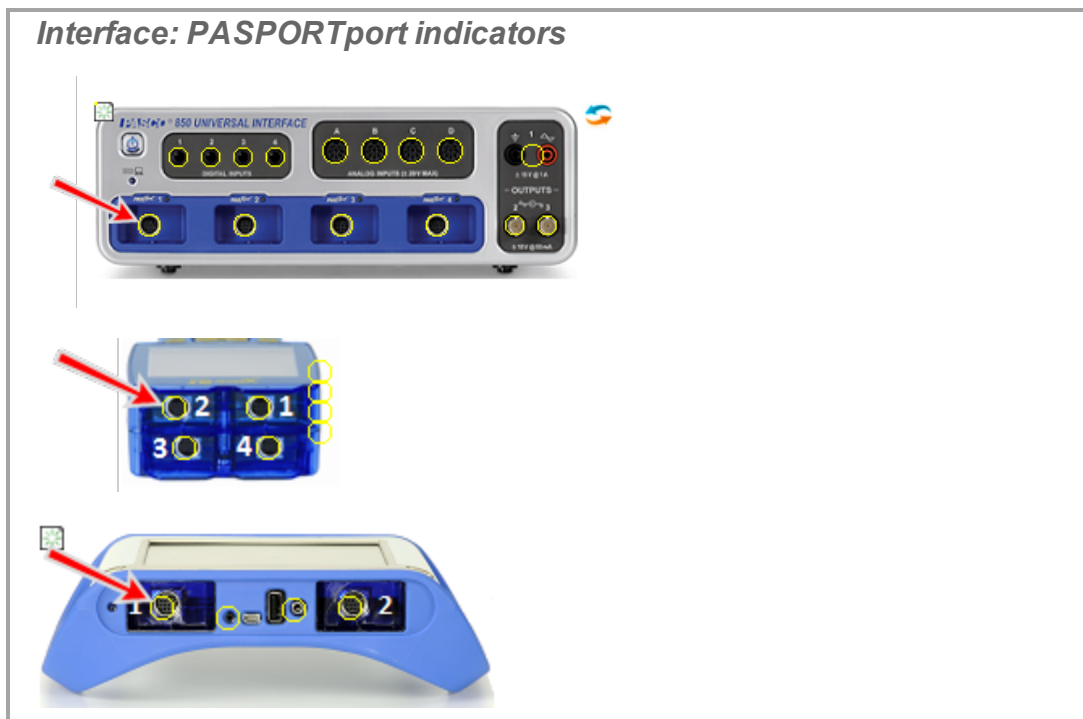
**Note:** Capstone automatically detects properly connected PASPORT sensors.

*Automatically detected PASPORT sensors*

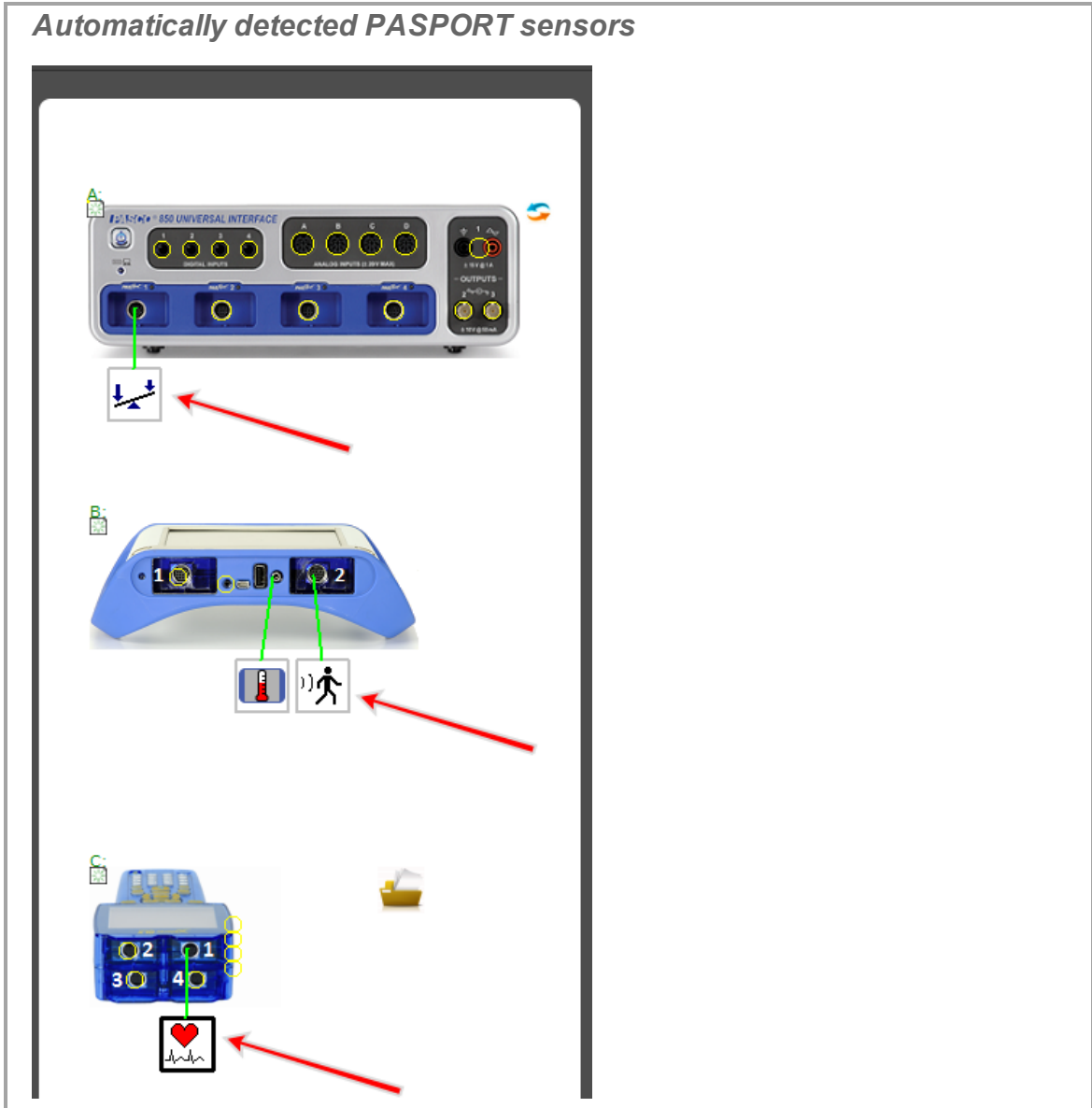


To manually add a PASPORT sensor:

- a. Click within the yellow circle of the port to add a sensor. 



- b. Select the sensor to add from the drop down list.


➔ The sensor icon appears for the port.

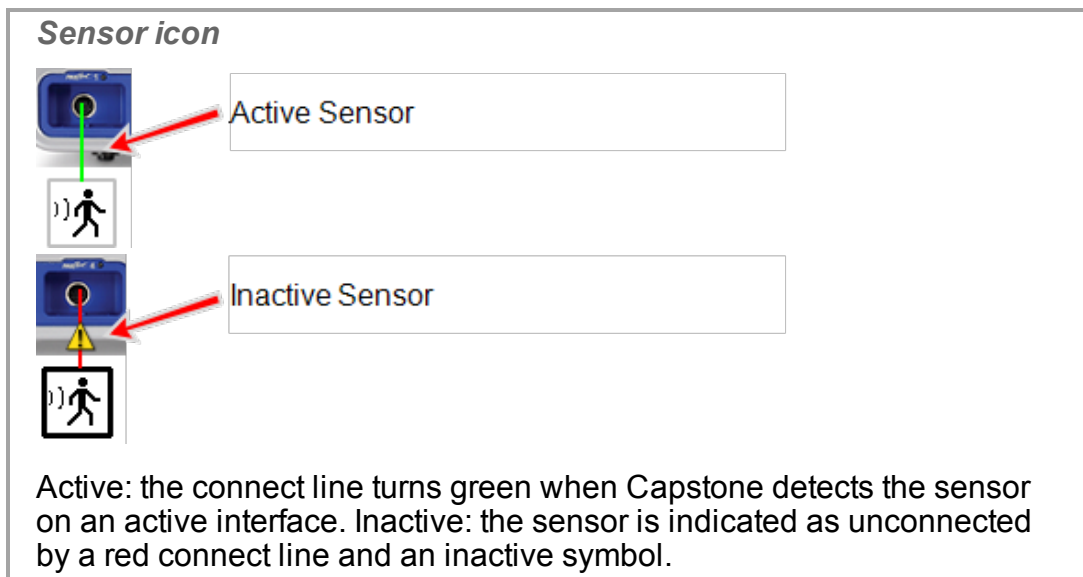


### Analog ScienceWorkshop sensor (Universal Interface)


**Note:** Capstone automatically detects Universal Interface (UI) sensors.

To add a *non-connected* sensor to Capstone:

- a. Click within the yellow circle of the port to add a sensor. 

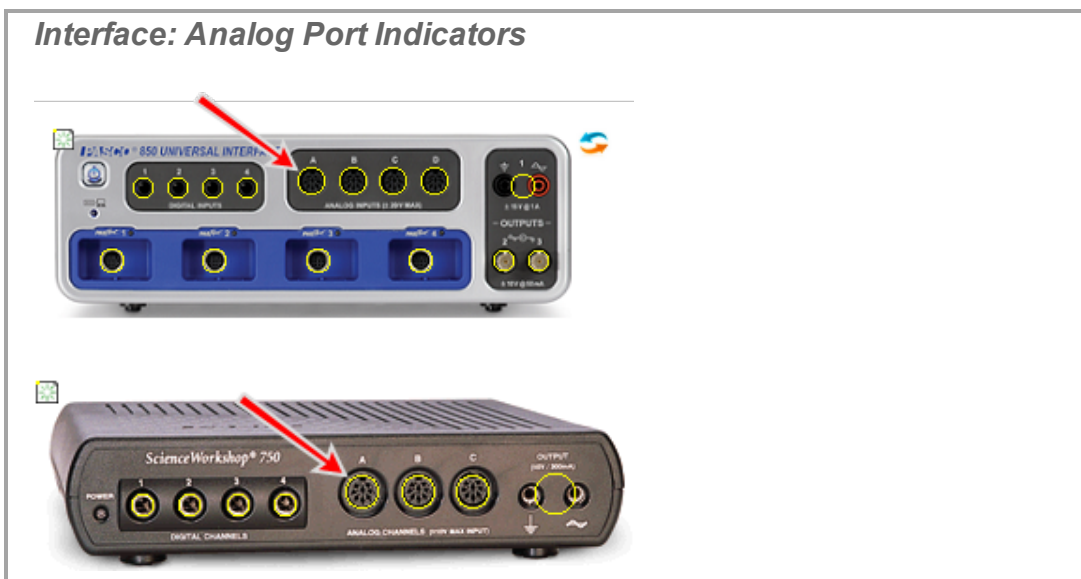


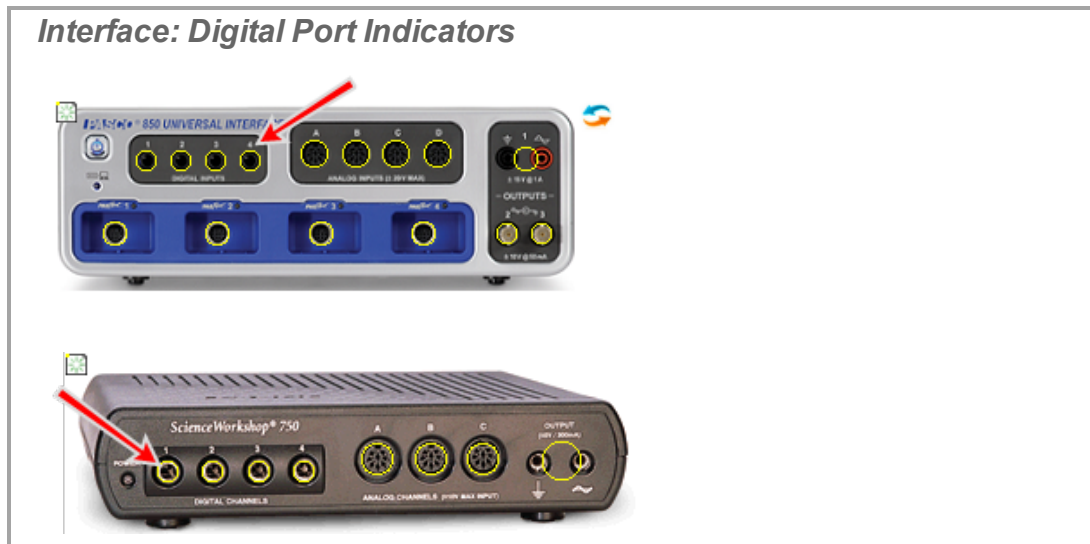
- b. Select the sensor to add from the drop down list.
  - ➡ The sensor icon appears for the port.

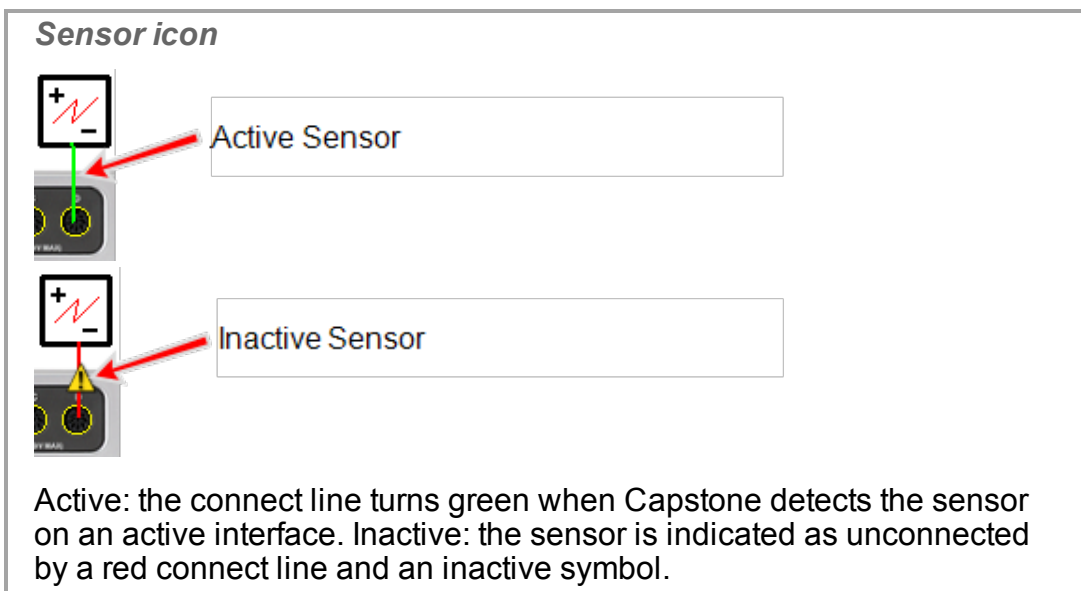

### Digital sensor (ScienceWorkshop)


**Note:** Capstone does not automatically detect most digital sensors.


- a. Click within the yellow circle of the port to add a sensor. 




- b. Select the sensor to add from the drop down list.

➔ The sensor icon appears for the port.

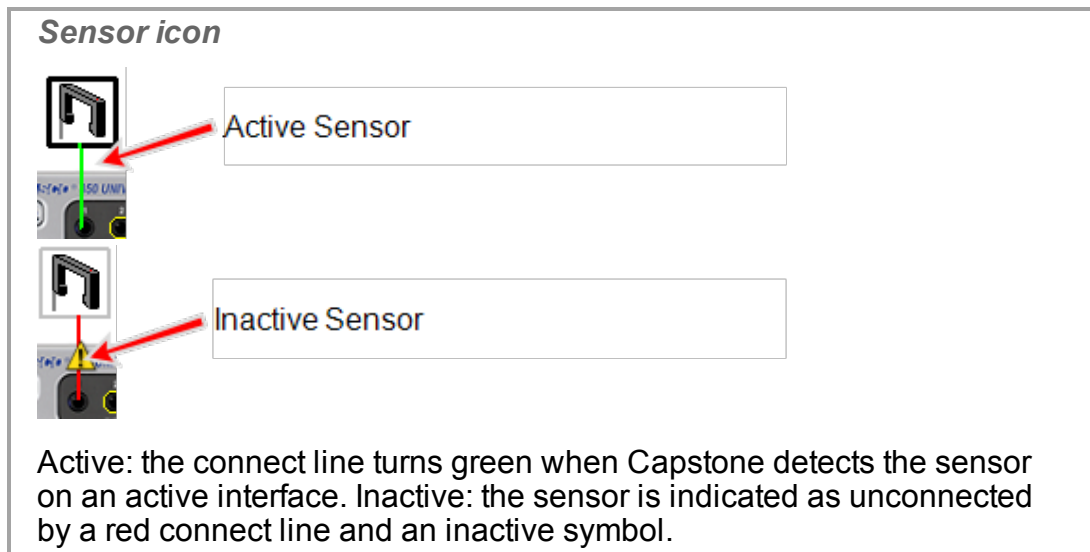


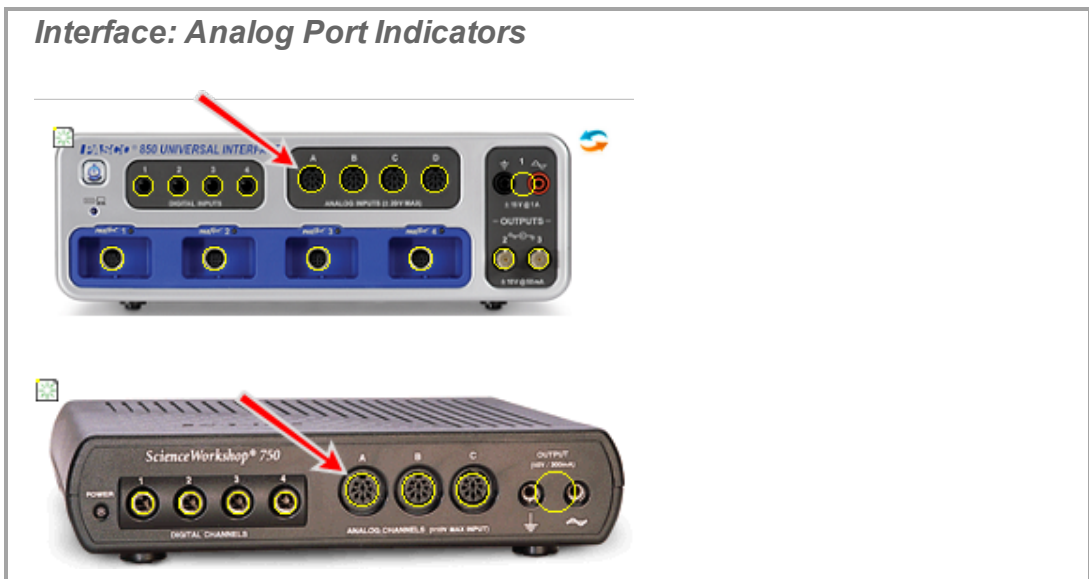
- c. Optional: to use the sensor as a timer, click **Timer Setup**  in the Tools Palette.

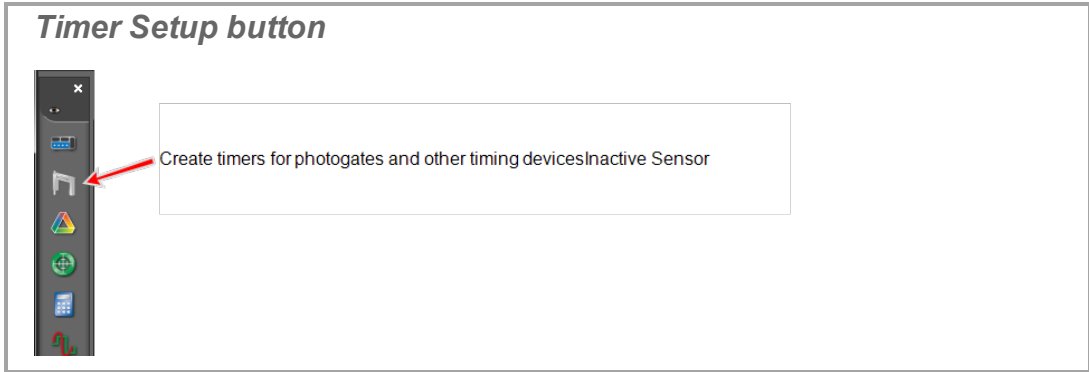

### Analog sensor (ScienceWorkshop)


**Note:** Capstone does not automatically detect most analog ScienceWorkshop sensors.

- a. Click within the yellow circle of the port to add a sensor. 



- b. Select the sensor to add from the drop down list.  
 ➔ The sensor icon appears for the port.

**Sensor icon**

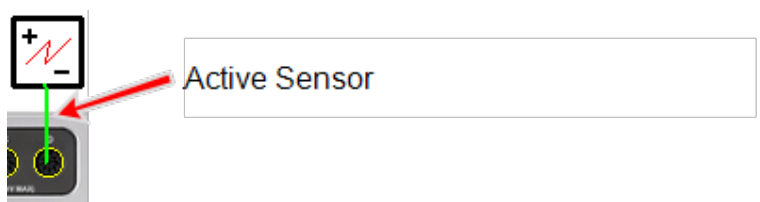

Active Sensor

Analog sensors do not toggle between inactive and active and display as active (green line) even when disconnected.


**Temperature sensor (SPARK SLS, SPARKlink, SPARKlink Air, and Xplorer GLX)**

**Note:** Capstone automatically detects temperature sensors connected to the internal temperature sensor port.

To manually add a temperature sensor:

- a. Click within the yellow circle of the temperature port to add a temperature sensor .


**Interface: Temperature port indicators**

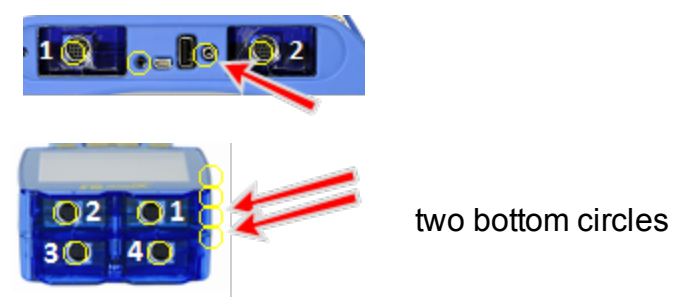

two bottom circles

- ➡ The sensor icon appears for the port.

**Sensor icon**

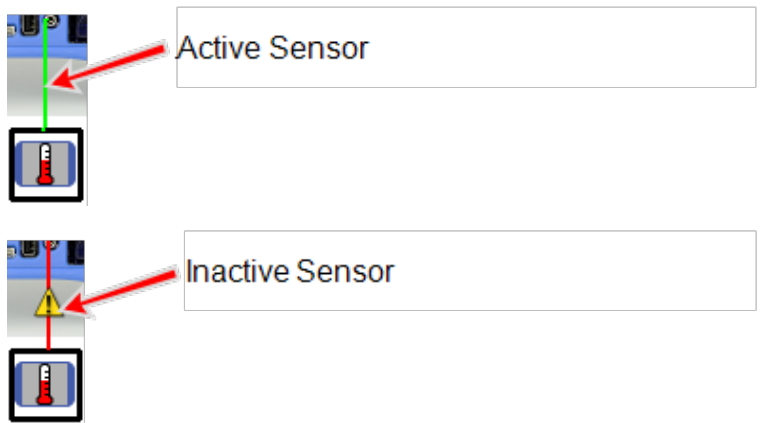

Active Sensor

Inactive Sensor

Active: the connect line turns green when Capstone detects the sensor on an active interface. Inactive: the sensor is indicated as unconnected by a red connect line and an inactive symbol.


### Voltage sensor (SPARK SLS, SPARKlink, SPARKlink Air, and Xplorer GLX)

**Note:** Capstone automatically detects sound sensors connected to the internal sound sensor port.

To manually add a voltage sensor:

- a. Click within the yellow circle of the port to add a voltage sensor.


**Interface: Voltage port indicators**

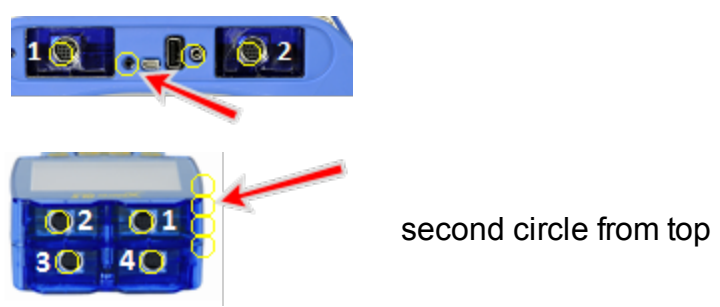

second circle from top

- ➔ The sensor icon appears for the port.

**Sensor icon**

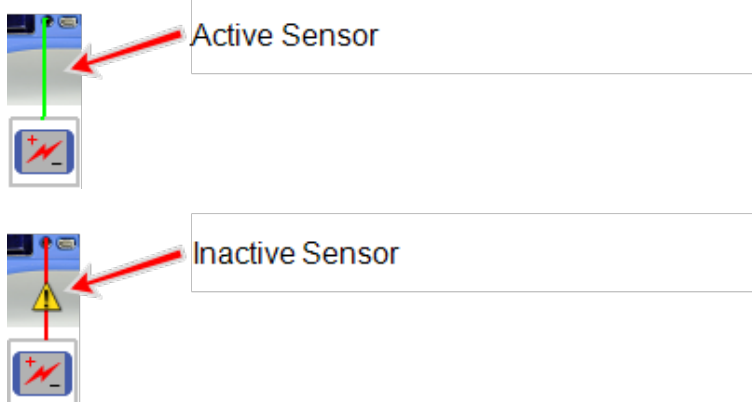

Active: the connect line turns green when Capstone detects the sensor on an active interface. Inactive: the sensor is indicated as unconnected by a red connect line and an inactive symbol.


### Sound sensor (Xplorer GLX)

**Note:** Capstone automatically detects voltage sensors connected to the internal voltage sensor port.

To manually add a sound sensor:

- a. Click within the yellow circle of the port to add a sound sensor .

**Interface: Sound port indicators**

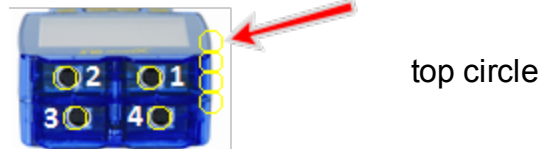

top circle

- b. Select the **GLX Sound Level Sensor** or the **GLX Sound Sensor**.
  - ➔ The sensor icon appears for the port.



**Sensor icon**

Active Sensor

Inactive Sensor

Active: the connect line turns green when Capstone detects the sensor on an active interface. Inactive: the sensor is indicated as unconnected by a red connect line and an inactive symbol.

➔ Next step: **"Adding a display ": 56**

### Connecting a sensor with the 850 Interface

1. Plug in PASPORT sensors, digital ScienceWorkshop sensors, or analog ScienceWorkshop (black) sensors.

**850 Universal Interface ports**


- 1 Digital ScienceWorkshop – 4 ports
- 2 Analog ScienceWorkshop – 4 ports
- 3 PASPORT – 4 ports **Tip<sup>1</sup>**

2. Add the sensor in Capstone:


#### PASPORT sensor


**Note:** Capstone automatically detects properly connected PASPORT sensors.

<sup>1</sup>PASPORT sensors are automatically detected.

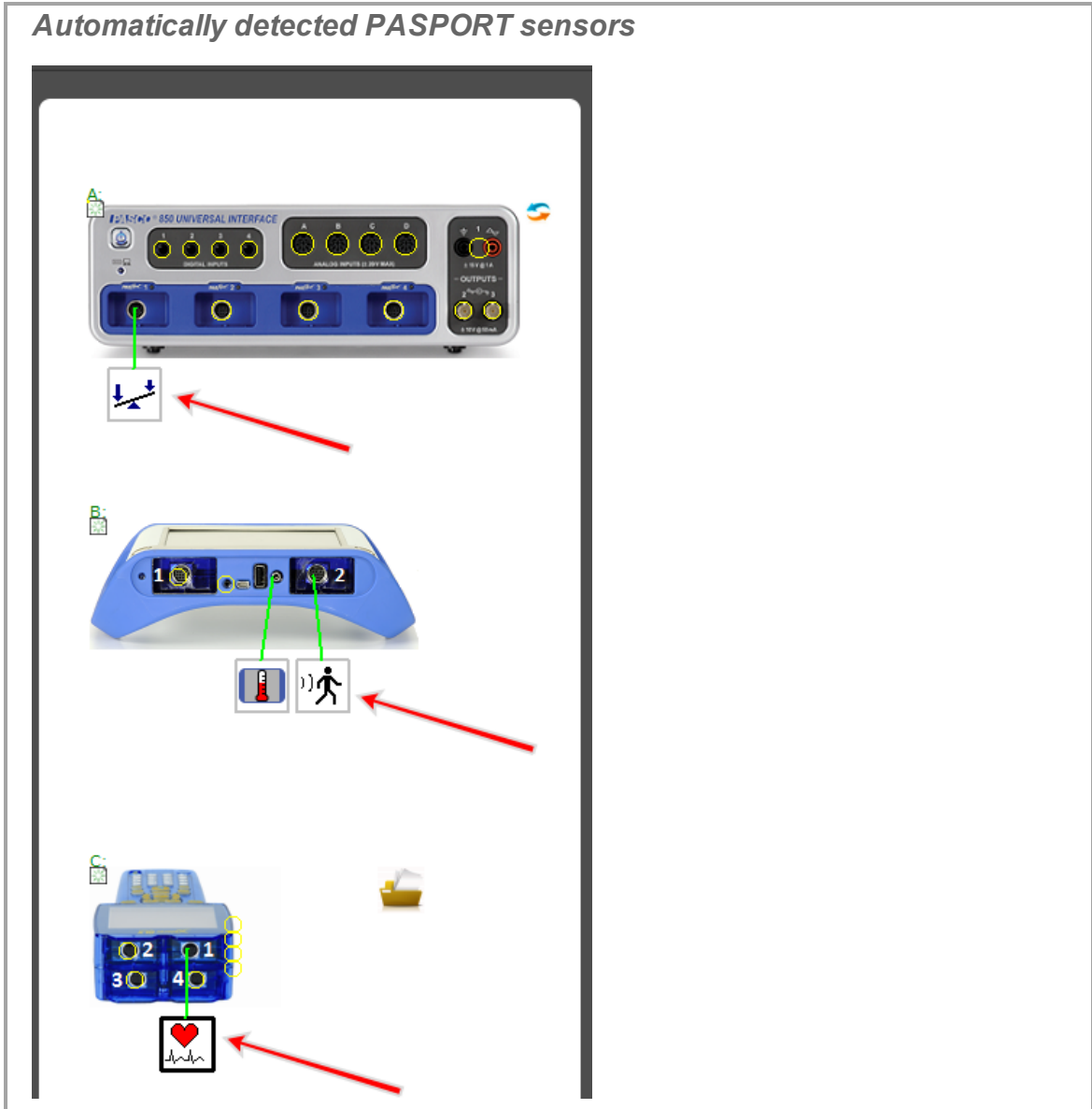


To manually add a PASPORT sensor:

- a. Click within the yellow circle of the port to add a sensor. 

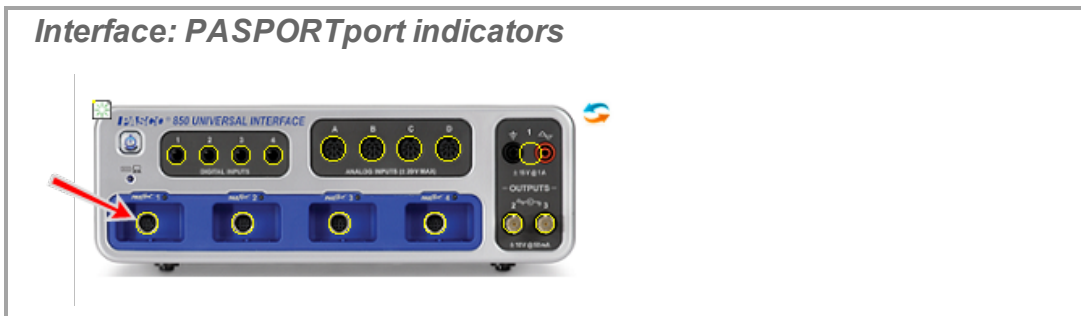

- b. Select the sensor to add from the drop down list.

➔ The sensor icon appears for the port.

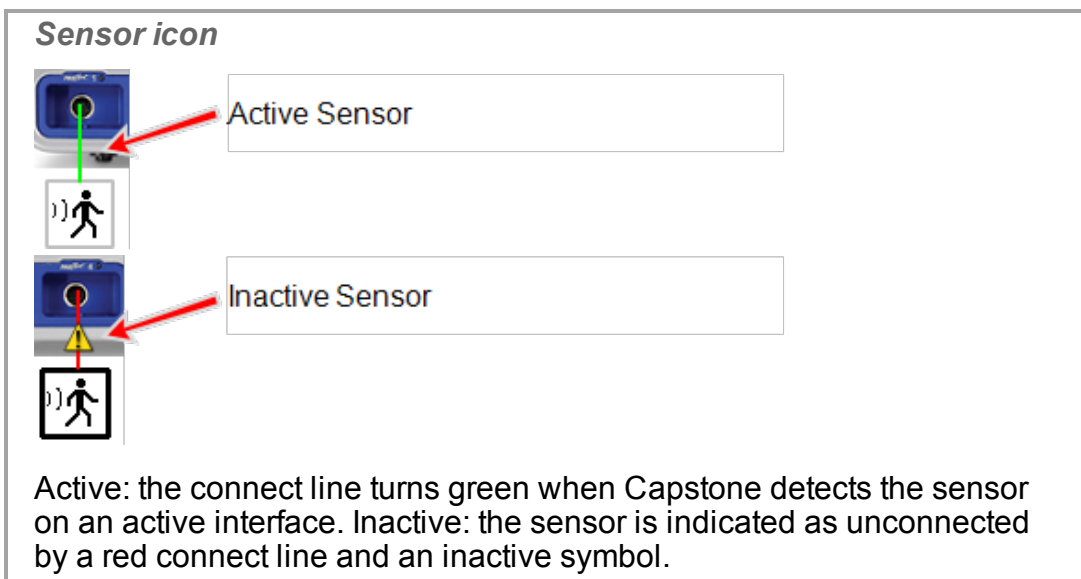

### Digital sensor (ScienceWorkshop)


**Note:** Capstone does not automatically detect most digital sensors.


- a. Click within the yellow circle of the port to add a sensor.




- b. Select the sensor to add from the drop down list.

➔ The sensor icon appears for the port.

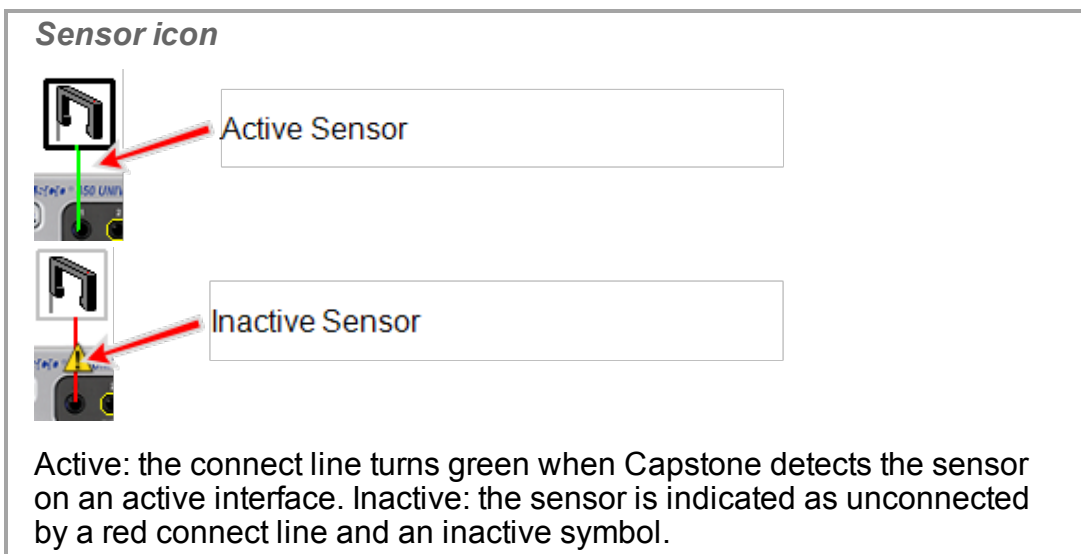


- c. Optional: to use the sensor as a timer, click **Timer Setup**  in the Tools Palette.

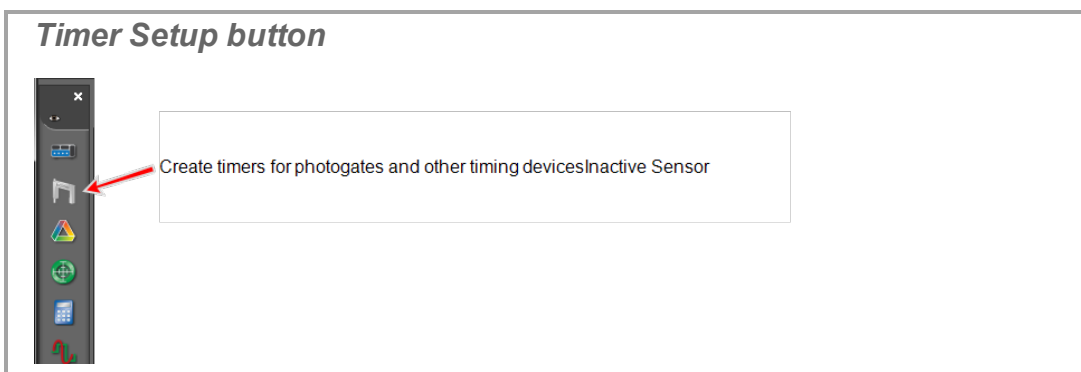


## Analog sensor (ScienceWorkshop)

**Note:** Capstone does not automatically detect most analog ScienceWorkshop sensors.

- a. Click within the yellow circle of the port to add a sensor. 



- b. Select the sensor to add from the drop down list.  
 ➔ The sensor icon appears for the port.

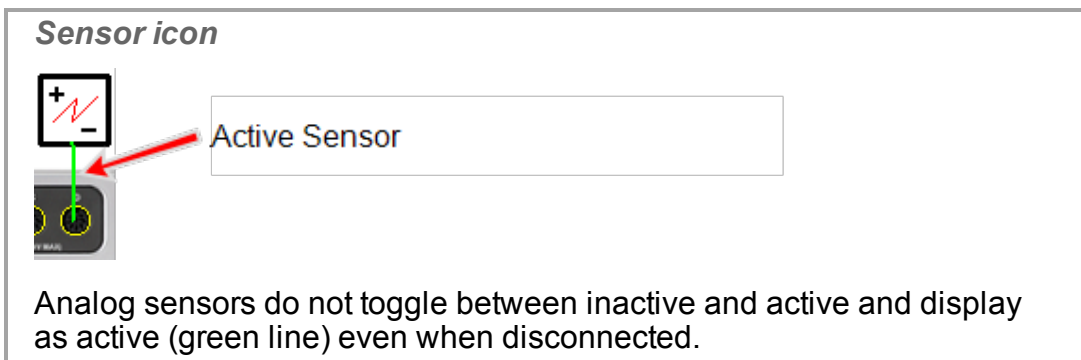


## Analog ScienceWorkshop sensor (Universal Interface)

**Note:** Capstone automatically detects Universal Interface (UI) sensors.


To add a *non-connected* sensor to Capstone:

- a. Click within the yellow circle of the port to add a sensor.

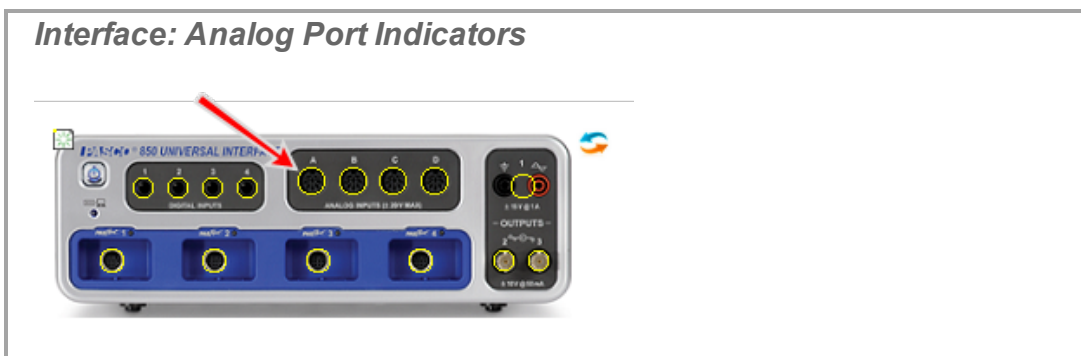

- b. Select the sensor to add from the drop down list.

➡ The sensor icon appears for the port.

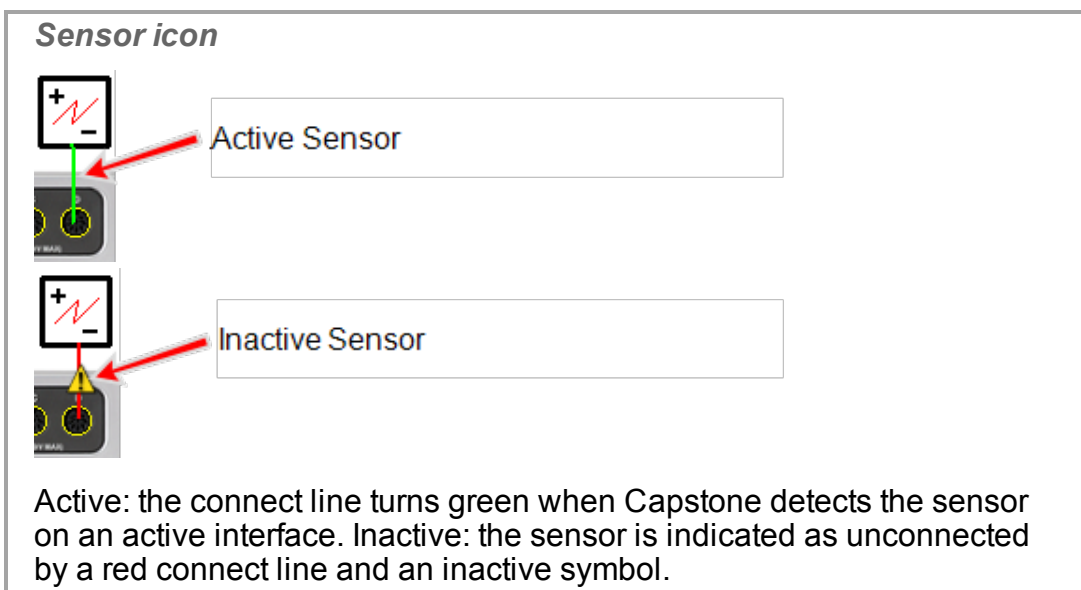

➡ Next step: **"Adding a display ": 56**

### Connecting a sensor with the 750 Interface

1. Plug digital ScienceWorkshop sensors or analog ScienceWorkshop (black) sensors.

750 Interface ports



2. Add the sensor in Capstone:

Digital sensor (ScienceWorkshop)

**Note:** Capstone does not automatically detect most digital sensors.

a. Click within the yellow circle of the port to add a sensor.


Interface: Digital Port Indicators




b. Select the sensor to add from the drop down list.


➡ The sensor icon appears for the port.

Sensor icon

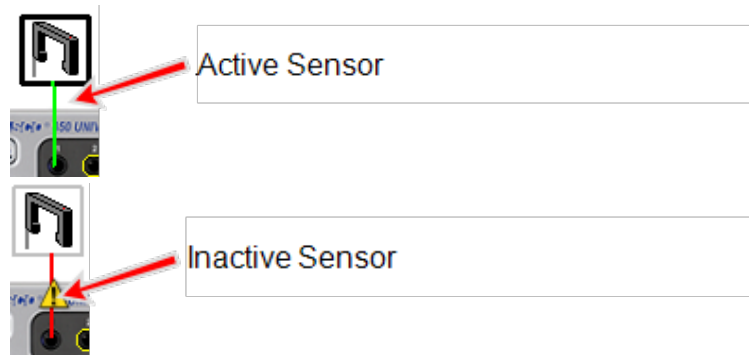



Active: the connect line turns green when Capstone detects the sensor on an active interface. Inactive: the sensor is indicated as unconnected by a red connect line and an inactive symbol.

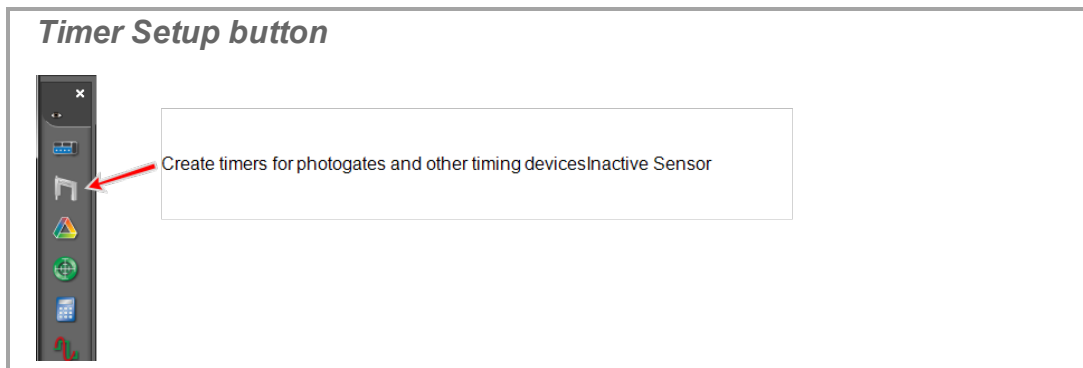- c. Optional: to use the sensor as a timer, click **Timer Setup**  in the Tools Palette.




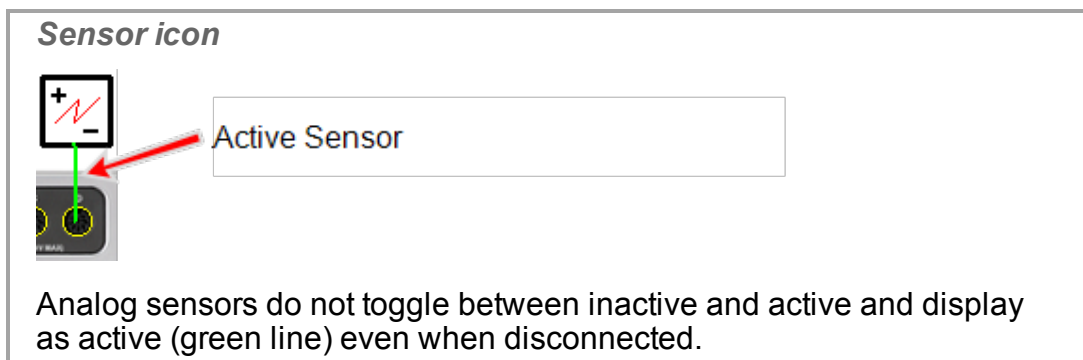
### Analog sensor (ScienceWorkshop)

**Note:** Capstone does not automatically detect most analog ScienceWorkshop sensors.

- a. Click within the yellow circle of the port to add a sensor.  



- b. Select the sensor to add from the drop down list.  
 The sensor icon appears for the port.




### Analog ScienceWorkshop sensor (Universal Interface)




**Note:** Capstone automatically detects Universal Interface (UI) sensors.

To add a *non-connected* sensor to Capstone:

- a. Click within the yellow circle of the port to add a sensor. 




- b. Select the sensor to add from the drop down list.
  - ➡ The sensor icon appears for the port.

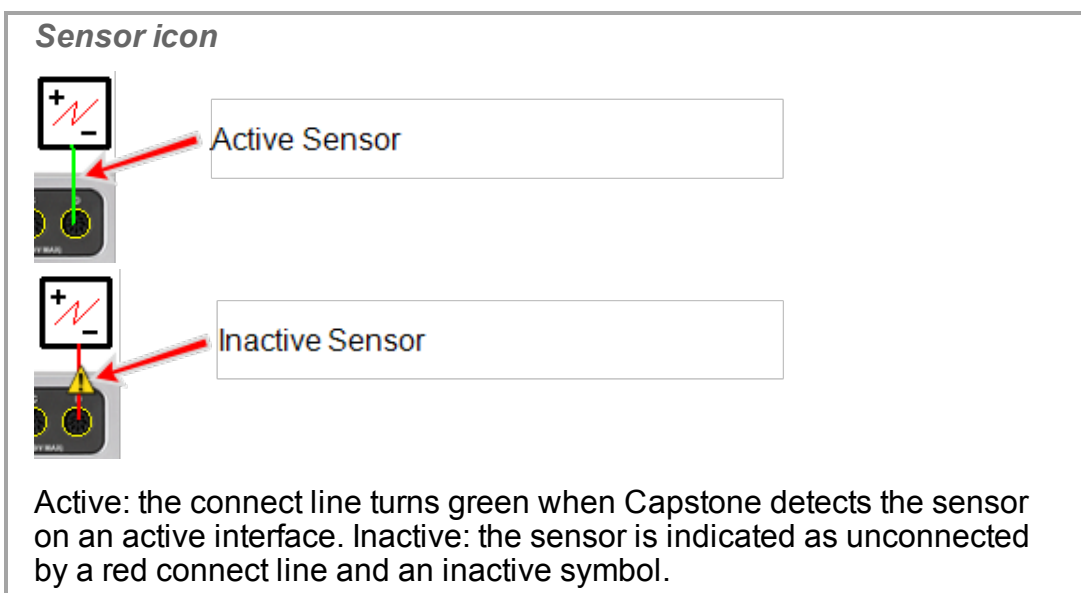

➡ Next step: **"Adding a display ": 56**

## Connecting a sensor with the SPARK SLS, SPARKlink, SPARKlink Air


1. Plug in PASPORT sensors, temperature probe, and/or voltage leads.

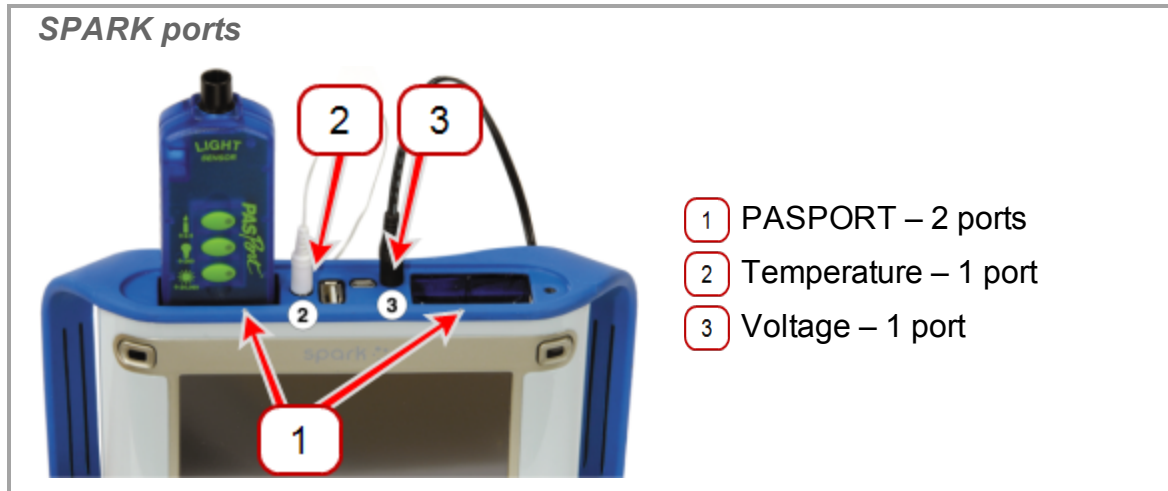

2. Press and hold the power button.
3. Configure the interface for Capstone. ([See "Connecting an interface": 17](#))
4. Add the sensor in Capstone:


### PASPORT sensor

**Note:** Capstone automatically detects properly connected PASPORT sensors.

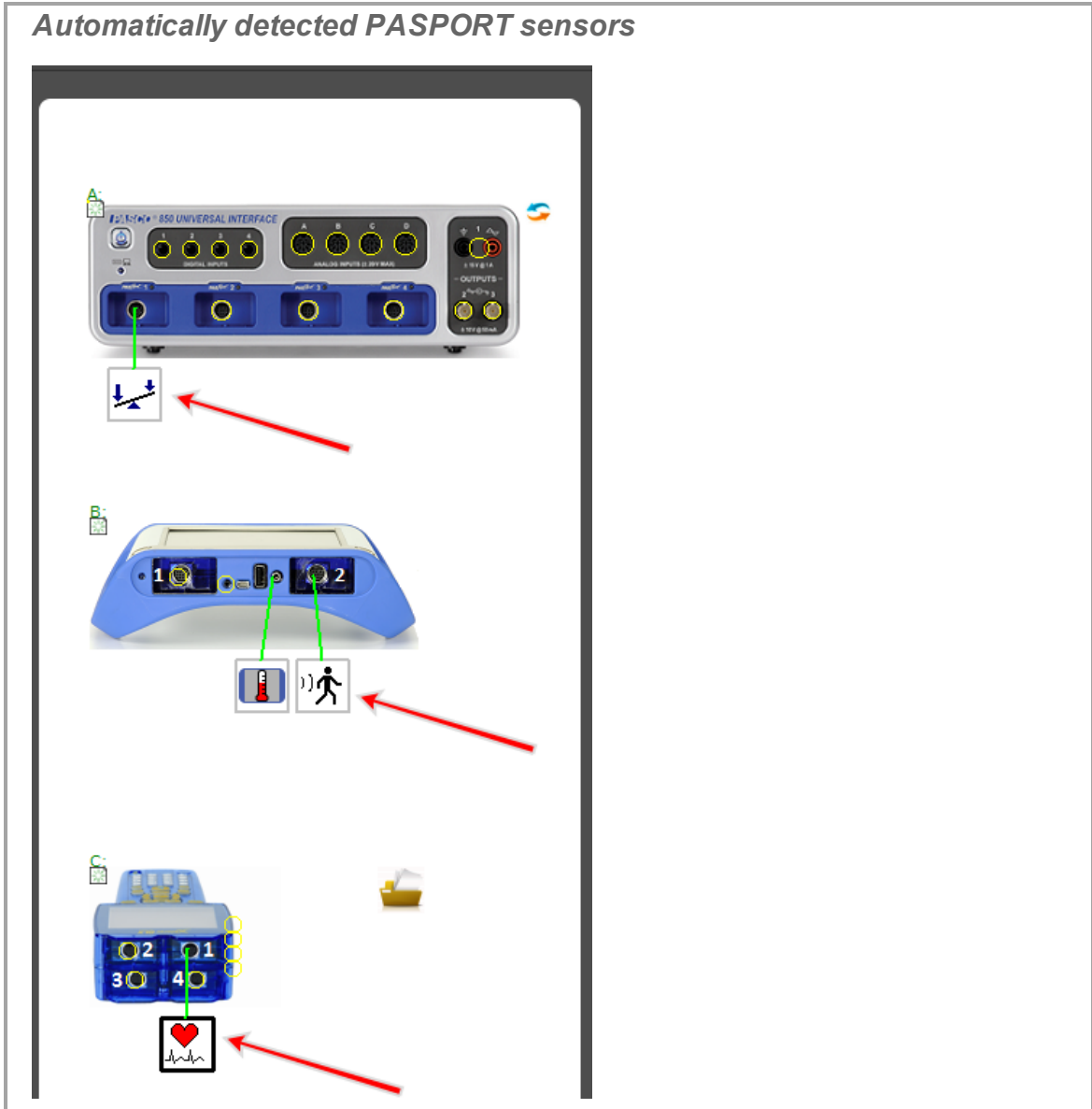


To manually add a PASPORT sensor:

- a. Click within the yellow circle of the port to add a sensor.

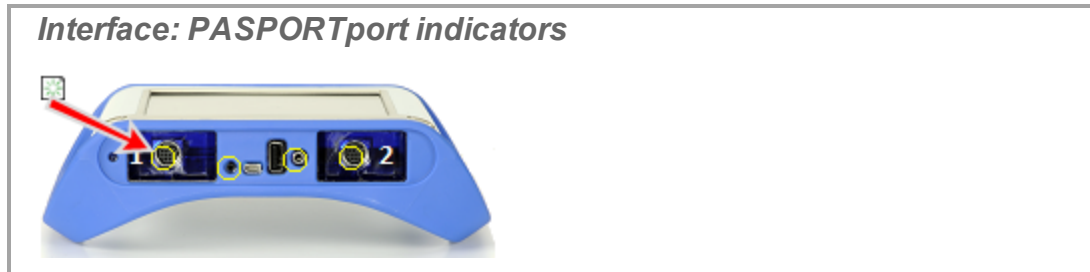

- b. Select the sensor to add from the drop down list.

➡ The sensor icon appears for the port.

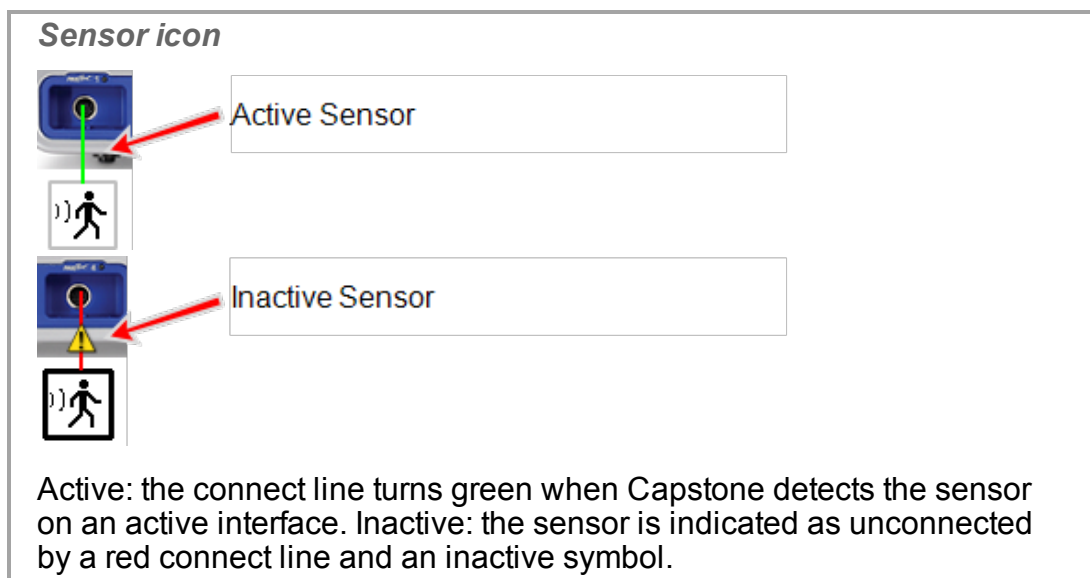


## Temperature sensor

**Note:** Capstone automatically detects temperature sensors connected to the internal temperature sensor port.


To manually add a temperature sensor:

- a. Click within the yellow circle of the temperature port to add a temperature sensor .

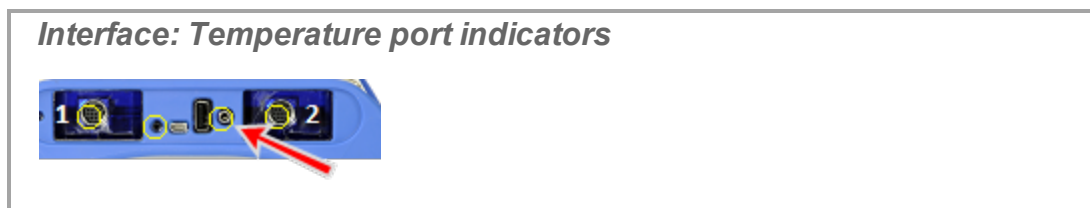

➔ The sensor icon appears for the port.

**Sensor icon**

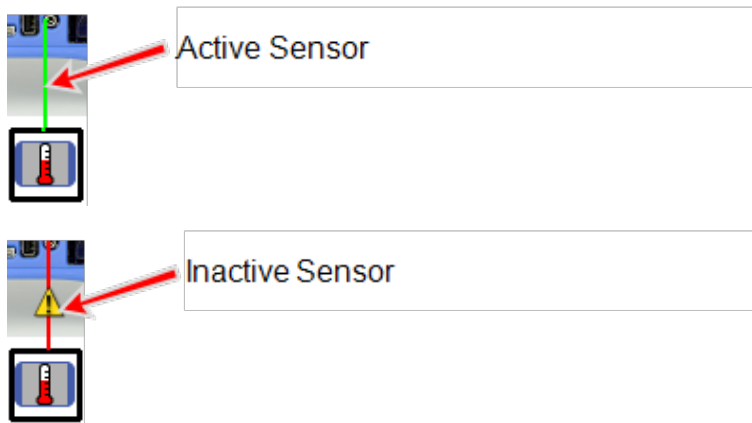

Active: the connect line turns green when Capstone detects the sensor on an active interface. Inactive: the sensor is indicated as unconnected by a red connect line and an inactive symbol.


### Voltage sensor

**Note:** Capstone automatically detects sound sensors connected to the internal sound sensor port.

To manually add a voltage sensor:


- a. Click within the yellow circle of the port to add a voltage sensor.

**Interface: Voltage port indicators**



➔ The sensor icon appears for the port.

*Sensor icon*

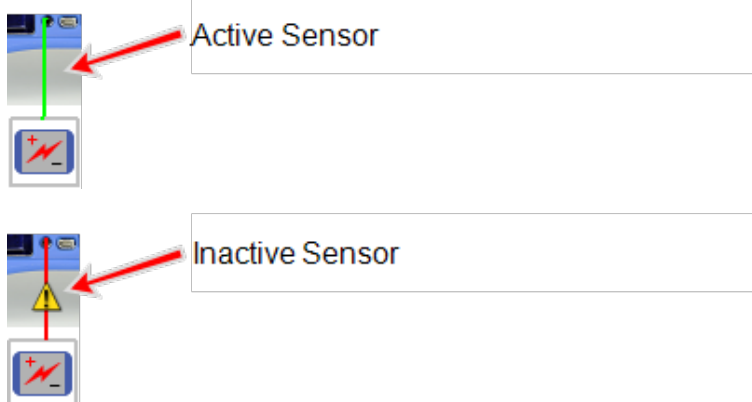

Active Sensor


Inactive Sensor

Active: the connect line turns green when Capstone detects the sensor on an active interface. Inactive: the sensor is indicated as unconnected by a red connect line and an inactive symbol.

➤ Next step: **"Adding a display ": 56**

## Connecting a sensor with the Xplorer GLX


1. Plug in PASPORT sensors, temperature probe, or voltage leads.

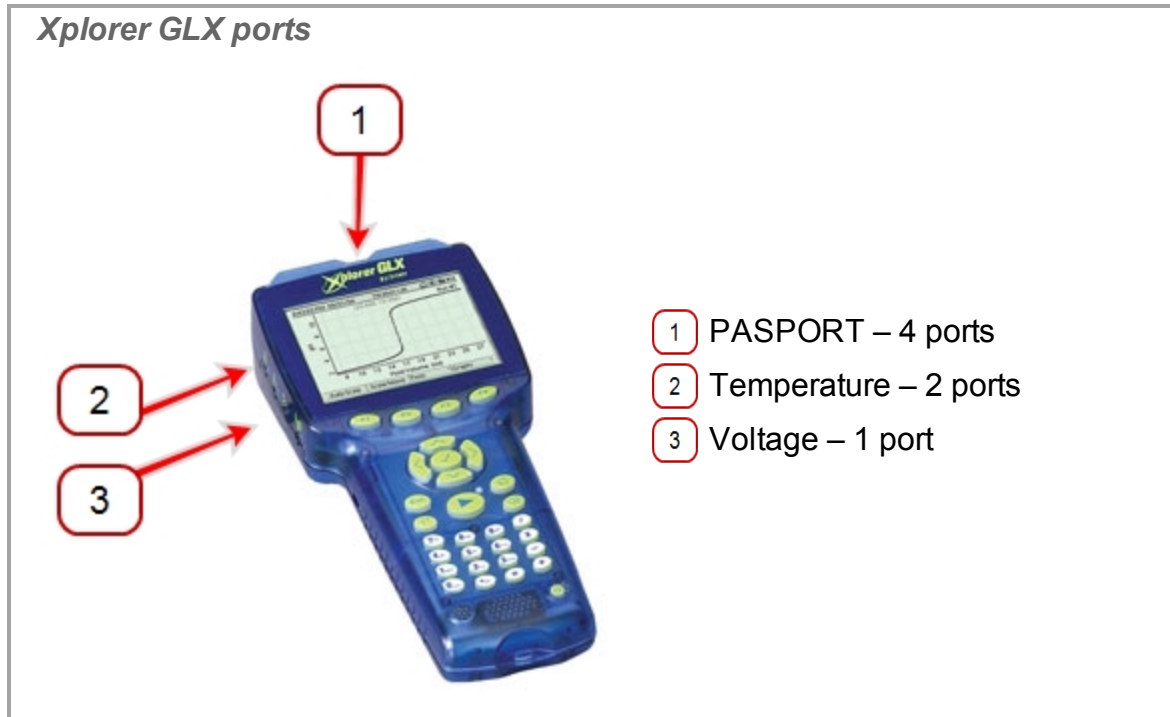

3. Press and hold the power button.
4. Configure the interface for Capstone. ([📎 See "Connecting an interface": 17](#))
5. Add the sensor in Capstone:

### PASPORT sensor

**Note:** Capstone automatically detects properly connected PASPORT sensors.

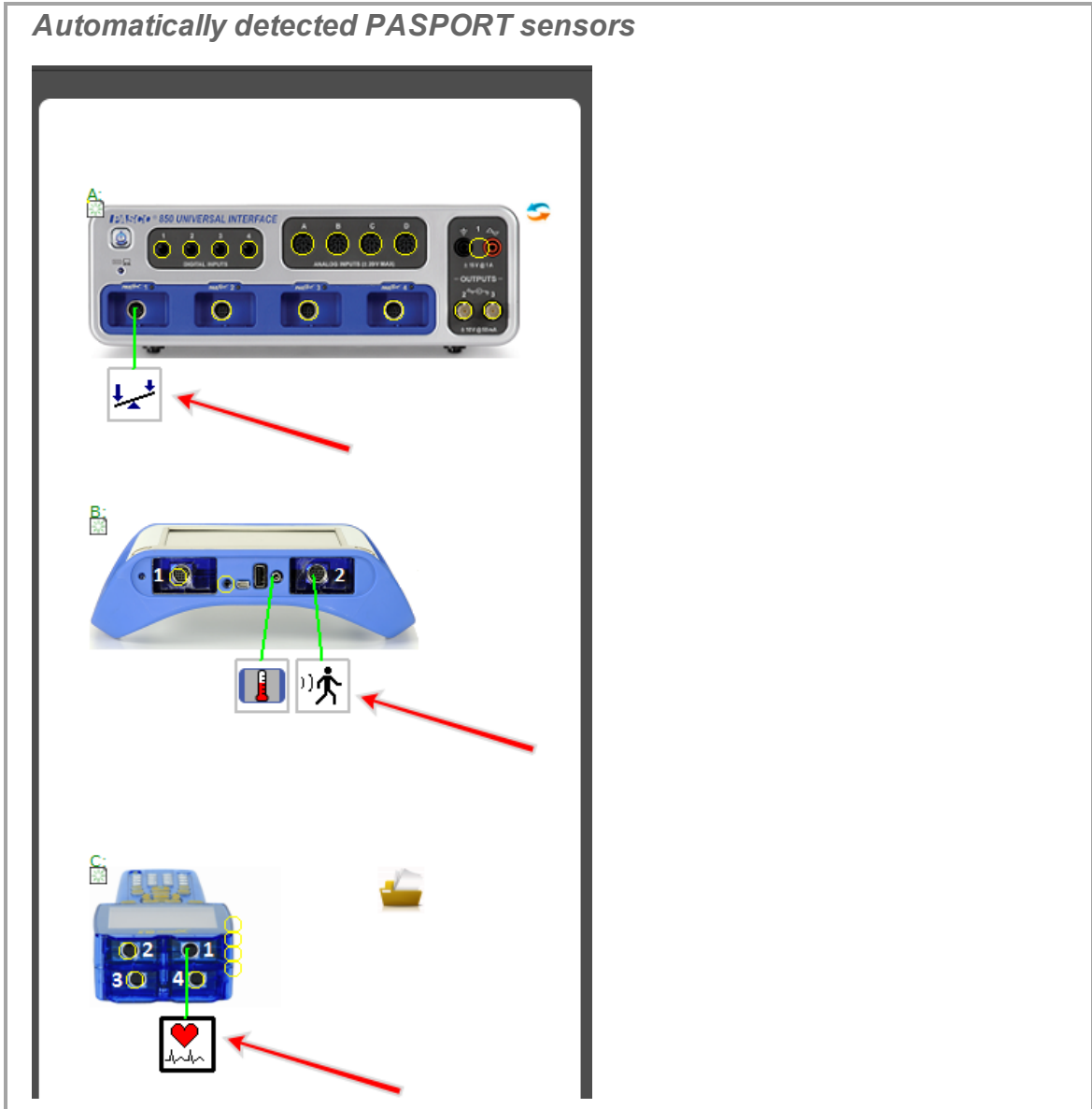

To manually add a PASPORT sensor:



- a. Click within the yellow circle of the port to add a sensor.




*Interface: PASPORTport indicators*



- b. Select the sensor to add from the drop down list.

➔ The sensor icon appears for the port.

*Sensor icon*

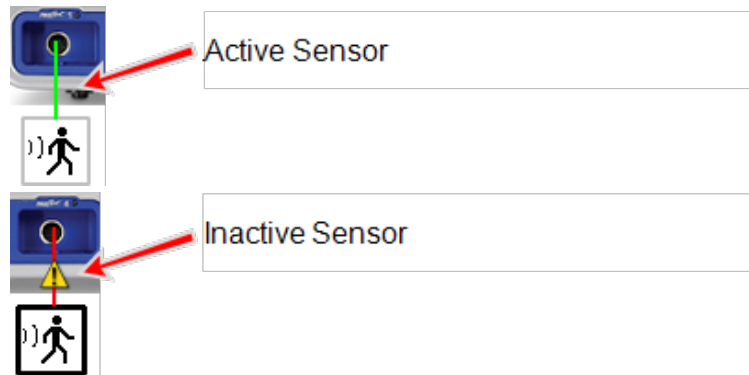

Active: the connect line turns green when Capstone detects the sensor on an active interface. Inactive: the sensor is indicated as unconnected by a red connect line and an inactive symbol.

## Temperature sensor

**Note:** Capstone automatically detects temperature sensors connected to the internal temperature sensor port.

To manually add a temperature sensor:


- a. Click within the yellow circle of the temperature port to add a temperature sensor .

*Interface: Temperature port indicators*



➔ The sensor icon appears for the port.

**Sensor icon**

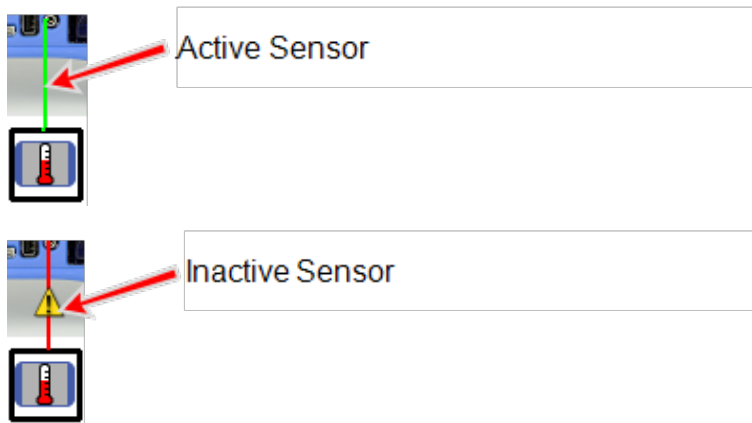

Active: the connect line turns green when Capstone detects the sensor on an active interface. Inactive: the sensor is indicated as unconnected by a red connect line and an inactive symbol.


### Voltage sensor

**Note:** Capstone automatically detects sound sensors connected to the internal sound sensor port.

To manually add a voltage sensor:

- a. Click within the yellow circle of the port to add a voltage sensor.


**Interface: Voltage port indicators**

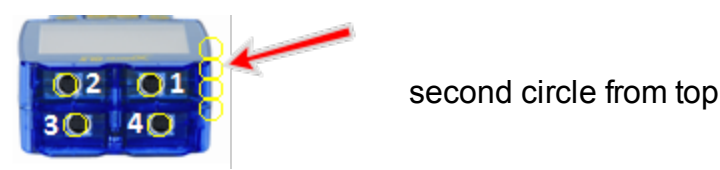

second circle from top

➔ The sensor icon appears for the port.

**Sensor icon**

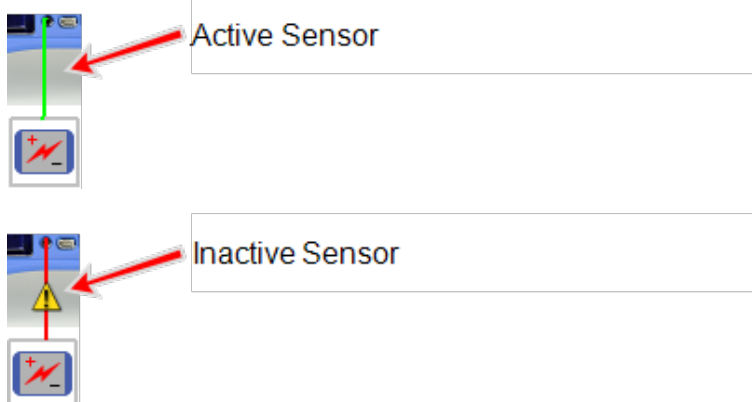

Active Sensor

Inactive Sensor

Active: the connect line turns green when Capstone detects the sensor on an active interface. Inactive: the sensor is indicated as unconnected by a red connect line and an inactive symbol.


## Sound sensor

**Note:** Capstone automatically detects voltage sensors connected to the internal voltage sensor port.

To manually add a sound sensor:


- a. Click within the yellow circle of the port to add a sound sensor .

**Interface: Sound port indicators**



top circle

- b. Select the **GLX Sound Level Sensor** or the **GLX Sound Sensor**.
  - ➔ The sensor icon appears for the port.

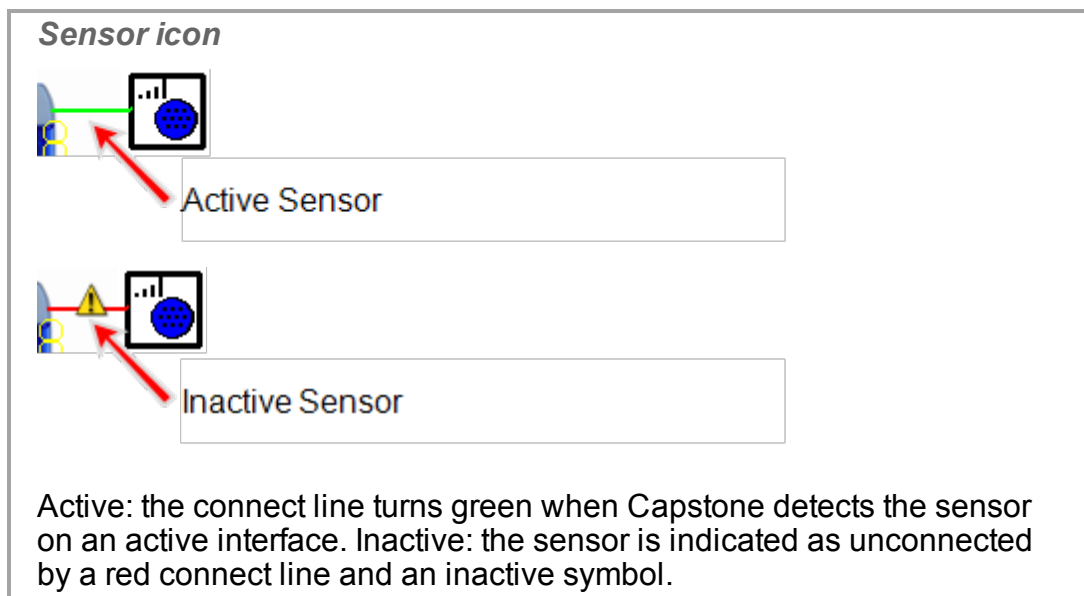

➤ Next step: **"Adding a display ": 56**

## Connecting a sensor with the Digital Adapter or Photogate Port

The Digital Adapter (PASCO part PS-2159) allows any ScienceWorkshop digital sensor sensors to be connected to a PASPORT interface to be used with Capstone. It also allows the use of the ScienceWorkshop Motion Sensor (CI-6742A) and the Rotary Motion Sensor (CI-6538).

The Photogate Port (PS-2123) is an obsolete adapter that supports switch-type sensors, but not the Rotary Motion Sensor or Motion Sensor.

### Procedure

1. Connect the adapter to your PASPORT interface.
2. Connect a digital sensor to the adapter.
  - ➡ You can connect a switch-type sensor to either port of the adapter.
  - ➡ If you are connecting a Motion Sensor or Rotary Motion Sensor, connect the yellow plug to Port 1 and the black plug to Port 2.
3. A list of digital sensors and sensor configurations appears.
4. Optionally, connect a second switch-type sensor to the other port of the adapter.
5. In the list, click one of the sensors or configurations to select it.

➤ Next step: **"Adding a display ": 56**

## Connecting a sensor with the Analog Adapter

The Analog Adapter (PASCO part PS-2158) allows analog ScienceWorkshop sensors to be connected to a PASPORT interface to be used with Capstone.

### Procedure

1. Connect the Analog Adapter to your PASPORT interface.
2. Connect an analog sensor to the adapter.
  - ➡ A list of sensors appears.
3. Click one of the sensors to select it.
4. Optionally, click the **Gain:** box and click a gain setting.
5. Click **OK**.

➡ Next step: "[Adding a display](#)": 56

## Connecting a sensor with the AirLink2

The AirLink2 sensor interface can communicate with computers using the USB connection or the via the bluetooth interface.

➡ Next step: "[Adding a display](#)": 56

## Adding a display


Capstone provides a highly flexible environment for workbook page layout.

Pages can be created using any combination of templates, drag and drop to targets, and moving and resizing display elements.



### Use a QuickStart template

1. Click one of the provided templates.  **Tip**<sup>1</sup>

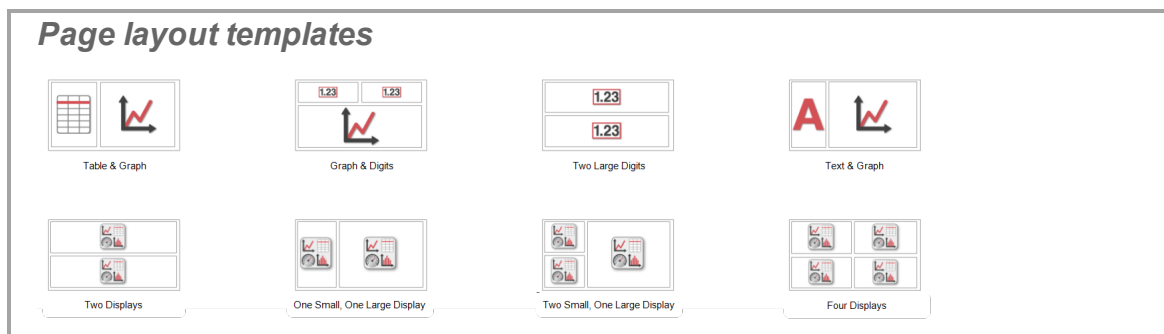
<sup>1</sup>You can add or remove display elements after you select a template.



2. The selected layout appears in the Display Area
3. Click **<Select Measurement>** for each display.
4. Select the measurement in the drop-down list.

**Note:** Sensors that are inactive (not ready to collect data) appear with .  
( See "Connecting an interface": 17 and "Connecting a sensor": 27)

## Use drag & drop and targets

1. Click one of the display icons in the Displays Palette.  
 See "Tour the Displays Palette ": 10 for a description of the available display types.
2. Drag the icon to the Display Area.
3. Use the Page layout targets to quickly format the page.

### Using Page layout targets


1. Drag a display icon to one of the five targets.
2. Move the cursor around open area of the page.
3. The target turns darker as you approach it and is selected when the border on the target appears.
4. Drop the icon on a selected target.

 Page targets divide the available area into the following five sections:  **Tip<sup>1</sup>**

**Note:** Page targets divide either an open area or an existing display.



<sup>1</sup>Use targets to quickly divide an existing display into two displays.

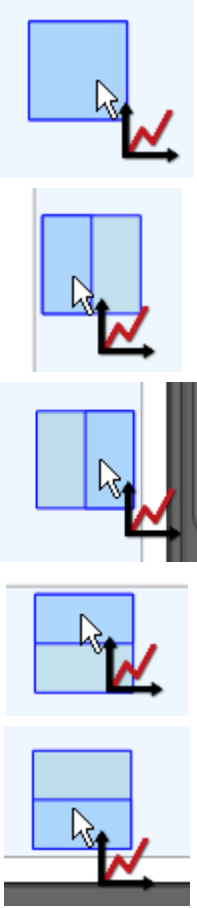**Page layout targets**



- entire available area (not used for splitting an existing display)
- left half (divided vertically—left half shows darker than right)
- right half (divided vertically—right half shows darker than left)
- top half (divided horizontally—top half shows darker than bottom)
- bottom half (divided horizontally—bottom half shows darker than top)

4. The selected layout appears in the Display Area **Tip<sup>1</sup>**
5. Add more displays noting the Page layout targets.
6. Click the **<Select Measurement>** button for each display.
7. Select the measurement in the drop-down list.

**Note:** Sensors that are inactive (not ready to collect data) appear with .  
( See "Connecting an interface": 17 and "Connecting a sensor": 27)

## Moving and resizing


Capstone uses handles to move and resize individual displays on a page. Capstone includes vertical and horizontal alignment borders for easy display alignment.

### Moving a display

1. Click on the display you wish to move to make it active.

<sup>1</sup>You can add or remove display elements after you select a template.


2. Move the cursor until the border darkens and the move (hand) symbol appears.

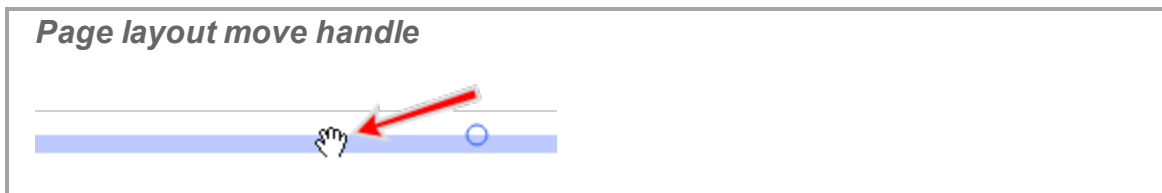


3. Click and drag the display to new location.


**Note:** The alignment guides appear (dotted lines) when the display border aligns with adjacent displays. 

## Resizing a display

1. Click on the display to move to make it active.
2. Move the cursor to one of the eight resize handles (circle ).

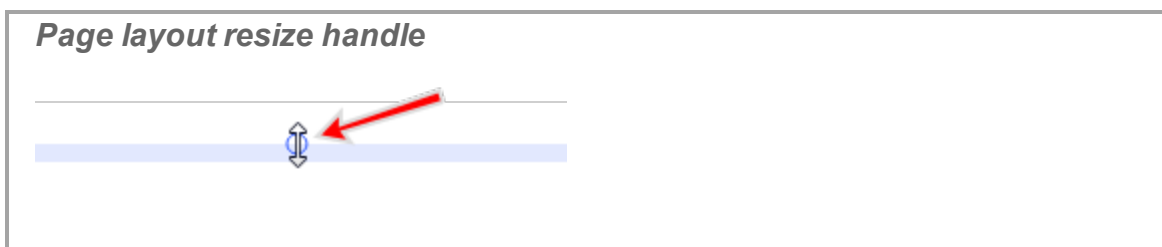


3. Click and drag the border to new position.

**Note:** The alignment guides appear (dotted lines) when the display border aligns with adjacent displays. 

## Realigning displays

1. Move the cursor between displays until the desired move border symbol appears.



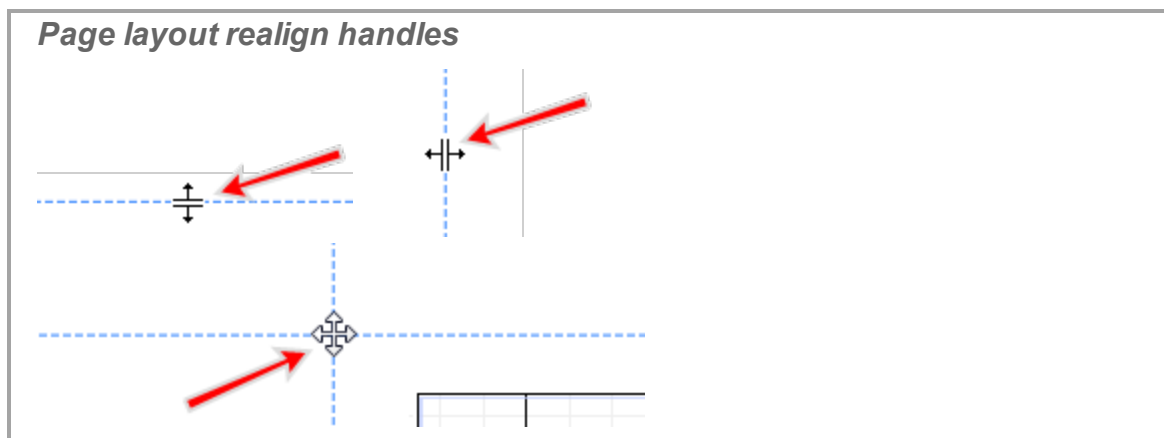

2. Click and drag the border to new position.

This action might affect multiple displays.

ⓘ It might be necessary to move displays to align to a common alignment border. (  
🔗 See "Moving a display": 58)

➡ Next step: **"Starting an experiment": 60**

## Starting an experiment

Capstone provides several controls in the Controls Palette for collecting data.

### Pre-work


1. Connect one or more interfaces. (🔗 See "Connecting an interface": 17)
2. Connect one or more sensors. (🔗 See "Connecting a sensor": 27)
3. Add one or displays. (🔗 See "Adding a display ": 56)





**Note:** Sensors that are inactive (not ready to collect data) appear with ⚠.  
(🔗 See "Connecting an interface": 17 and "Connecting a sensor": 27)

### Start data collection with default settings

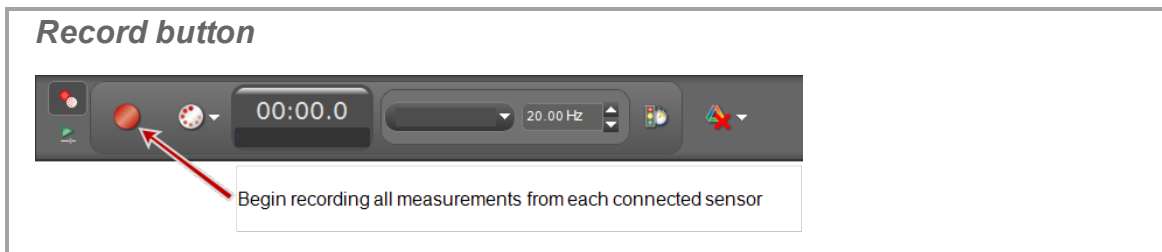**Note:** Measurements in displays that are inactive (not ready to collect data) appear with ⚠. Additionally, the data record button turns gray  if there is not at least one active interface and sensor.  
(🔗 See "Connecting an interface": 17 and "Connecting a sensor": 27)





1. Click **Record**  in the Controls Palette.




- ➔ Capstone begins recording all available data.
  - ➔ **Record**  toggles to **Stop** .
2. Click **Stop**  in the Controls Palette to stop data collection.
    - ➔ Collected data appear in all displays of the current page.
    - ➔ The data run is listed in the legend for each display.
    - ➔ Recorded data appear on the **Sensor Data Summary** tab () in **Data Summary**, and is selectable from any data display with a default name **Run #%1**, where %1 is an automatically generated run number.

 See "[How do I record data in PASCO Capstone?](#) ": 73 for details.

## Removing data runs

- Delete Last Run: click **Delete Last Run**  in the Controls Palette .
- Delete Selected Runs: click the down arrow next to **Delete Last Run**  and select a specific run.
- Delete All Runs: click the down arrow next to **Delete Last Run**  and select **Delete All Runs**.  **Tip<sup>1</sup>**

**Note:** The run data is removed from all displays, legends, and the data summary. Click the **Undo** button  at the top to recover the data.

 Next step (optional): [Downloading experiments](#)

## Downloading Capstone experiments

PASCO provides several ready-made experiments for Capstone.  **Tip<sup>2</sup>**

<sup>1</sup>You can delete all data for the run and repeat data collection.

<sup>2</sup>The lab might contain multiple pages including: ◇ introduction ◇ theory ◇ setup & procedures ◇ analysis ◇ conclusions

1. Go to the Capstone experiments web page: <http://www.pasco.com/family/capstone-experiments/index.cfm>
2. Select a subject area, for example **Thermodynamics**.
3. Select the **Free Experiments Files** tab.
4. Scroll through the list of ready-made experiments and click to download the experiment files.
5. Copy the files from your download folder to your Capstone experiments folder.
  - ➡ The download might contain the following:
    - Capstone application files; these have a ".cap" file extension
    - Editable lab guides
    - Images
6. Open the \*.cap file from Capstone.
  - a. Click **File > Open Experiment...**
  - b. Navigate to the experiment file.
  - c. Click **Open**.

## 3 How do I...

### Installation and Update

"How do I automate a network installation?": 66

"How do I force an update to an interface?": 66

### Experiment Files

"How do I download and open PASCO ready-made experiments?": 67

"How do I save my experiment?": 68

"How do I export all data?": 68

"How do I open DataStudio experiments and data?": 68

"How do I open Xplorer GLX, SPARK SLS, and other non-Capstone experiments and data?": 69

"How do I import data from Microsoft Excel\*? ": 69

"How do I export data to Microsoft Excel\*?": 70

"How do I import Xplorer GLX data?": 71

### Data and Measurements

"How do I record data in PASCO Capstone? ": 73

"How do I delete data runs?": 79

"How do I choose one or more runs to display?": 79

"How do I identify a measurement when I have multiple of the same type?": 80

"How do I change the numerical format of data?": 81

"How do I change the numerical format of statistics?": 83

"How do I exclude, delete, and restore selected data?": 84

"How do I edit a measurement?": 87

"How do I create a pure mathematical function?": 87

### Workbooks

"How do I show, hide, and lock the palettes?": 89

"How do I change what display toolbar buttons are visible?": 90

### Sensors

"How do I zero a measurement?": 92

"How do I switch the sign on a sensor?": 93

"How do I calibrate a sensor?": 94

"How do I calibrate a motion sensor?": 95

"How do I calibrate a drop counter?": 96

"How do I calibrate a pH sensor?": 97

## Calculations

"How do I create a calculation? ": 99

"How do I use my calculation?": 102

"How do I select a measurement in the calculator?": 103

"How do I see the result of a calculation?": 103

"How do I reorder calculations?": 105

## Plots

"How do I plot multiple measurements on separate plot areas with shared X-axis?": 107

"How do I zoom to a selection?": 108

"How do I pan and scale?": 110

"How do I select active data?": 111

"How do I plot multiple measurements on same plot with multiple Y-axes?": 113

"How do I plot multiple measurements on same plot with same Y-axes?": 114

"How do I apply a curve fit to a portion of the data?": 117

"How do I change curve fit parameters?": 118

"How do I adjust the weights for the weighted linear fit?": 119

"How do I lock and unlock the axes of a graph?": 121

"How do I change layer order of a display?": 123

## Tables

"How do I reorder columns in a table?": 124

"How do I use a table to build data in Keep Mode?": 124

## Text Boxes

"How do I enter subscripts, superscripts, Greek letters, and symbols?": 127

"How do I enter an equation into a text display?": 127

## Scope

"How do I compare scope traces from the same sensor when monitoring?": 130

"How do I select the scope trigger?": 131

"How do I pre-trigger?": 134

"How do I freeze and record the scope display?": 135

## Video

"How do I add a video source?": 137

"How do I analyze a video?": 138

"How do I manually synchronize video with data?": 140

"How do I adjust the frame rate of a recorded video?": 141

"How do I adjust the resolution of a video?": 141

"How do I make files with videos smaller?": 141

"How do I troubleshoot video problems?": 142

## How do I... Installation and Update

"How do I automate a network installation?": 66

"How do I force an update to an interface?": 66

### How do I automate a network installation?

You must have a valid site license to perform a network installation. PASCO Capstone uses an [InstallShield](#) installer, which packages the [Windows Installer](#) (msi) into an executable setup file: `PASCO_Capstone_Setup.exe`. Both the InstallShield installer and the packaged msi can be controlled using command line switches: See [InstallShield Note 1039936](#) for details on the InstallShield switches and [Microsoft Note 227091](#) for information on the msi switches.

Below is an example of a silent installation:

1. Copy the installation file `PASCO_Capstone_Setup.exe` that you have either downloaded or obtained from a CD to the root directory of your hard drive "C:".
2. Go to **Start > Run...**
3. Type `C:/PASCO_Capstone_Setup.exe /s /v"/qb"`. (/s = installer runs silently, /v = passes the parameter that follows immediately after in quotes to the msi installer, /qb = msi runs with no modal dialog boxes.)
4. Run the software.
5. Enter your license key when prompted.
6. After the license key is accepted, copy `ResourceFiles.rc` located at `C:\Documents and Settings\All Users\Application Data\PASCO Scientific\ResourceFiles.rc` (Windows XP) or `C:\Program Data\PASCO Scientific\ResourceFiles.rc` (Window 7) to the same location on the other computers by means of a [batch copy](#) command or other mechanism.

Contact "[Teacher and Technical Support](#)": [ii](#) for instructions to install on Windows 8 and MAC.

### How do I force an update to an interface?

PASCO Capstone prompts to update the interface if a new firmware version exists.

Follow these steps if an interface requires re-installing existing firmware to correct firmware corruption issues:

1. Unplug the interface from the USB port or turn it off.
2. Run PASCO Capstone.
3. Hold down **CTRL+U** while connecting or turning on the interface.

## How do I... Experiment Files

"How do I download and open PASCO ready-made experiments?": 67

"How do I save my experiment?": 68

"How do I export all data?": 68

"How do I open DataStudio experiments and data?": 68

"How do I open Xplorer GLX, SPARK SLS, and other non-Capstone experiments and data?": 69

"How do I import data from Microsoft Excel\*? ": 69

"How do I export data to Microsoft Excel\*?": 70

"How do I import Xplorer GLX data?": 71

### How do I download and open PASCO ready-made experiments?

PASCO provides several ready-made experiments for Capstone.  **Tip**<sup>1</sup>

1. Go to the Capstone experiments web page: <http://www.pasco.com/family/capstone-experiments/index.cfm>
2. Select a subject area, for example **Thermodynamics**.
3. Select the **Free Experiments Files** tab.
4. Scroll through the list of ready-made experiments and click to download the experiment files.
5. Copy the files from your download folder to your Capstone experiments folder.
  - ➡ The download might contain the following:
    - Capstone application files; these have a ".cap" file extension
    - Editable lab guides
    - Images
6. Open the \*.cap file from Capstone.
  - a. Click **File > Open Experiment...**
  - b. Navigate to the experiment file.
  - c. Click **Open**.

<sup>1</sup>The lab might contain multiple pages including: ◇ introduction ◇ theory ◇ setup & procedures ◇ analysis ◇ conclusions



## How do I save my experiment?

Click **File** > **Save Experiment**.

All of the data, displays, toolbar states and formatting are saved.


**Note:** Deleted data and undo history are not saved.

## How do I export all data?


1. Click **File** > **Export Data...**
2. Select a file type: **Text (Tab delimited) (\*.txt)**; **CSV (Comma delimited) (\*.csv)**.
3. Navigate to the location to save the file.
4. Click **Save**.

## How do I open DataStudio experiments and data?

PASCO Capstone can open and convert most DataStudio 1.9 files to Capstone format.

1. Click **File** > **Open Experiment...**
2. Click on the file type pull-down menu, and select: **DataStudio 1.9 files (\*.ds)**
3. Click **Open**.
  - ➔ Capstone places *each* display from the DataStudio on separate Capstone work-book pages.
  - ➔ Each display is sized to the full page.
  - ➔ Each page (tab) is named with the display name.
4. Recommended: save the experiment as a Capstone file. ( See "[How do I save my experiment?](#)": 68)
5. Re-layout the Capstone pages as desired.

To combine multiple displays on a single page:

  - a. Resize and move the display that will contain multiple displays to create a space for additional display(s). ( See "[Moving and resizing](#)": 58)
  - b. Click on page tab containing display to move.
  - c. Resize the display smaller than full page.
  - d. Select the display border, right click and select Cut.

- e. Click on the page, right-click and select Paste.
- f. Resize and move as desired.

## How do I open Xplorer GLX, SPARK SLS, and other non-Capstone experiments and data?

### Xplorer GLX

PASCO Capstone can open and convert most DataStudio 1.9 and Xplorer GLX files to Capstone format.

(For files resident on the Xplorer GLX: see ["How do I import Xplorer GLX data?": 71](#))

1. Click **File > Open Experiment...**
2. Click on the file type pull-down menu, and select: **GLX Files (\*.glx)**
3. Click **Open**.
4. Recommended: save the experiment as a Capstone file. ([See "How do I save my experiment?": 68](#))

### Xplorer and 500 Dataloggers

TBD

### SPARK SLS Experiments


TBD

## How do I import data from Microsoft Excel\*?

### Tip<sup>1</sup>

1. Open Microsoft Excel with the file containing your data.
2. Highlight the data within one column.
3. Press **CTRL+C** to copy the data.
4. Open PASCO Capstone.
5. Create a Table. ([See "Adding a display ": 56](#))
6. Click on **<Select Measurement>** in a column.

<sup>1</sup>Capstone supports export as tab-delimited text, which can be imported into almost any spreadsheet or data analysis program. Note that Capstone is a complete data collection, visualization, and analysis tool; therefore, it is generally easier to analyze the data in Capstone than to export to another software program for analysis.

7. Select **Create New > User-Entered Data**.
8. Optional: Rename the measurement and give it appropriate units.
9. Click in the cell of the **User-Entered Data** column where you would like to have the data pasted.
10. Press **CTRL+V** to paste.
11. Repeat for additional columns of data.
  - ➔ Each column of data is a separate data set.
  - ➔ Imported data are listed in the **User-Entered Data Summary** tab () of **Data Summary**.


## How do I export data to Microsoft Excel\*?

### Tip<sup>1</sup>

#### Export All Data

1. Select **File > Export Data...**
2. Select **CSV (comma delimited)** as the file type.
3. Click **Save**.

#### Export Selected Data

1. Create a Table. ( See "Adding a display ": 56)
2. Select the data that you would like to copy:
  - All data: click anywhere in the table and press **CTRL+A**.
  - All of the data in a column: click anywhere on the heading of the column where there is no text.
  - Visible part of a table: click and drag to select data.
3. Press **CTRL+C** to copy data to clipboard.
4. Open Microsoft Excel\*.
5. Click in a cell where you would like the data to appear.
6. Press **CTRL+V** to paste data.

<sup>1</sup>Capstone supports export as tab-delimited text, which can be imported into almost any spreadsheet or data analysis program. Note that Capstone is a complete data collection, visualization, and analysis tool; therefore, it is generally easier to analyze the data in Capstone than to export to another software program for analysis.


**Note:** You can define the format in Excel to include significant digits. (Format as **Number** with a defined number of decimal places.)

## How do I import Xplorer GLX data?


**Note:** For Xplorer GLX files on the host computer: see "[How do I open DataStudio experiments and data?](#)": 68

To transfer experiments (including data files) from the Xplorer GLX to Capstone:

1. Connect the Xplorer GLX to Capstone.
2. Click **Hardware Setup** in the Tools Palette: .

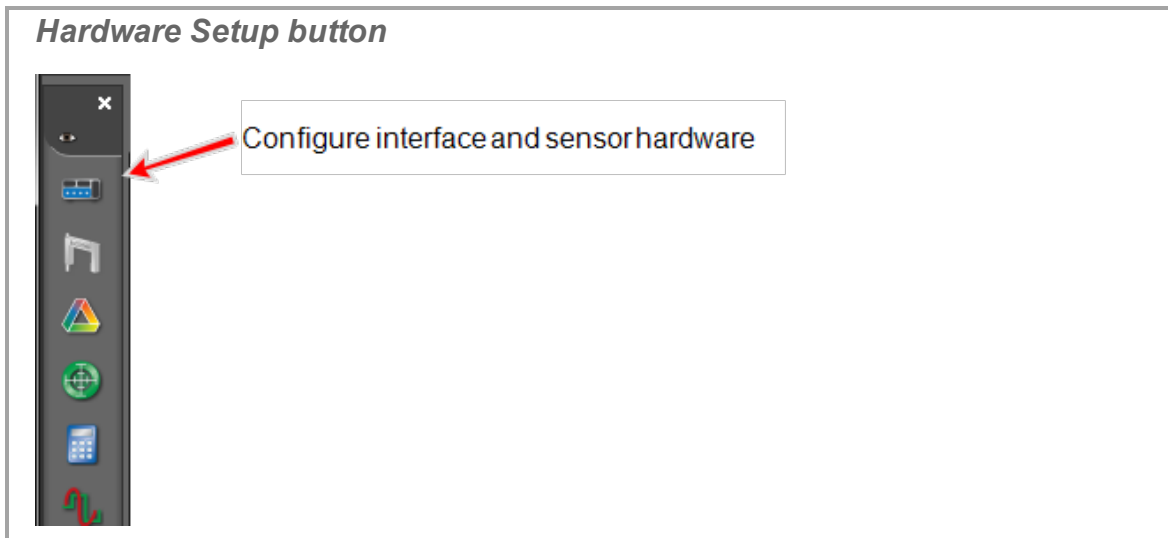

3. Click the yellow file folder icon next to the Xplorer GLX: .

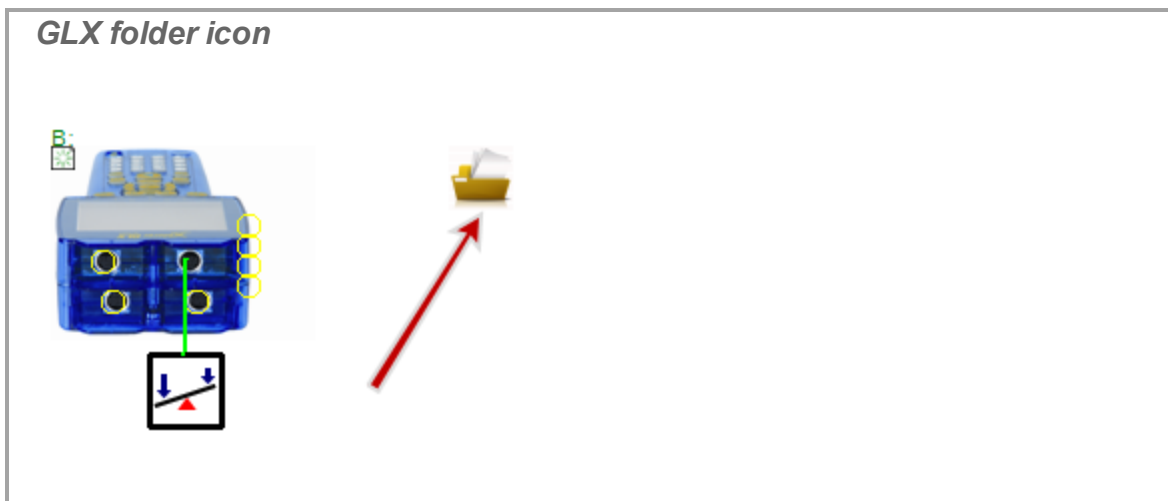

4. Click **Download All**.
5. Navigate a folder and click **Select Folder**.
6. Click **File > Open Experiment...**
7. Select **GLX Files (\*.glx)** from the file type drop down.
8. Navigate to the saved file .
9. Click **Open**.

Capstone opens the experiment and data for analysis and further data acquisition.

## How do I... Data and Measurements

"How do I record data in PASCO Capstone? ": 73

"How do I delete data runs?": 79

"How do I choose one or more runs to display?": 79

"How do I identify a measurement when I have multiple of the same type?": 80

"How do I change the numerical format of data?": 81

"How do I change the numerical format of statistics?": 83

"How do I exclude, delete, and restore selected data?": 84


"How do I edit a measurement?": 87




"How do I create a pure mathematical function?": 87

### How do I record data in PASCO Capstone?

There are three data recording modes in PASCO Capstone. Data recording is controlled by the settings on the record tab of the Controls Palette.

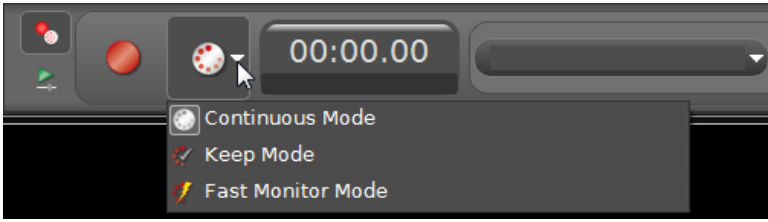*Record modes*





	<b>Continuous Mode</b>	Set to record data at the current sample rate settings
	<b>Keep Mode</b>	Set to keep data points when commanded
	<b>Fast Monitor Mode</b>	Set to acquire data without recording; required for high speed sampling in Scope and FFT displays

### Detailed Explanations


## Continuous Mode *Set to record data at the current sample rate settings*


### Behavior

- All measurement **Data**<sup>1</sup> continuously records as soon as **Record**  is clicked in the Controls Palette until the **Stop**  is clicked (or the hardware is no longer able to process the data—at which time data collection will stop with an error message).
- Data records at the specified sample rate(s).


 See "[What is Continuous Mode?](#)": 168


### Procedure

1. Select  **Continuous Mode** in the Controls Palette.
2. Adjust the sample rates of each of the measurements:
  - ➡ Ensure sensors are sufficiently responsive for desired data while low enough so that the computer system can keep up with data sampling and does not generate excessive file sizes.

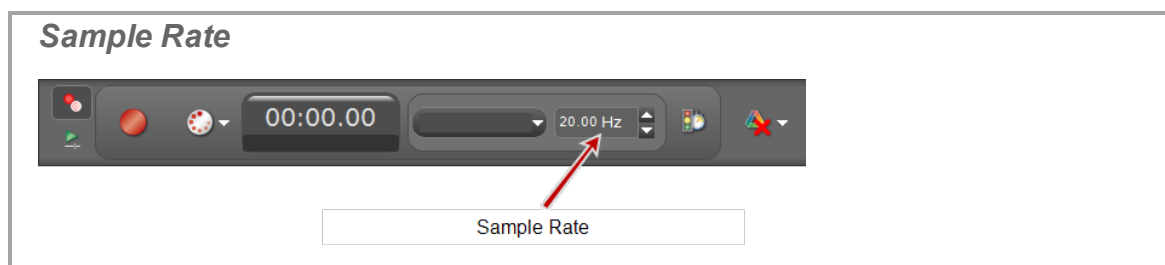

4. Click **Record**  in the Controls Palette.

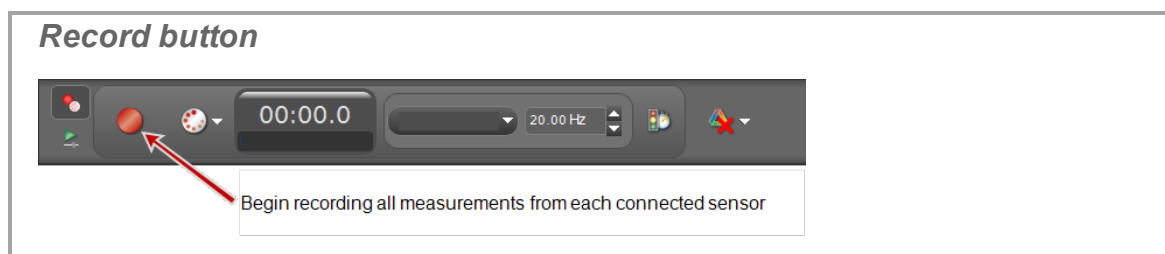

5. Click **Stop**  in the Controls Palette to stop data collection.
  - ➡ Collected data appear in all displays of the current page.

<sup>1</sup>A data set including measurement and user-defined data from sensors and time data.

- ➔ The data run is listed in the legend for each display.
- ➔ Recorded data appear on the **Sensor Data Summary** tab (📎) in **Data Summary**, and is selectable from any data display with a default name **Run #%1**, where %1 is an automatically generated run number.


 **Keep Mode** *Set to keep data points when commanded*


**Behavior**

- Manual recording mode.
- The measurement sample rate(s) determine the rate at which data samples are *previewed* (not recorded).
- **Data**<sup>1</sup> records when and only when **Keep Sample** ✓ is clicked in the Controls Palette while monitoring data.
- One instance of all data is recorded each time **Keep Sample** ✓ is clicked.


**Procedure**

 Tip: You can use a table with **Keep Mode** to build data for specific points. (📎 See "[How do I use a table to build data in Keep Mode?](#)": 124)\_

1. Select  **Keep Mode** in the Controls Palette.
2. Adjust the sample rates of each of the measurements so that the sensors are sufficiently responsive for desired data.

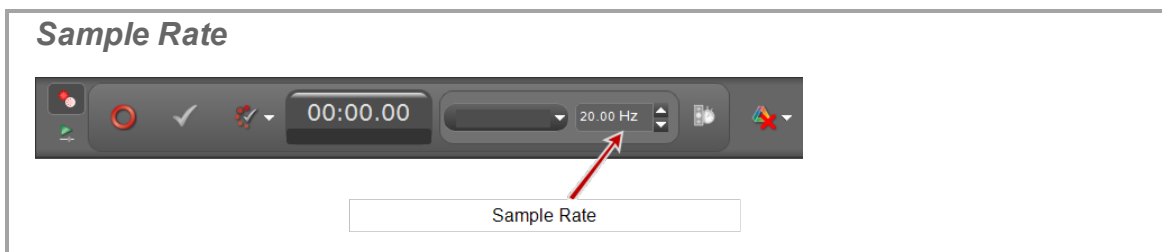



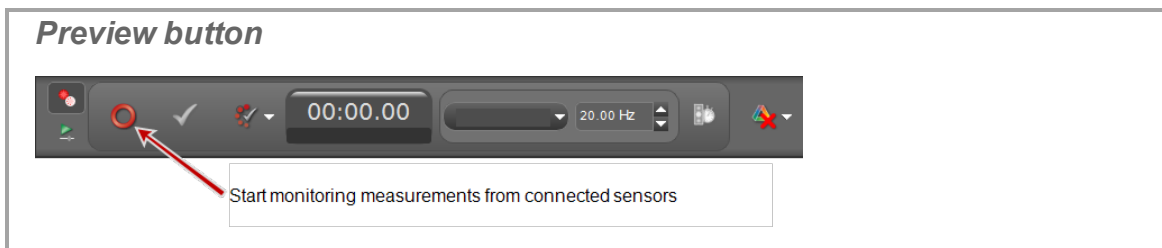


3. Click **Preview**  in the Controls Palette.






<sup>1</sup>A data set including measurement and user-defined data from sensors and time data.



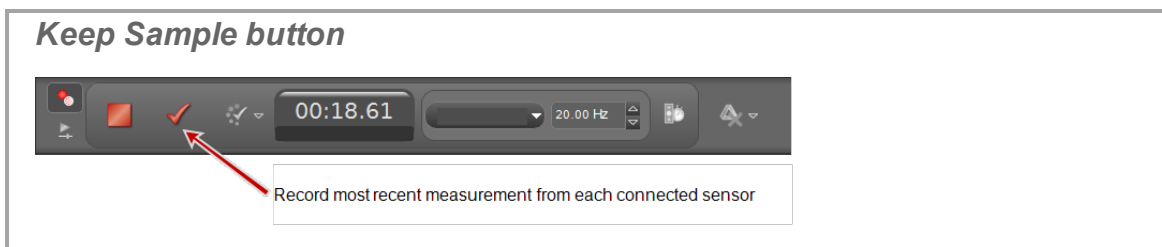

- ➔ Capstone begins monitoring data and displays previews.
  - ➔ **Preview**  toggles to **Stop** .
  - ➔ **Keep Sample**  button is active.
4. Click **Keep Sample**  to record the data currently monitored.
- ➔ One instance of all data is recorded.



5. Repeatedly click **Keep Sample**  to record data.
- ➔ One instance of all data is recorded each time **Keep Sample**  is clicked.
6. Optional: Replace values in a table. (See ["Replacing a current value in a table": 125.](#))




**Note:** You can replace data in tables while in preview mode for the run (before clicking **Stop** .

See ["How do I edit a measurement?": 87](#) to edit data after you click **Stop** .

7. Click **Stop**  in the Controls Palette to stop observing the data.
- ➔ Collected data appear in all displays of the current page.
  - ➔ The data run is listed in the legend for each display.
  - ➔ Recorded data appear on the **Sensor Data Summary** tab () in **Data Summary**, and is selectable from any data display with a default name **Run #%1**, where **%1** is an automatically generated run number.


**Fast Monitor Mode** Set to acquire data without recording; required for high speed sampling in Scope and FFT displays

**Behavior**

- Intended for use with workbooks pages containing the FFT or Scope displays.
- Designed for viewing periodic data or triggered transient events.
- **Monitor**  displays data *without* recording.
- **Snapshot**  records the monitored data.
  - Only **Data**<sup>1</sup> that is currently visible in the display is recorded each time Snapshot  is clicked.
  - Unsaved buffer data is also recorded when closing Capstone.
- Can repeatedly monitor and record data points.



 See "What is Fast Monitor Mode?": 169

**Procedure**

1. Select  **Fast Monitor Mode** in the Controls Palette.
2. Adjust the sample rates of each of the measurements so that the sensors are sufficiently responsive for desired data.




3. Select one or more scope or FFT displays. ( See "Adding a display ": 56)


**Note:** To record a *single event*, use the scope display and click the single-shot button . ( See "What does the Single Shot button do?": 178)


4. Click **Monitor**  in the Controls Palette.


<sup>1</sup>A data set including measurement and user-defined data from sensors and time data.

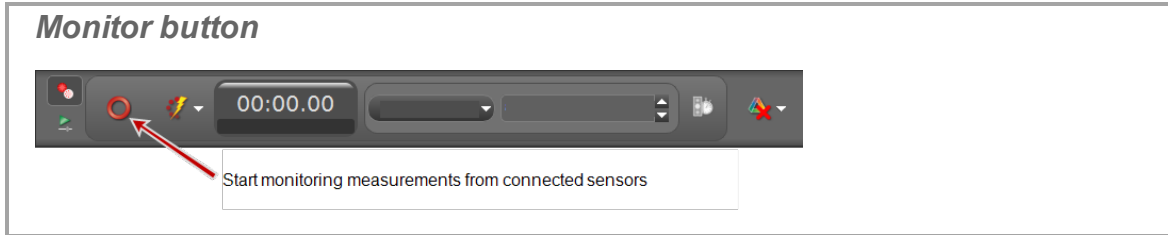

5. Click **Stop**  in the Controls Palette to stop observing the data.

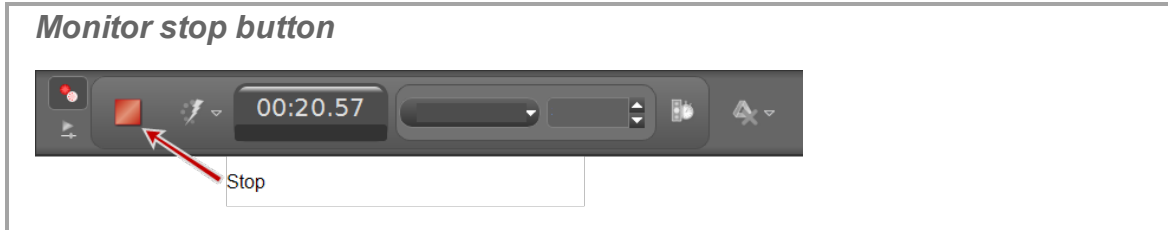

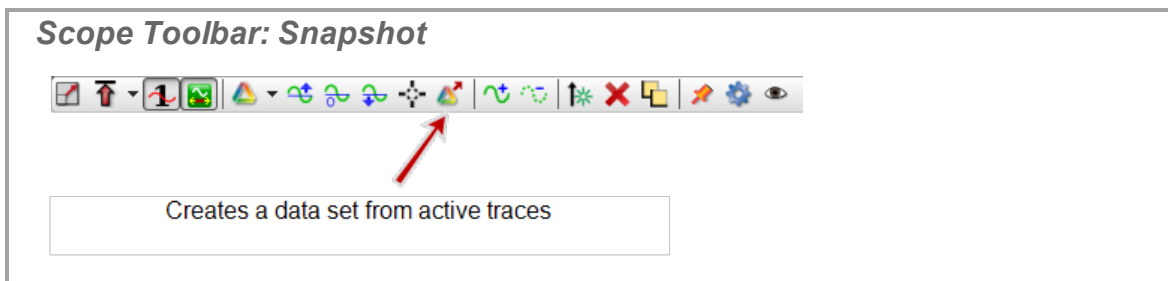

6. Scale the data to fit the data to record within the displays. ([See "How do I pan and scale?": 110](#))
7. Click **Snapshot**  in the scope toolbar.









- Capstone records the data.


**Notes:**

Only the data that is visible in the plot area is recorded as a run when **Snapshot**  is clicked.

Recorded data appear on the sensor tab in the Data Summary tab () and is selectable from any data display with a default name **Monitor Run--ScopeSnapshot%1** where **%1** is an automatically generated run number.

## How do I delete data runs?

- Delete Last Run: click **Delete Last Run**  in the Controls Palette .
- Delete Selected Runs: click the down arrow next to **Delete Last Run**   and select a specific run.
- Delete All Runs: click the down arrow next to **Delete Last Run**   and select **Delete All Runs**.  **Tip<sup>1</sup>**


**Note:** The run data is removed from all displays, legends, and the data summary. Click the **Undo** button  at the top to recover the data.

## How do I choose one or more runs to display?

By default, PASCO Capstone displays the last run recorded or the currently selected measurement in Replay mode.

To select a different run to display:

1. Click the pull-down arrow (▼) next to **Run Selection**   in the toolbar.

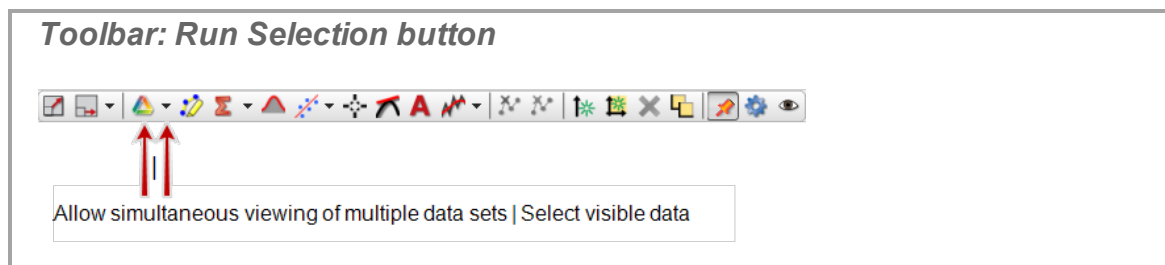


2. Select the run to display.

To display multiple runs:

<sup>1</sup>You can delete all data for the run and repeat data collection.

1. Click to depress **Run Selection**  in the toolbar.

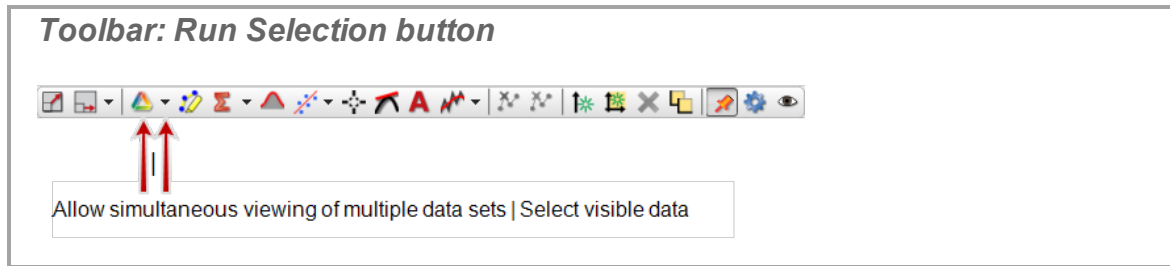

Selected runs display simultaneous with the button is depressed.

2. Click the pull-down arrow (▼) next to **Run Selection**   in the toolbar.
3. Select the one or more runs to display.

## How do I identify a measurement when I have multiple of the same type?

### One instance of a measurement

Format for a single instance of a specific measurement:

*Measurement Type (units)*

**Example: Force (N)**

### Multiple instances of measurement—one interface connected

Format for multiple instances of a specific measure with one interface connected:

*Measurement Type, Port (units)*

**Example: Force, Ch P1 (N)**

**Where:**

Measurement Type = Force

Port on interface = P1 (Port identifier on the interface "P1" → PASPORT sensor number 1)

### Multiple instances of measurement—multiple interfaces connected

Format for multiple instances of a specific measure with multiple interfaces connected:

*Measurement Type, Channel:Port (units)*

**Example: Force, Ch A:P1 (N)**


**Where:**

Measurement Type = Force

Channel = A (Interface identifier "A")


Port = P1 (Port identifier on the interface "P1" → PASPORT sensor number 1)

**Setup with multiple interfaces**

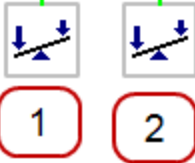

**Data Summary**


- ① Force, Ch A:P1 (N)
- ② Force, Ch A:P2 (N)
- ③ Force, Ch B:1 (N)

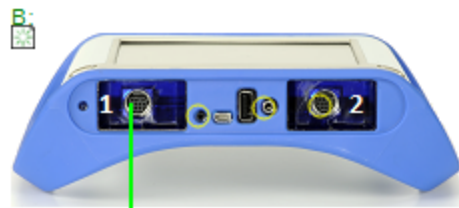



**How do I change the numerical format of data?**

**Note:** Changing data precision correspondingly changes the precision of statistics that use the data.

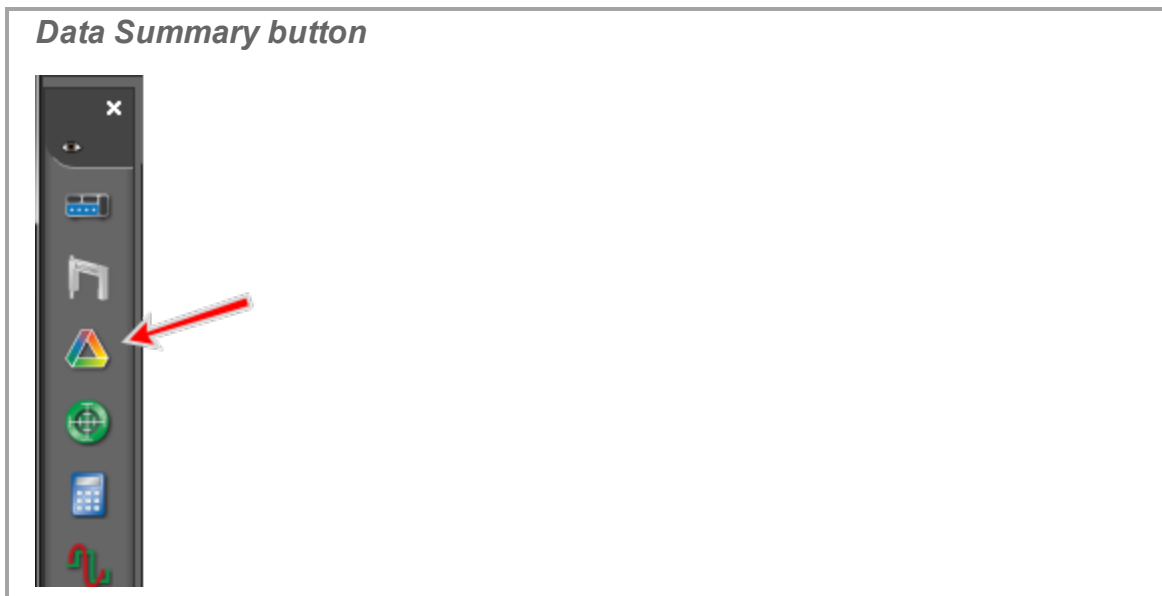1. Click **Data Summary**  in the Tools Palette.



2. Click the **Sensor Data** tab .
3. Click the measurement to change.
4. Click **Properties**  that appears to the right of the measurement.  
The **Properties** window opens.
5. Click **Numerical Format**.
6. Set the desired numerical format:

Numerical Format	Description
<b>Number of Decimal Places</b>	Set the number of digits that will appear to the right of the decimal.
<b>Apply Scientific Notation Thresholds</b>	<input checked="" type="checkbox"/> Select to enable Scientific Notation Thresholds
<b>Scientific Notation Thresholds</b>	Determines when the unit range transitions into scientific notation.
<b>Lower Exponent Threshold ≤ Upper Exponent Threshold ≥</b>	Apply limits for which the unit range transitions into scientific notation.
<b>Apply Zero Level Thresholds</b>	<input checked="" type="checkbox"/> Select to enable Apply Zero Level Thresholds
<b>Apply Zero Level Thresholds</b>	Apply the limit under which the number is considered to be zero.


7. Click **OK**.



## How do I change the numerical format of statistics?

**Note:** Changing data statistics precision correspondingly changes the precision of data.

To change the numerical format of the statistics, change the numerical format of the data on which the statistics are based.

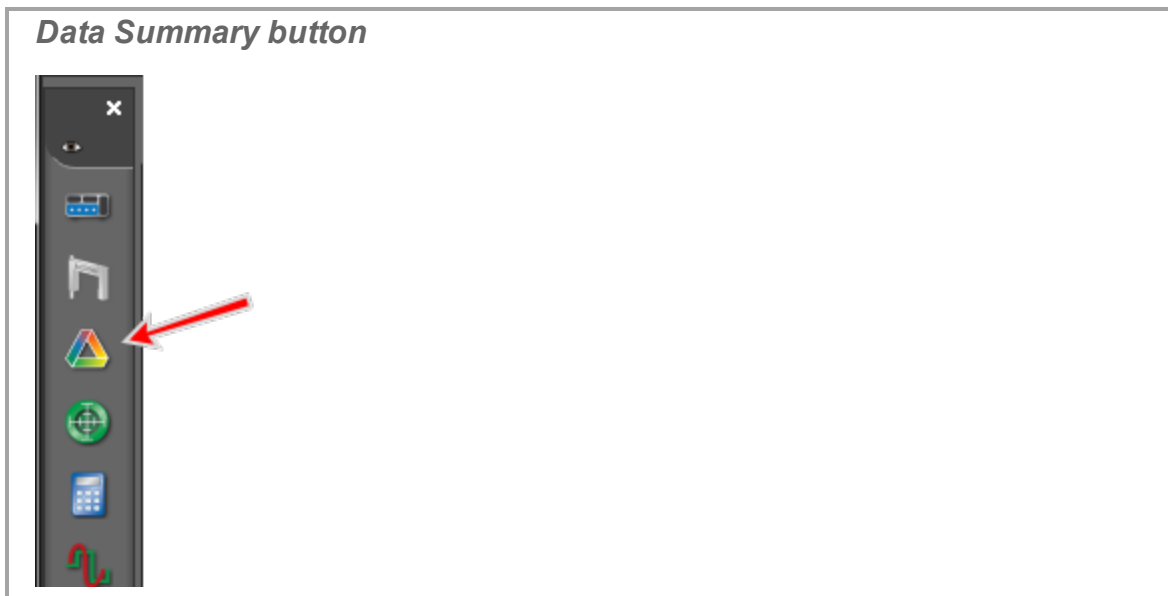1. Click **Data Summary**  in the Tools Palette.



2. Click the **Sensor Data** tab .
3. Click the measurement to change.
4. Click **Properties**  that appears to the right of the measurement.  
The **Properties** window opens.
5. Click **Numerical Format**.



6. Set the desired numerical format:


Numerical Format	Description
Number of Decimal Places	Set the number of digits that will appear to the right of the decimal.
Apply Scientific Notation Thresholds	<input checked="" type="checkbox"/> Select to enable Scientific Notation Thresholds
Scientific Notation Thresholds	Determines when the unit range transitions into scientific notation.
Lower Exponent Threshold $\leq$ Upper Exponent Threshold $\geq$	Apply limits for which the unit range transitions into scientific notation.
Apply Zero Level Thresholds	<input checked="" type="checkbox"/> Select to enable Apply Zero Level Thresholds
Apply Zero Level Thresholds	Apply the limit under which the number is considered to be zero.

7. Click **OK**.


## How do I exclude, delete, and restore selected data?

**Note:** You can create User-Entered Data to copy and edit/delete data in the list. (See "How do I edit a measurement?": 87)

### Table

- Highlight cells (click one cell or click and drag across multiple cells  **Tip<sup>1</sup>**).
- Exclude Selected Data:** Right-click and select **Exclude Data**.
- Delete Selected Data:** Right-click and select **Delete Cell Data**.

**Note:** Excluded data are masked in the data set and are not used by Capstone for data operations including display and calculations. Excluded data can be restored at any time.

**Note:** Deleted data are removed from the data set and the data cannot be restored at a later time. You can immediately undo a delete action by clicking **Undo** .

<sup>1</sup>Click the table heading to highlight all data in the table column.

*Deleted and Excluded Data in a Table*


0.900		0.30
0.950		<del>0.26</del>
1.000		<del>0.25</del>
1.050		<del>0.19</del>
1.100		<del>0.18</del>
1.150	2	<del>0.18</del>
1.200		<del>0.18</del>
1.250		<del>0.21</del>
1.300		<del>0.28</del>
1.350		0.35
1	1.400	0.42
	2.200	0.42

1 Deleted data are removed. Note the gap in time data.

2 Excluded data display with strikethrough.

## Graph


1. Click **Highlight range of points in active data**  in the toolbar.

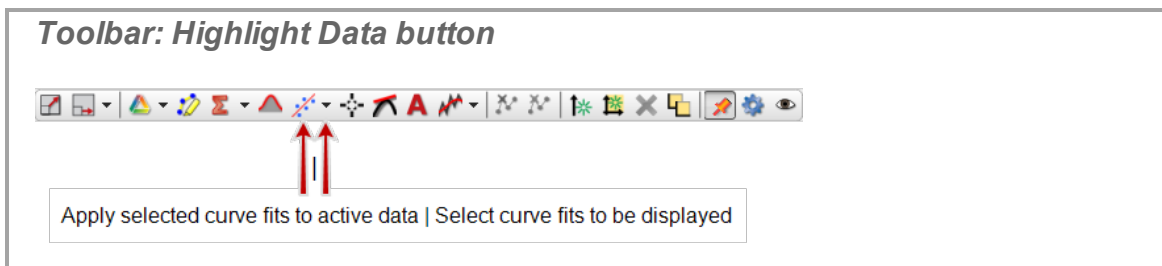


2. Move and resize the selection rectangle as needed.
3. **Exclude Selected Data:** Right-click and select **Exclusions > Exclude Highlighted Data Points**.
4. **Delete Selected Data:** Right-click and select **Deletions > Delete Highlighted Data Points**.
5. **Delete Excluded Data:** Right-click and select **Deletions > Delete ALL Excluded Data Points**.

**Note:** Excluded data are masked in the data set and are not used by Capstone for data operations including display and calculations. Excluded data can be restored at any time.

**Note:** Deleted data are removed from the data set and the data cannot be

restored at a later time. You can immediately undo a delete action by clicking **Undo** .

*Deleted and Excluded Data in a Plot*

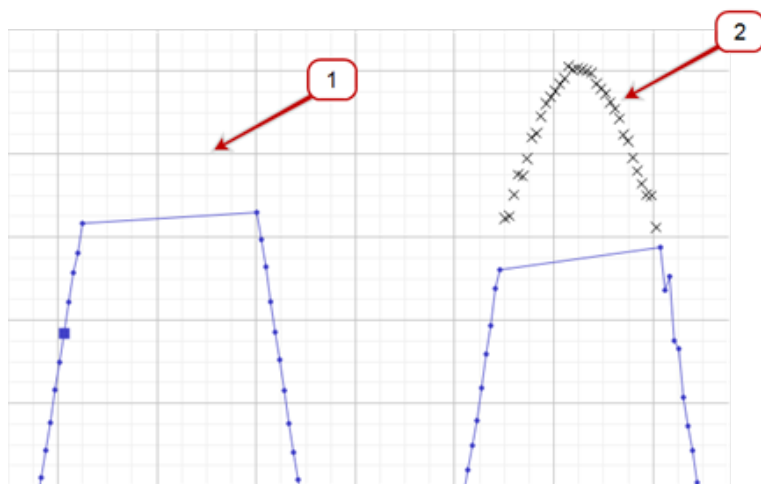


1 Deleted data are removed.

2 Excluded data display with "X" marks.


**To restore excluded data**

**Table**

1. Highlight excluded cells.  **Tip<sup>1</sup>**
2. Right-click and select **Restored Excluded Data**

**Graph**

1. Click **Highlight range of points in active data**  in the toolbar.

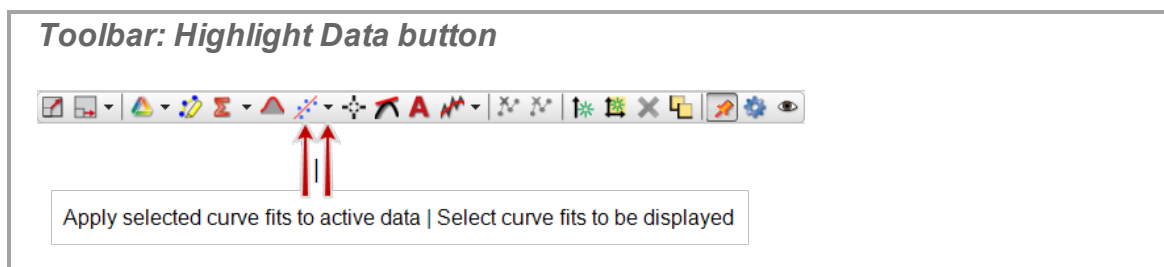

2. Move and resize the selection rectangle as needed to highlight excluded data points.
3. **Restore Selected Data:** Right-click and select **Exclusions > Restore Excluded Data Points Within Highlighter's Bounds**.
4. **Restore All Data:** Right-click and select **Exclusions > Restore ALL Excluded Data Points**.

<sup>1</sup>Click the table heading to highlight all data in the table column.

## How do I edit a measurement?

Measurements cannot be edited directly in PASCO Capstone. You can create User-Entered Data to copy and edit/delete data in the list.

**Note:** You can use Exclude Data to exclude selected data from analysis and Displays. (🔗 See "How do I exclude, delete, and restore selected data?": 84)

1. Create a Table. (🔗 See "Adding a display ": 56)
2. Click **<Select Measurement>** in the first column.
3. Select the measurement you would like edit.
4. Click **<Select Measurement>** in the second column.
5. Select **Create New > User-Entered Data**.
6. Click on the top of the first column.
7. Right click and select **Copy Values** (or use **CTRL+C**).
8. Click in the first cell of the **User Data %1** (where %1 is serial number) column.
9. Right-click and select **Paste Values** (or use **CTRL+V**).
10. Edit data in cells:
  - a. Change values in cells as needed.
  - b. Highlight cells and right-click to select: **Clear Values**, **Insert Cells**, **Delete Cell Data**, or **Exclude Data** as needed.
11. Click on **User Data %1** in the column heading, select rename, and enter a meaningful name and units.
 

➔ Imported data are listed in the **User-Entered Data Summary** tab (🔗) of **Data Summary**.

## How do I create a pure mathematical function?

Graph a purely mathematical function by defining a **model<sup>1</sup>** and then creating a calculation on the model within the calculator.

### Procedure

1. Create a model in the calculator.
  - a. Open the Calculator. (🔗 See "How do I create a calculation? ": 99)
  - b. Right click on a new calculation line and select **Insert Model Range**.


<sup>1</sup>A sequence of values within a range and a defined number of steps or step size.

The **Edit Model Range Properties** window opens.

- c. Set the **Start Value**: of the range.
- d. Set the **End Value**: of the range.
- e. Select option:

Option	Action	Example Format:
<b>Step Count:</b>	Set the number of values between the Start Value and End Value.	{0 .. 10;100} Start Value: 0; End Value: 10; Step Count: 100
<b>Step Size:</b>	Set the increment size between values in the range. A smaller step size will produce more values in the range and a larger step size will produce fewer values in the range.	{0 .. 10;+1} Start Value: 0; End Value: 10; Step Size: 1

2. Click **OK**.
3. Insert an equation using the Model on a new calculation line. ([See "How do I create a calculation?": 99](#))

Calculations appear on the **Calculation Data Summary** tab () in **Data Summary**, and are selectable from any data display. ([See "How do I use my calculation?": 102](#))

### Example in Calculator

$y = \{0 \dots 10;100\}$

$z = \sin(y)^2$

$w = \text{integral}([y(\text{units})], y)$

**Where:**

y is a model

z and w are functions based on model y.


## How do I... Workbooks

"How do I show, hide, and lock the palettes?": 89

"How do I change what display toolbar buttons are visible?": 90

### How do I show, hide, and lock the palettes?

1. Right-click anywhere in the gray area surrounding a workbook page.

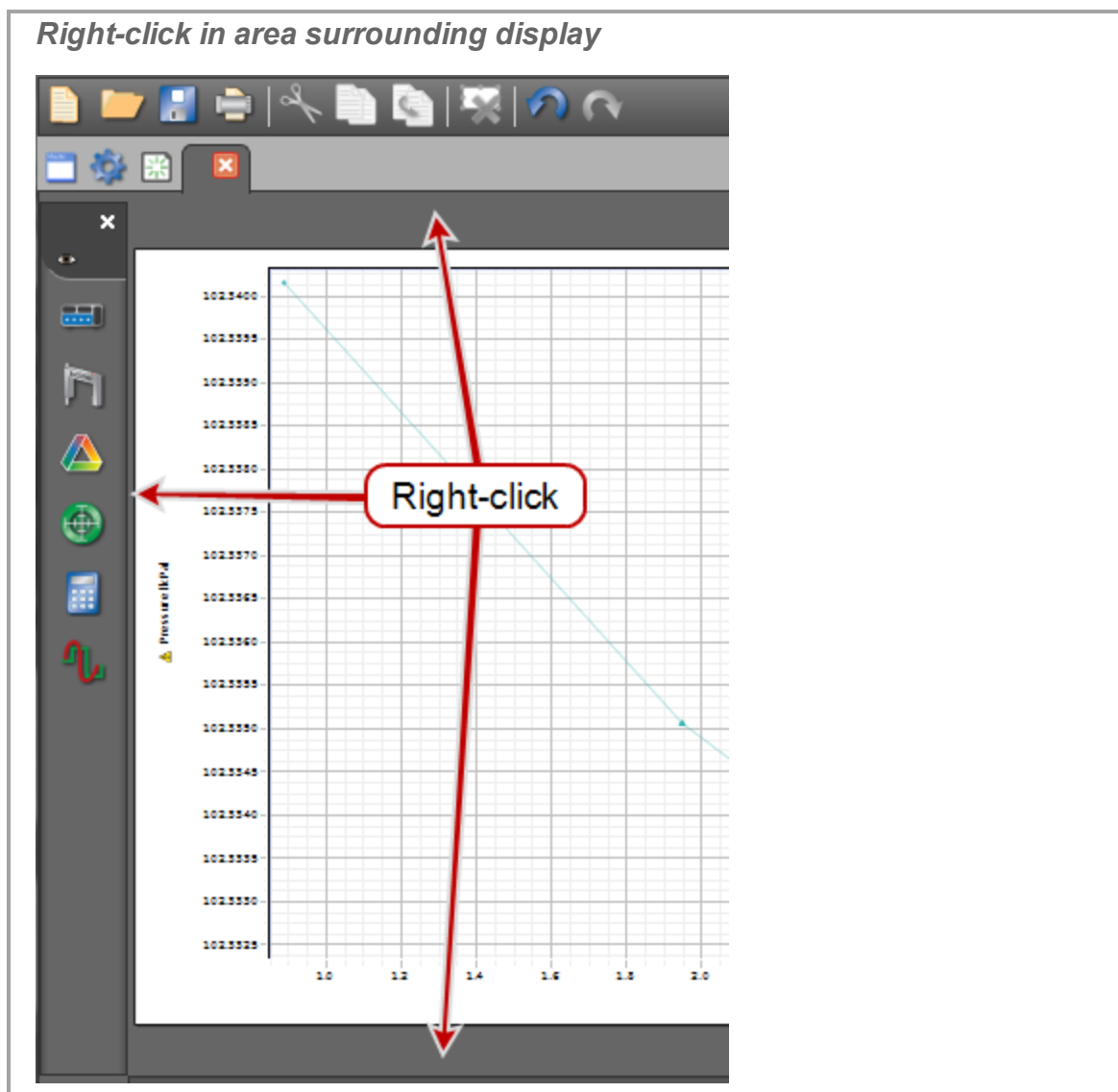

2. Show/hide: click the palette that you would like to show (☑) or hide (☐).


3. Lock/unlock: click **Lock Page** to lock (☑) or unlock page (☐).


➔ Locked pages:

- Displays cannot be added or removed.
- Text Entry Box elements are active and accept user entries.
- Displays can be scaled and panned.


## How do I change what display toolbar buttons are visible?

1. Click on **Visibility Menu**  in the display Toolbar.

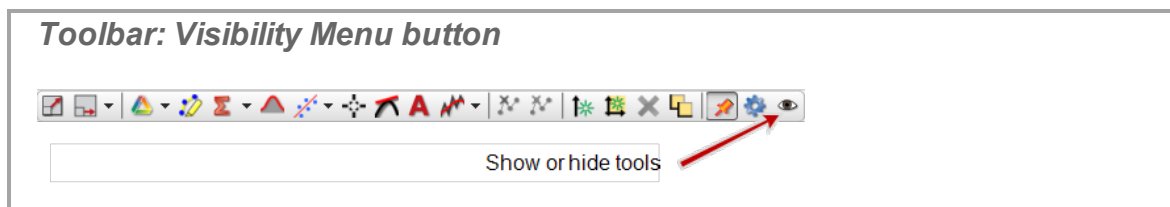

➔ If the **Visibility Menu** button  is not available: right-click anywhere in the gray

area surrounding a workbook page and deselect **Lock Page**.

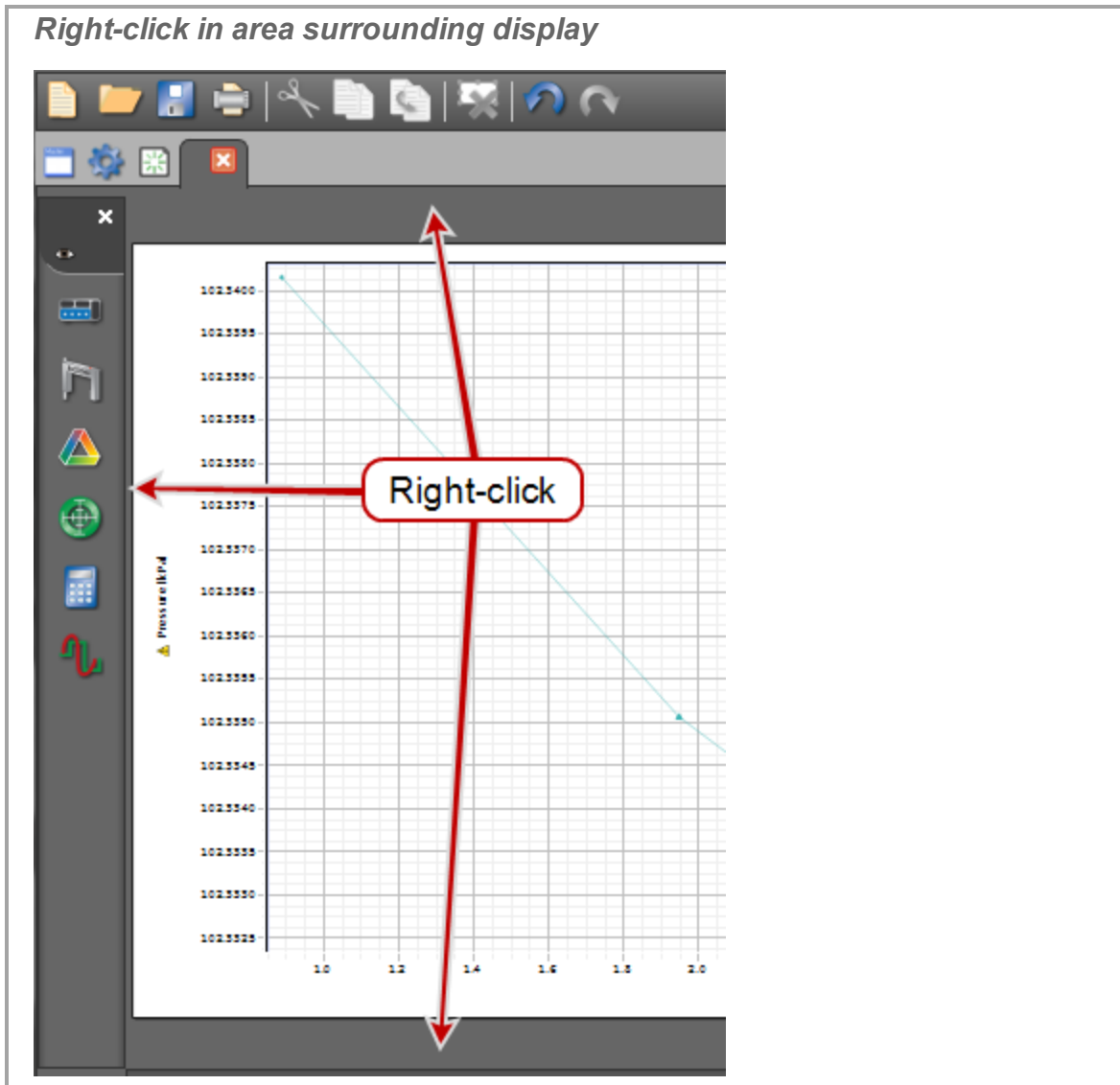

2. Click the buttons that you would like to show (☑) or hide (☐) on the toolbar.



## How do I... Sensors

"How do I zero a measurement?": 92

"How do I switch the sign on a sensor?": 93

"How do I calibrate a sensor?": 94

"How do I calibrate a motion sensor?": 95


"How do I calibrate a drop counter?": 96


"How do I calibrate a pH sensor?": 97


### How do I zero a measurement?

**Note:** If a sensor is selected in the Controls Palette that supports zeroing, the Controls Palette includes a **Zero Sensor Now** button .

#### Procedure

1. Connect the sensor to PASCO Capstone. (See "Connecting a sensor": 27)
2. Click **Hardware Setup** in the Tools Palette: .

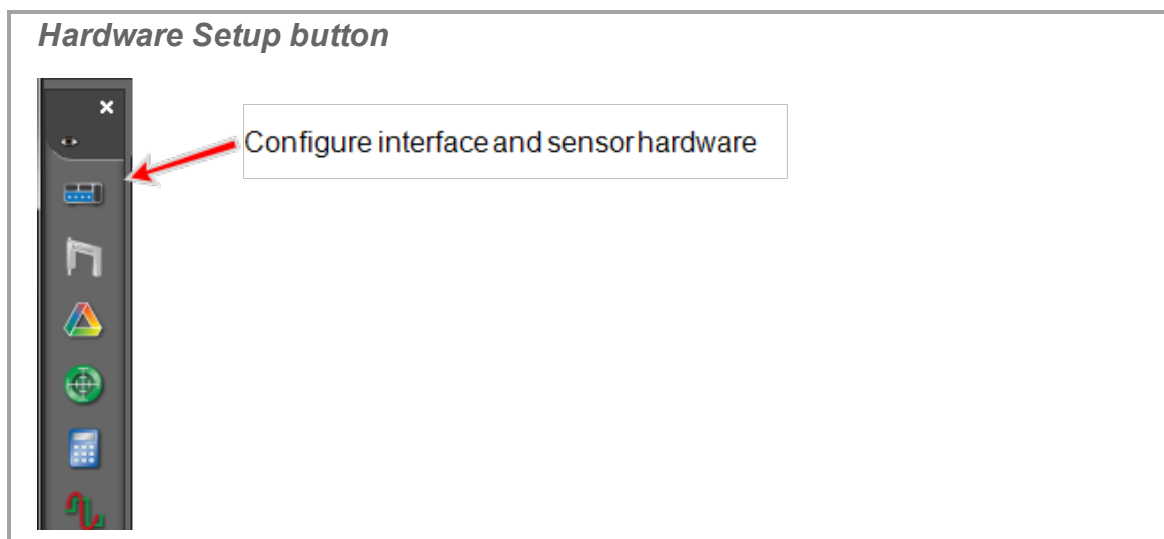

3. Click on the sensor to zero.
4. Click **Properties**  in the lower right corner of the Hardware Setup window.
  - ➔ The **Properties** window opens.
5. Select option:


➔ These options will not appear for sensors that do not have these capabilities.


Option	Description
<b>Zero Sensor Measurements at Start</b>	Zero this sensor's measurements each time a new run is recorded.
<b>Zero Sensor Now</b>	Zero this sensor's measurements when the button is clicked.

6. Click **OK**.

## How do I switch the sign on a sensor?

### Method 1: configuring sensor (that supports sign change)

1. Click Hardware Setup in the Tools Palette: .

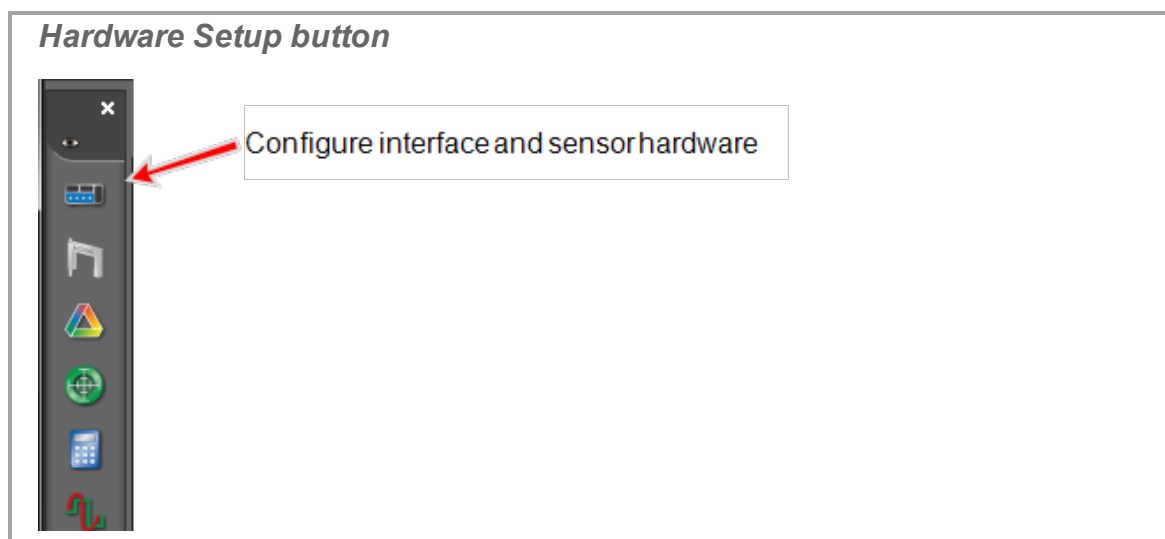

2. Click on the sensor to Change Sign.

3. Click **Properties**  in the lower right corner of the Hardware Setup window.

4. The **Properties** window opens.

5. Click **Change Sign** .

➔ This option will not appear for sensors that do not have this capability.

6. Click **OK**.


➔ The sensor automatically switches the sign when recording data.


### Method 2: using QuickCalc (any sensor)

1. Create a Table. ([See "Adding a display": 56](#))

2. Click **<Select Measurement>** and select the measurement to change sign.

3. Click **QuickCalc**  in the toolbar.

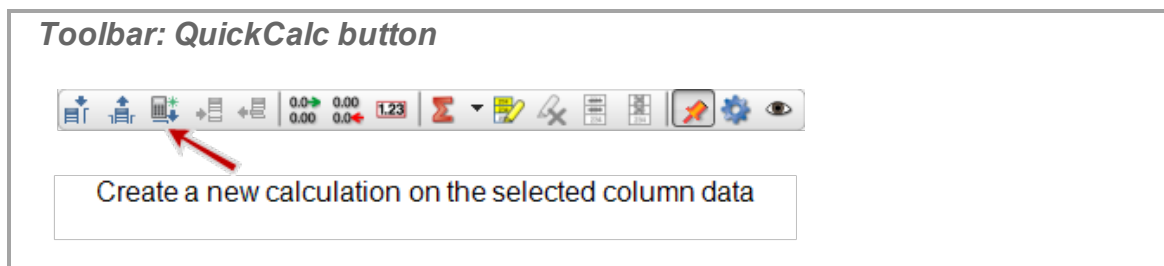


- ➔ A new table column is added with a default calculation that includes the selected measurement.
4. Rename the calculation to something meaningful.
  5. Add a negative sign in the calculation before the measurement.  
Example: **ForceA=- [ Force (N) ]**
  6. Press **Enter**.  
➔ The new column is the opposite sign of the original measurement data.
  7. Click on the units in the new column and change to something meaningful.  
➔ Calculations appear on the **Calculation Data Summary** tab () in **Data Summary**, and are selectable from any data display. (🔗 See "How do I use my calculation?": 102)

### Method 3: using Calculator (any sensor)

Use the Calculator to create a calculation that changes the sign. (🔗 See "How do I create a calculation? ": 99)


Example: **ForceA=- [ Force (N) ]**


- ➔ Calculations appear on the **Calculation Data Summary** tab () in **Data Summary**, and are selectable from any data display. (🔗 See "How do I use my calculation?": 102)

## How do I calibrate a sensor?

PASCO Capstone includes calibration wizards for many sensors.

### Procedure

1. Connect the sensor to PASCO Capstone. (🔗 See "Connecting a sensor": 27)
2. Click **Calibration**  in the Tools Palette.

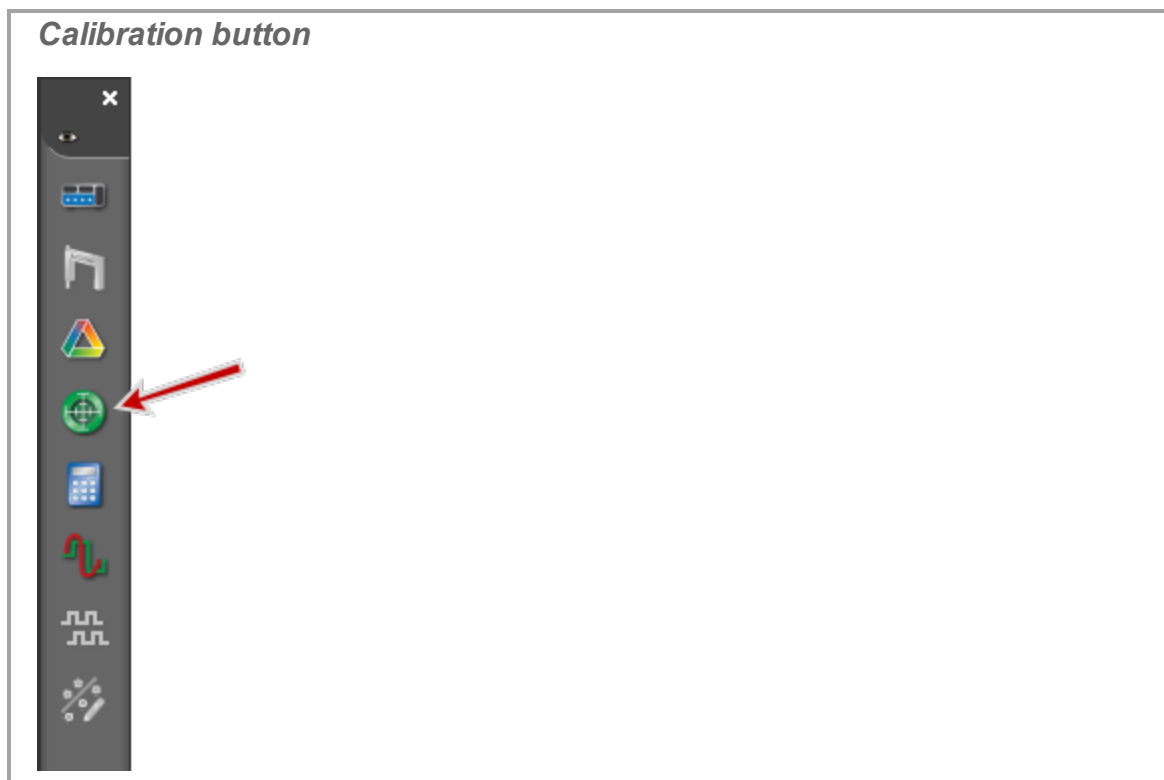

3. Follow the instructions within Capstone.

➔ Sensors that do not have a wizard in Capstone will not be listed.

For detailed procedures, see:

"How do I calibrate a motion sensor?": 95


"How do I calibrate a pH sensor?": 97



"How do I calibrate a drop counter?": 96

"Can multiple measurements be calibrated simultaneously?": 175

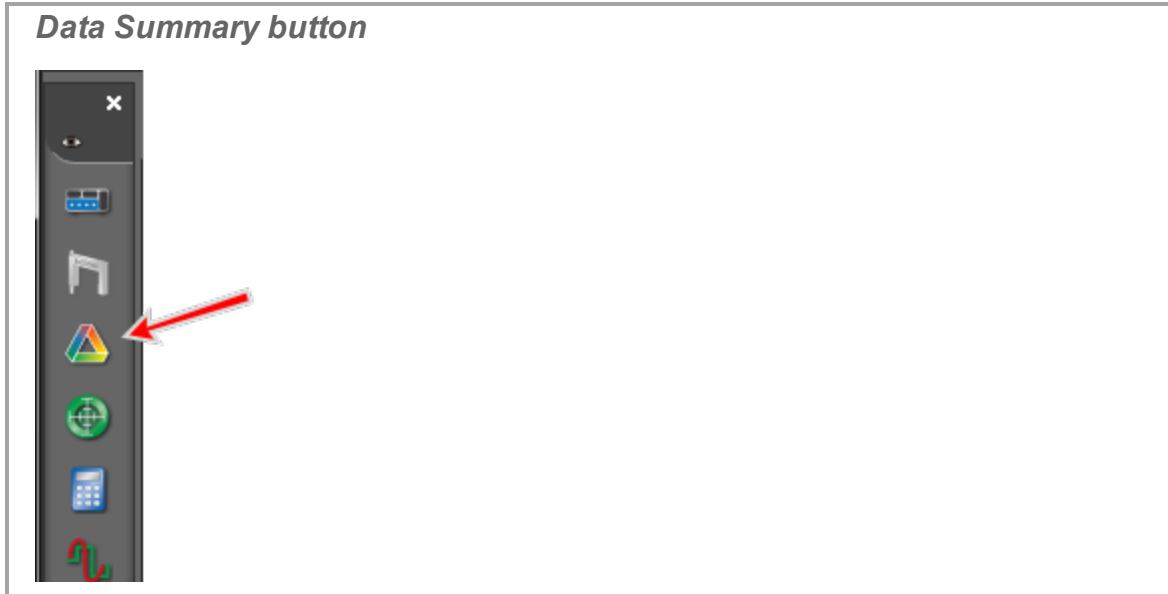

## How do I calibrate a motion sensor?

1. Connect the Motion Sensor to PASCO Capstone. See "[Connecting a sensor](#)": 27.
2. Place a marker at one meter from the face of the motion sensor.
3. Record the distance to the marker reported by the motion sensor.
4. Click **Data Summary** in the Tools Palette.





5. Click the **Sensor Data Summary** tab .
6. Click on the **Motion Sensor** properties icon .
7. Divide the value in the **Speed of Sound (m/s)** field by the recorded distance in meters (from step 3).
8. Enter the calculated value (from step 7) in the **Speed of Sound (m/s)** field.
9. Click **OK**.

## How do I calibrate a drop counter?

1. Connect the Drop Counter to PASCO Capstone. ( See "[Connecting a sensor](#)": 27)
2. Measure the volume from the Drop Counter for 10 drops.
3. Calculate the volume per drop (divide volume from step 2 by 10).
4. Click **Data Summary**  in the Tools Palette.



5. Click the **Sensor Data Summary** tab .
6. Click on the **Drop Counter** properties icon .
7. Enter the calculated value (from step 3) in the **Drop Volume (mL)** field.
8. Click **OK**.



## How do I calibrate a pH sensor?


**Note:** Simultaneous calibration is not recommended for pH, ISE, ORP, Conductivity, stainless-steel temperature sensors without insulating sleeves, or Dissolved Oxygen sensors because these sensors will not read correctly if the probes are placed in the same solution simultaneously.

### Equipment Required

Two aqueous pH buffer standards that bracket the pH values of interest.

### Procedure

1. Connect the pH Sensor to PASCO Capstone. ( See "[Connecting a sensor](#)": 27)
2. Click **Calibration**  in the Tools Palette.

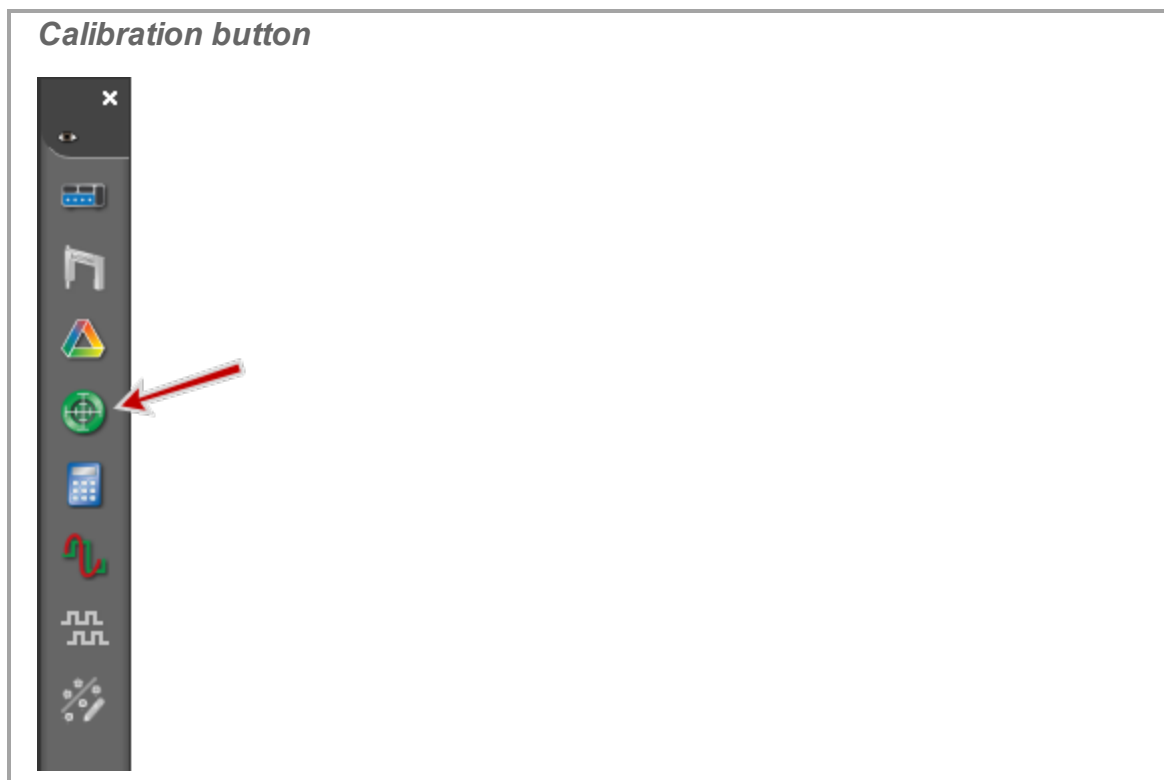

3. Choose the probes you would like to calibrate now: **pHMeasurement**
4. Click **Next**.
5. Choose the type of calibration you would like to perform: **Two Standards (2 point)**.
6. Click **Next**.
7. Enter the first known pH buffer value in the **Standard Value** text box.
8. Rinse and dry the pH Sensor probe.
9. Place the probe in the corresponding buffer solution.
10. Stir the probe until the **Current Value** stabilizes.
11. Click the **Set Current Value to Standard Value** button.
12. Click **Next**.
13. Enter the second known pH buffer value in the **Standard Value** text box.
14. Rinse and dry off the probe.
15. Place the probe in the second buffer solution.
16. Stir the probe until the value **Current Value** stabilizes.
17. Click the **Set Current Value to Standard Value** button.
18. Click **Next**.
19. Click **Finish**.

Repeat the calibration procedure for other pH sensors using different probes.

## How do I... Calculations

"How do I create a calculation? ": 99



"How do I use my calculation?": 102

"How do I select a measurement in the calculator?": 103


"How do I see the result of a calculation?": 103

"How do I reorder calculations?": 105

### How do I create a calculation?

**Note:** Calculations appear on the **Calculation Data Summary** tab () in **Data Summary**, and are selectable from any data display. ( See "How do I use my calculation?": 102)


#### Tips

- Capstone supports simple to highly-complex calculations using a combination of math functions, measurements, user-defined data, constants, and text labels including symbols.
- Use either the keyboard or the **Calculator** keypad for numbers and simple operators.
- Add parenthesis to explicitly define order of operation.
- Multiplication is implied with adjacent arguments. Insert the **\*** symbol to explicitly denote multiplication.
- Type: **[** to quickly insert measurements (or click **Insert Data** .

#### Procedure

1. Click **Calculator**  in the Tools Palette.



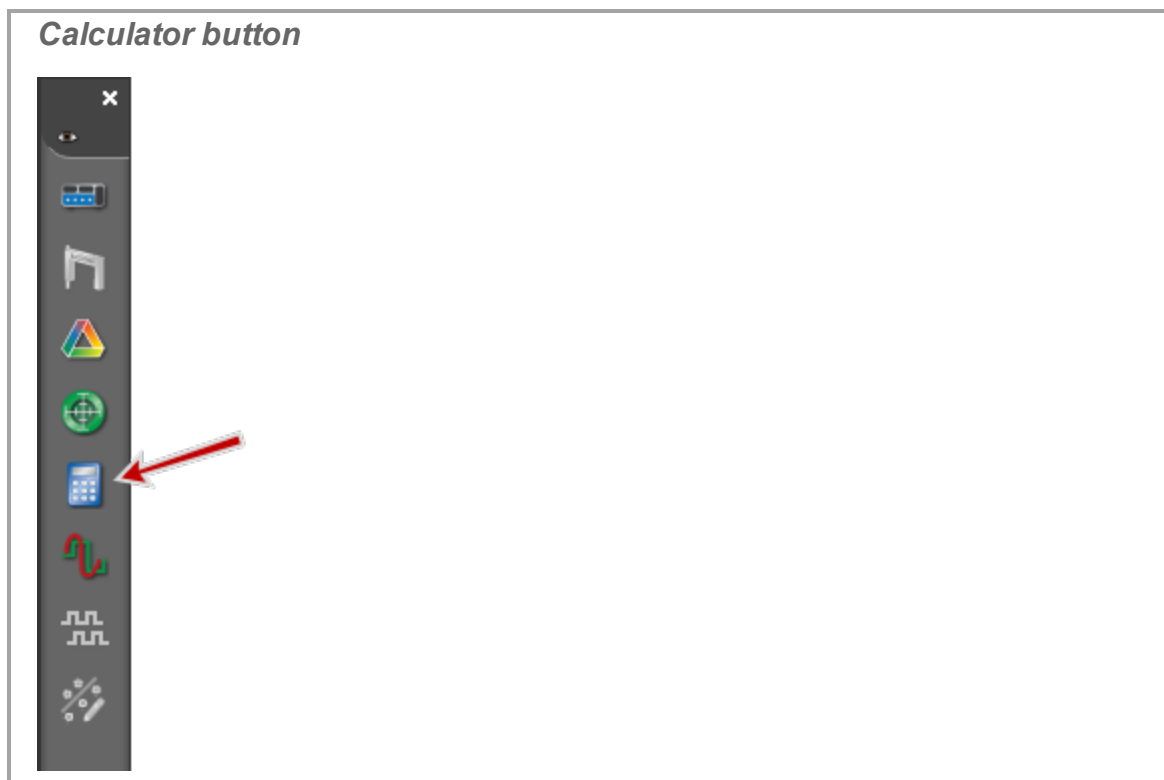

➔ The **Calculator** dialog opens.

2. Use a combination of [measurements](#), text, calculator functions, and numerals to create a calculation.

➔ The calculation is displayed **Calculations** list.

➔ Invalid or incomplete calculations are indicated with red font.

[See "How do I select a measurement in the calculator?": 103](#)

[See "How do I see the result of a calculation?": 103](#)

[See "How do I reorder calculations?": 105](#)

[See "How do I use my calculation?": 102](#)

**Example Procedure:**  $x = \frac{1}{2}at^2 + vt + x_0$

Define the following equation:

$$x = \frac{1}{2}at^2 + vt + x_0$$


**Where:**



x = Position (calculated)



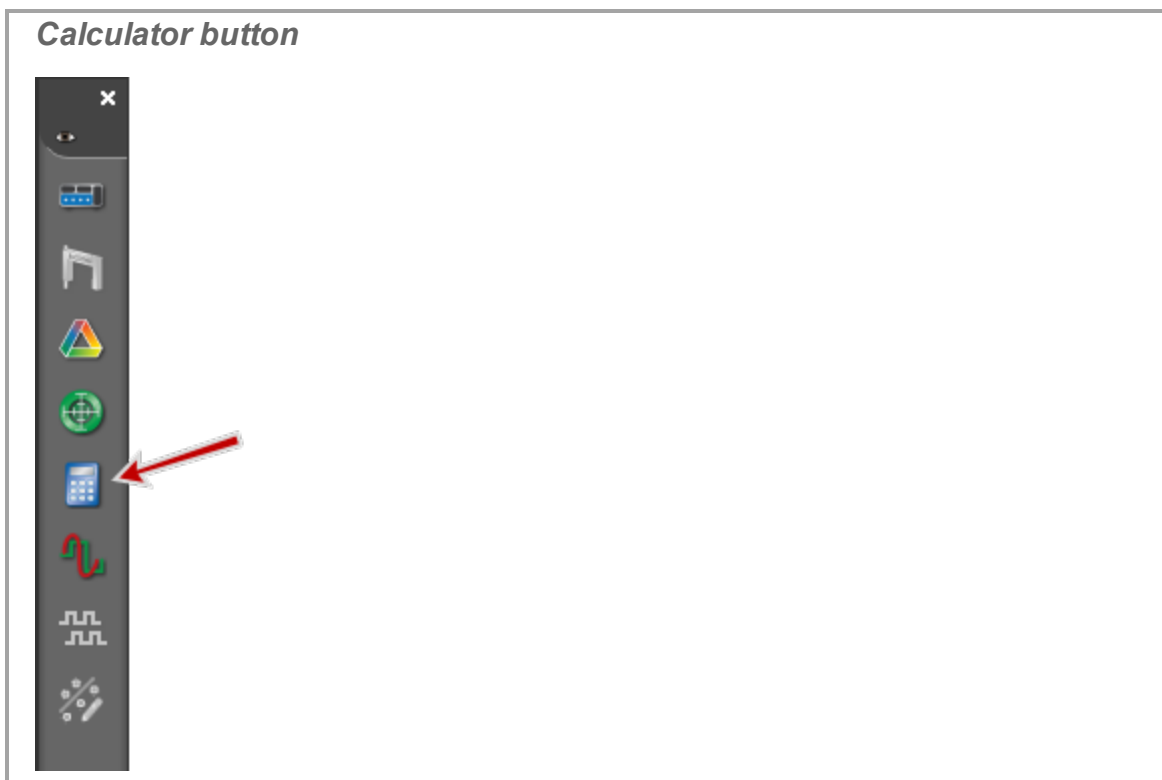

a = Acceleration (measured)

- v = Velocity (measured)
- t = Time (measured)
- $x_0$  = initial position (user-created constant)

1. Click **Calculator**  in the Tools Palette.



- ➔ The **Calculator** dialog opens.
2. Click in a new line of **Calculations**.
3. Create the constant  $x_0$ :
  - a. Type: **x**
  - b. Click  **$x_0$** , select **0**
  - c. Type: **=**
  - d. Type: **0**
4. Click in a new line of **Calculations**
5. Type: **x=1/2 (**
6. Type: **[** (or click **Insert Data** ).
  - ➔ A pop-up list opens.
7. Select the appropriate Acceleration measurement from the drop down list.
8. Type: **[** (or click **Insert Data** .

9. Select the Time measurement from the drop down list.
10. Click  $x^2$ .
11. Type: + (or click + )
12. Type: [ (or click **Insert Data** ).
13. Select the appropriate Velocity measurement from the drop down list.
14. Type [ (or click **Insert Data** ).
15. Select the **Time** measurement from the drop down list.
16. Type: + (or click + )
17. Type: [ (or click **Insert Data** ).
18. Select the **Constants** (under **Equations/Constants**) measurement from the drop down list.
19. Select  $x_0$ .
20. Click in the **Units** field.
21. Type: m and select  $m/s^2$  from the list.
22. Press **Enter** or **Accept** .

## Example Calculations

$Constant1=2*3$



$Calc1=[Acceleration(m/s^2)]^2$


$r^2=sin([Angle(rad)])$

$\phi=integral([Force(N)],[Time(s)])$

$Resultant\ Acceleration=sqrt(( [Acceleration-x(m/s^2)]^2+[Acceleration-y(m/s^2)]^2+[Acceleration-z(m/s^2)]^2))$

## How do I use my calculation?


**Note:** Calculations appear on the **Calculation Data Summary** tab () in **Data Summary**, and are selectable from any data display. ( See "How do I use my calculation?": 102)

1. Create a calculation. ( See "How do I create a calculation? ": 99)
2. Click **<Select Measurement>** in any display.

3. Select the calculation or constant to use under **Equations/Constants**.
4. Select the run to display. (📎 See "How do I choose one or more runs to display?": 79)

## How do I select a measurement in the calculator?

### Method 1

1. Position cursor for placement of measurement.
2. Click **Insert Data** .
  - ➡ A pop-up list opens.
3. Select the measurement from the list.


### Method 2

1. Right-click anywhere within the calculation to place the measurement.
  - ➡ A pop-up menu opens.
2. Select **Insert Data** from the menu.
3. Select the measurement from the list.

### Method 3

1. Position cursor for placement of measurement.
2. Enter a left-bracket “[ ”.
  - ➡ A pop-up list opens.
3. Select from the measurement from the list.


## How do I see the result of a calculation?

**Note:** Calculations appear on the **Calculation Data Summary** tab () in **Data Summary**, and are selectable from any data display. (📎 See "How do I use my calculation?": 102)

## Constants Only

If the calculation uses constants (no measurements) such that the calculation evaluates to a single number, then the result of the calculation is displayed in the result status line.

*Result status line*

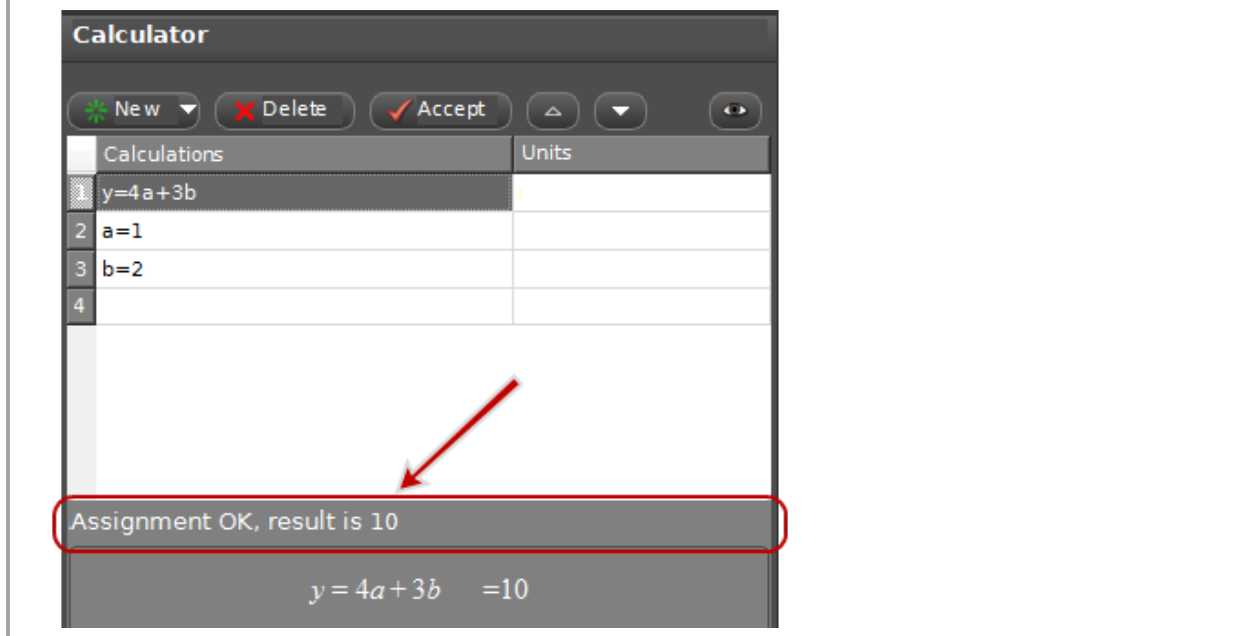

Constants can also be viewed in a table or other display, see below.

### Includes Variables (measurements)

The results are calculated using MathML and are associated with data runs. Calculated data can be viewed in a table or other display.

1. Create a calculation. (📎 See "How do I create a calculation? ": 99)
2. Click **<Select Measurement>** in any display.
3. Select the calculation or constant to use under **Equations/Constants**.
4. Select the run to display. (📎 See "How do I choose one or more runs to display?": 79)

### Example Calculations

$$\text{Constant1}=2*3$$

$$\text{Calc1}=[\text{Acceleration (m/s}^2)]^2$$

$$\Gamma^2=\sin ([\text{Angle (rad)}])$$

$$\Phi=\text{integral} ([\text{Force (N)}], [\text{Time (s)}])$$


Resultant Acceleration=sqrt(([Acceleration- x(m/s<sup>2</sup>)]<sup>2</sup>+ [Acceleration- y(m/s<sup>2</sup>)]<sup>2</sup>+ [Acceleration- z(m/s<sup>2</sup>)]<sup>2</sup>))

## How do I reorder calculations?

### Method 1

1. Click and drag the calculation that you would like to move.
2. When the colored insert line is at the desired insertion point, release the mouse button.

### Method 2

1. Click the calculation that you would like to move to select it.
2. Use the arrow keys  in the **Calculator** window.

## How do I... Plots

"How do I plot multiple measurements on separate plot areas with shared X-axis?": 107

"How do I zoom to a selection?": 108

"How do I pan and scale?": 110

"How do I select active data?": 111

"How do I plot multiple measurements on same plot with multiple Y-axes?": 113

"How do I plot multiple measurements on same plot with same Y-axes?": 114

"How do I apply a curve fit to a portion of the data?": 117

"How do I change curve fit parameters?": 118

"How do I adjust the weights for the weighted linear fit?": 119


"How do I lock and unlock the axes of a graph?": 121

"How do I change layer order of a display?": 123

## How do I plot multiple measurements on separate plot areas with shared X-axis?

### Attributes of multiple plot areas

- 1 The plots share the same x-axis measurement.
- 2 The y-axes are independent.
- 3 The plots from the different measurements do not overlap.

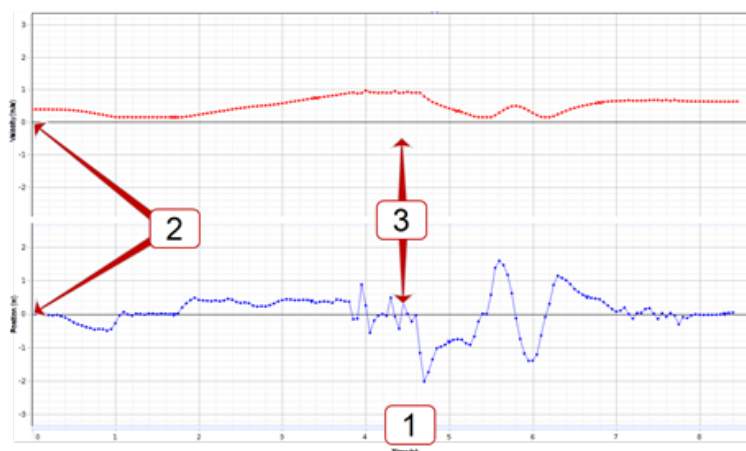


[See "How do I plot multiple measurements on same plot with multiple Y-axes?": 113](#)

[See "How do I plot multiple measurements on same plot with same Y-axes?": 114](#)

### Procedure

1. Click **Add new plot area to the Graph display**  in the toolbar.

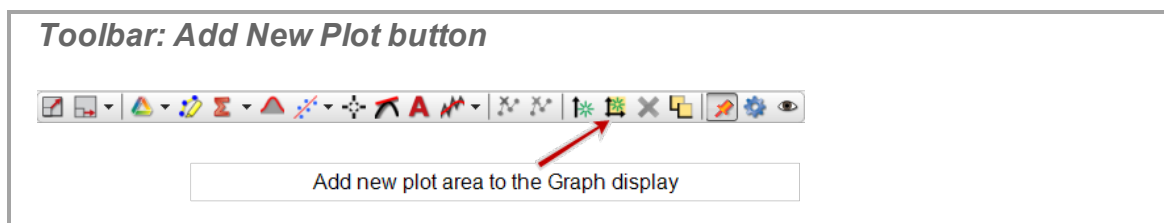

2. Click **<Select Measurement>** on the plot and choose the measurement for the vertical axis.



- To delete a plot area and its corresponding plots, click on the plot area and click **Remove active element, axis, or plot area** ✖ .

**Note:** You will not be allowed to have fewer than one plot area. Use **Delete Selected Display** ✖ in main toolbar to remove the display. (Scope display does not allow multiple plots.)

## How do I zoom to a selection?

### Zoom in

- If multiple runs are visible, click on the symbol corresponding to the plot to zoom in the legend.

**Legend: selecting the active plot**

Symbol and color in legend matches plot symbol and color. Click legend

- to make active. (The active plot is shaded in legend and darkest in display.)
- Click in the upper left corner of the legend to clear all active runs.


- Click **Data Highlighter**  in the toolbar.

**Toolbar: Highlight Data button**

Apply selected curve fits to active data | Select curve fits to be displayed

- ➔ A rectangle appears on the graph that is the same color as the target plot.


3. Click and drag anywhere within the highlighter to move it.
4. Click and drag the handles to resize the highlighter rectangle to the area to zoom.
  - ➔ Selected data is highlighted in yellow.
5. Click **Scale to Fit** in the toolbar.

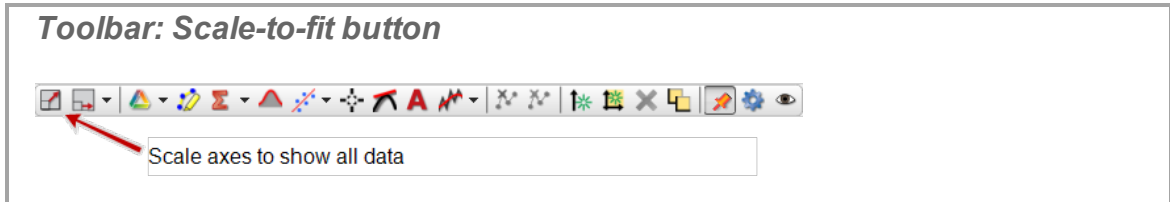

- ➔ The plot zooms to the selected section.
6. Optional: Remove the highlight rectangle:
    - a. Click the highlight rectangle
    - b. Press **Delete** to remove the rectangle


### Zoom out

1. Remove the highlight rectangle:
  - a. Click the highlight rectangle
  - b. Press **Delete** to remove the rectangle
2. Click **Scale to Fit** in the toolbar.

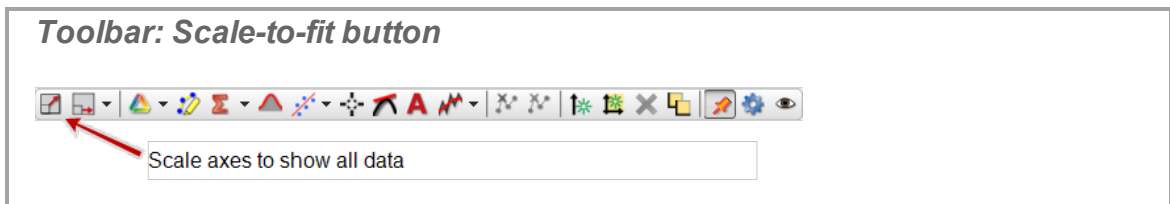

- ➔ The plot zooms out to include all data.

## How do I pan and scale?

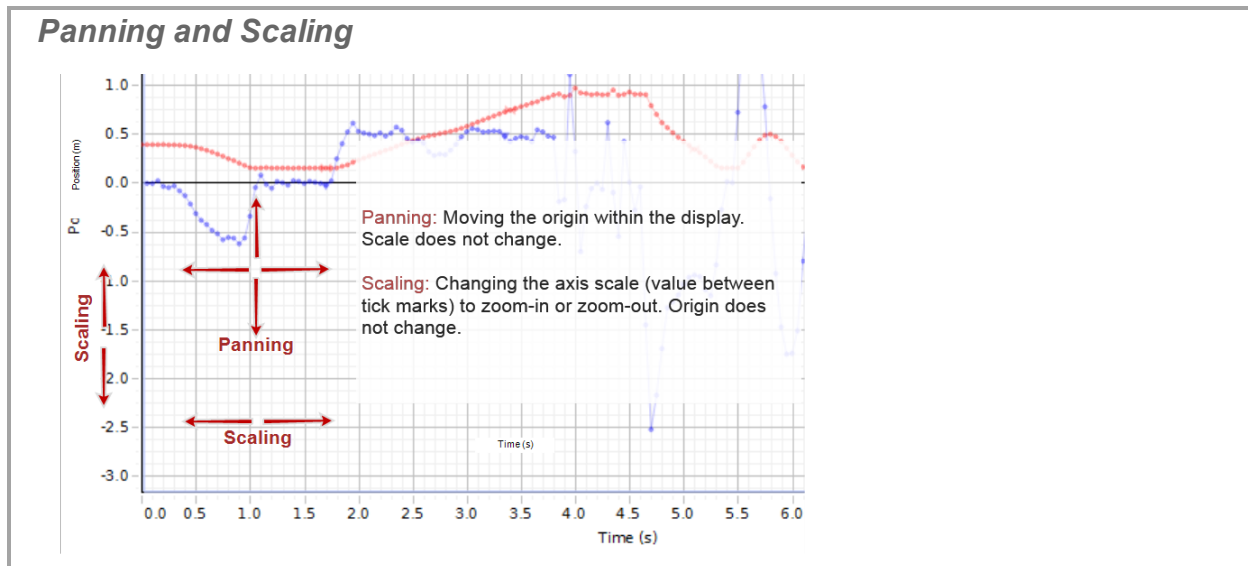

**Note:** If the plot does not pan or scale: see "How do I lock and unlock the axes of a graph?": 121.

### **Panning** *Moving the origin within the display. Scale does not change.*

1. Click within the plot area (except for the legend) and drag the plot in any direction.
2. Release the mouse button when the origin is in the desired location.

### **Scaling** *Changing the axis scale (value between tick marks) to zoom-in or zoom-out. Origin does not change.*

#### **Click and Drag Method**

1. Click on an axis to scale.
2. Drag away from origin to decrease scale (zoom-in).
3. Drag toward the origin to increase scale (zoom-out).

#### **Scroll Wheel Method**

1. Move the cursor to a location on an axis to scale.
2. Scroll back to decrease scale (zoom-in).
3. Scroll forward to increase scale (zoom-out).

↻ Scaling is performed on the axis at the point of the cursor.

💡 Tip<sup>1</sup>

### Zooming around a point *Changing the scale and origin around a single point.*

1. Move the cursor to a location on a point to remain stationary (except for the legend).
2. Scroll back to decrease scale (zoom-in).
3. Scroll forward to increase scale (zoom-out).

### How do I select active data?

By default, the last selected measurement of the last run is the active run.

To make another plot active:

- Click on the symbol corresponding to that plot in the legend (1 in the example)).

**Legend: selecting the active plot**

Symbol and color in legend matches plot symbol and color. Click legend (1) to make active. (The active plot is shaded in legend and darkest in display.)

(2) Click in the upper left corner of the legend to clear all active runs.

The image shows a plot window with a grid. Three data series are plotted: a blue line with circles, an orange line with triangles, and a pink line with squares. A legend is located in the upper right corner of the plot area. The legend contains three entries, each with a colored symbol and a text label. The first entry has a pink triangle and is highlighted with a red box labeled '1'. A red arrow points from this box to the pink triangle in the legend. The second entry has an orange triangle. The third entry has a blue triangle. A red box labeled '2' is in the upper left corner of the legend, with a red arrow pointing to it.

To clear all active runs:


<sup>1</sup>◇ Place the cursor at the origin to maintain the origin point or use click and drag method. ◇ Place the cursor at the intersection of two axis to scale simultaneously.

- Click in the upper left corner of the legend (2) in the example).

**Legend: selecting the active plot**

1 Symbol and color in legend matches plot symbol and color. Click legend to make active. (The active plot is shaded in legend and darkest in display.)

2 Click in the upper left corner of the legend to clear all active runs.

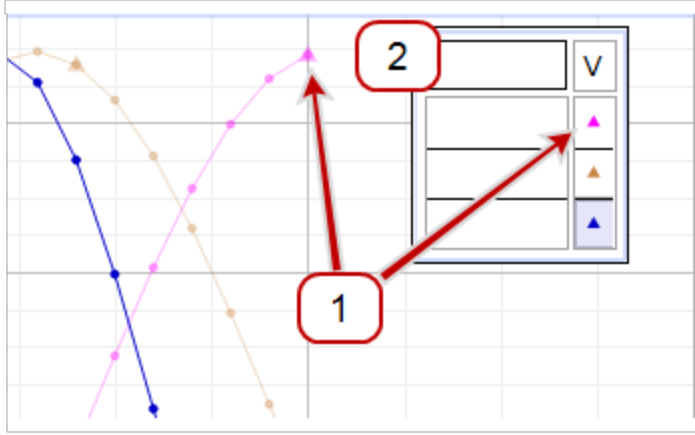


**Toolbars (including data functions):**

- ➡ Toolbars are active when a single display is active, and will only apply to the active display.
- ➡ For displays with multiple plots, the toolbar actions apply to the active plot within the display.

## How do I plot multiple measurements on same plot with multiple Y-axes?

### Attributes of multiple Y-axes


- 1 Measurements scale independently( except for the rescaling required to make the alignment of the tick mark values on each side of the graph match up).
- 2 The order of the axes may be changed once multiple axes are created.
- 3 Plots from different measurements use the same plot area and may overlap.

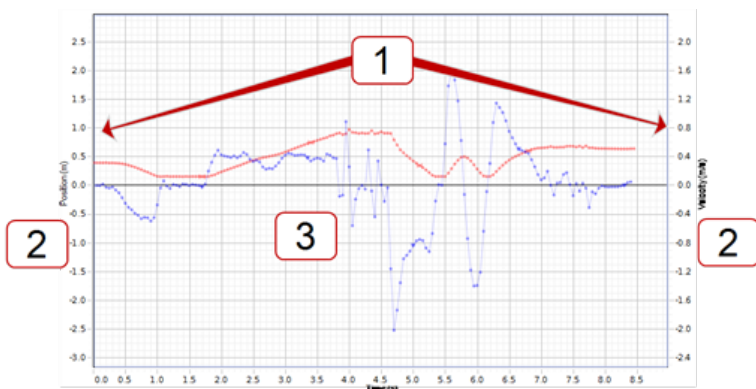


See "How do I plot multiple measurements on separate plot areas with shared X-axis?": 107

See "How do I plot multiple measurements on same plot with same Y-axes?": 114


### Procedure


1. Create a graph display. (See "Adding a display ": 56)
2. Click **Add new y-axis to active plot area** \*. (Click multiple times to add multiple axes.)

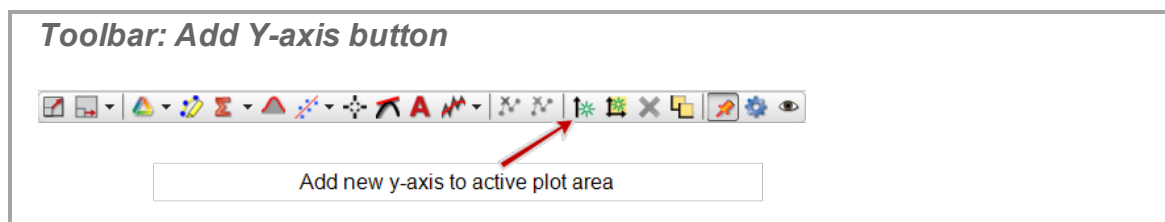

3. Click **<Select Measurement>** and choose the measurements for each vertical axis.
4. Optional: reorder axes:

- a. Click **Allow rearrangement of axes and plot areas**  in the toolbar.

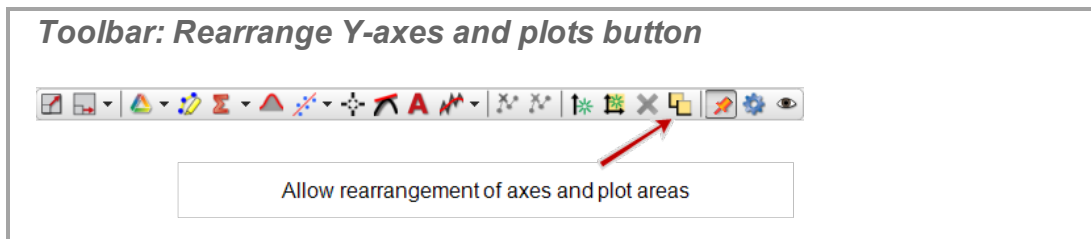


- b. Drag and drop axes one at a time.
- ➡ Axes can reside on the left, right, or a combination of both sides of the display.
- c. Click **Allow rearrangement of axes and plot areas** .
- ➡ Exits re-arrangement mode.

### Remove an axis

Click on axis to remove. (Must not be in re-arrangement mode.)

Click **Remove active element, axis, or plot area** .




**Note:** You will not be allowed to have fewer than one plot area. Use **Delete Selected Display**  in main toolbar to remove the display. (Scope display does not allow multiple plots.)

### How do I plot multiple measurements on same plot with same Y-axes?

PASCO Capstone allows one measurement on the x (independent variable) axis and multiple measurements on the y (independent variable) axis. Y-axis labels can be placed on either side. If the measurements share the same measurement type, then the measurements can be plotted with a shared Y-axis.

**Attributes of multiple measurements with a shared Y-axis**


- 1 Measurements must share the same measurement type. The shared generic Measurement Type name is displayed on the Y-axis.
- 2 Measurements scale together.
- 3 Plots from different measurements use the same plot area and may overlap.

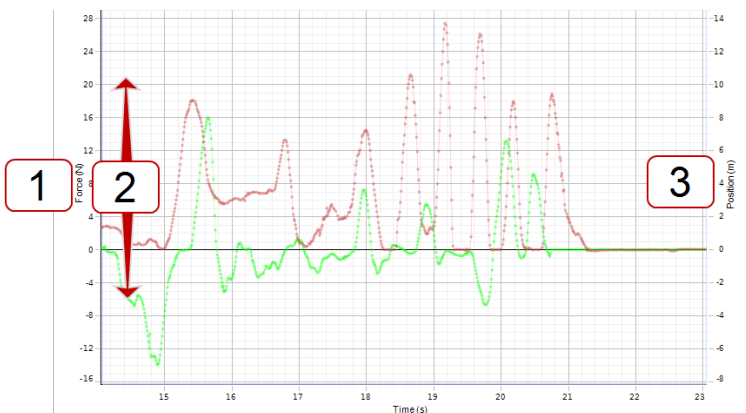

[See "How do I plot multiple measurements on separate plot areas with shared X-axis?": 107](#)

[See "How do I plot multiple measurements on same plot with multiple Y-axes?": 113](#)


**Procedure**

1. Create a graph display. ([See "Adding a display ": 56](#))
2. Click **<Select Measurement>** and select one of the measurements to display on the vertical Y-axis.
3. Click **Data Summary**  in the Tools Palette.




**Data Summary button**



4. Click the **Sensor Data** tab .
  5. Click and drag another measurement of the *same measurement type* to the graph display.
  6. Repeat for additional sensors of same type.
- ➔ The shared generic Measurement Type name is displayed on the Y-axis.

**Multiple-Y-axes**

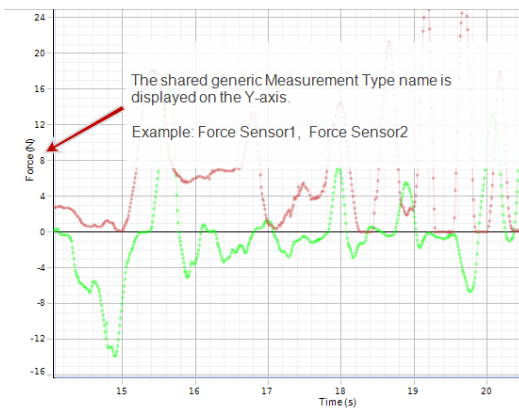

## How do I apply a curve fit to a portion of the data?

1. Select the plot to fit by clicking on its corresponding symbol in the legend.


**Legend: selecting the active plot**

- 1 Symbol and color in legend matches plot symbol and color. Click legend to make active. (The active plot is shaded in legend and darkest in display.)
- 2 Click in the upper left corner of the legend to clear all active runs.

2. Click **Data Highlighter**  in the toolbar.

**Toolbar: Highlight Data button**

Apply selected curve fits to active data | Select curve fits to be displayed

- ➡ A rectangle appears on the graph that is the same color as the target plot.
3. Click and drag anywhere within the highlighter to move it.
4. Click and drag the handles to resize the highlighter rectangle.
- ➡ Selected data is highlighted in yellow.
5. Click the down arrow (▼) on **Curve Fit** .
6. Select the curve fit that you wish to apply to the selected data from the pull-down menu.

**Note:** Move the highlighter rectangle to see fits for different regions. (Capstone will not simultaneously perform curve fits on two or more different regions of the same data set.)

## How do I change curve fit parameters?

PASCO Capstone supports modifying curve fit parameters for high-precision fits. (Initial curve fits are usually automatic and accurate.)

### Procedure


1. If not visible, show the **Curve Fit Editor**  in the Toolbar. (See "How do I show, hide, and lock the palettes?": 89)





2. Select the plot to fit by clicking on its corresponding symbol in the legend.

**Legend: selecting the active plot**

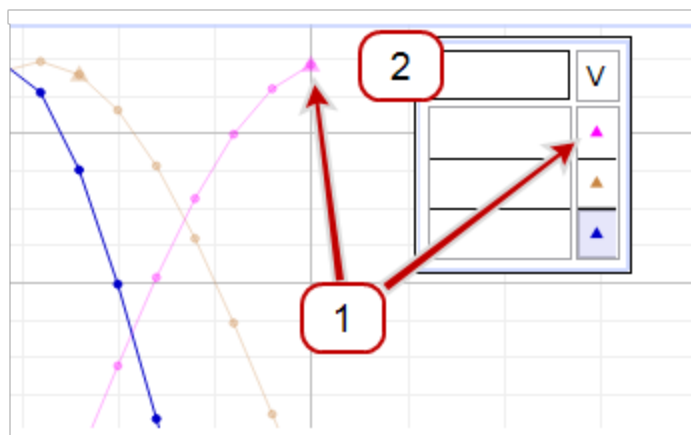- 1 Symbol and color in legend matches plot symbol and color. Click legend to make active. (The active plot is shaded in legend and darkest in display.)
- 2 Click in the upper left corner of the legend to clear all active runs.






- 3. Click **Curve Fit Editor**  in the Toolbar.
- 4. Enter the value to modify in the **Initial Guess** box of the corresponding parameter.
- 5. Click the **Lock** check box () next to parameter.
  - ↻ Locks the manual entry. Capstone will not overwrite.
- 6. Repeat for each parameter to modify.
- 7. Click the **Update Fit** button.
  - ↻ The curve fit in the graph uses the manually entered values.
  - ↻ Repeat steps to iteratively improve the curve fit.
  - ↻ To reset to default settings and unlock parameter values, click **Reset**.
- 8. Click **Curve Fit Editor**  in the Toolbar to close the editor.

### How do I adjust the weights for the weighted linear fit?

The weights for the weighted linear fit come from the errors in the data set. To define the error associated with each data value:

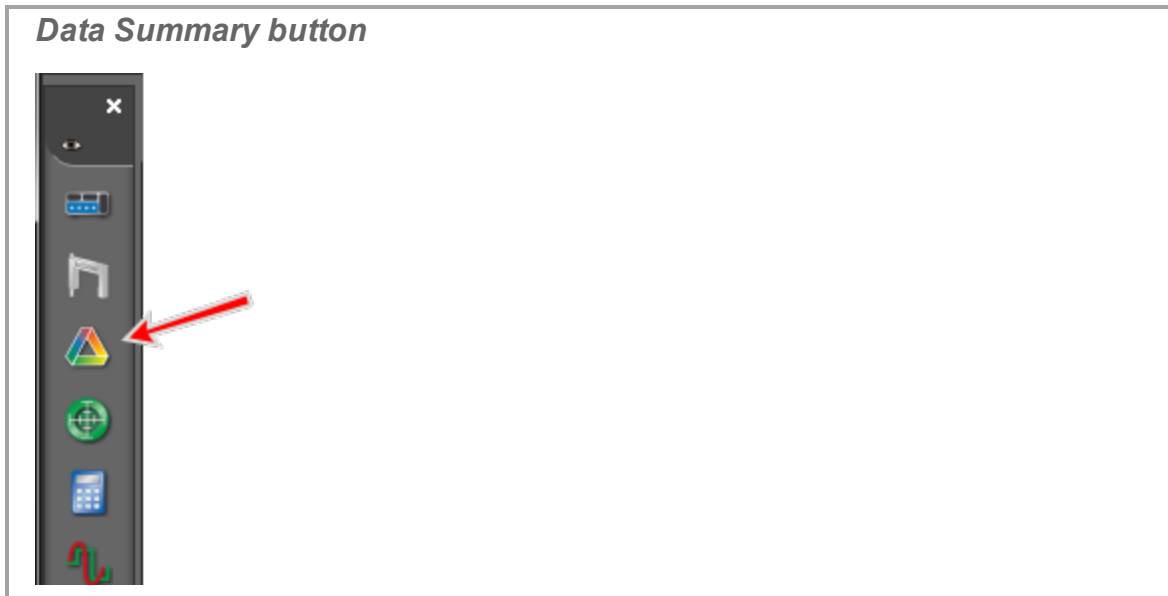1. Click **Data Summary**  in the Tools Palette.




2. Click the **Sensor Data** tab .
3. Click the measurement to change.
4. Click **Properties**  that appears to the right of the measurement.
  - ➔ The **Properties** window opens.
5. Click the **Errors** section.
6. Define the error type and values.

Error Type	Description
<b>Default Absolute Error</b>	Specifies the default absolute error value for this measurement and is used when Default Errors are in effect.
<b>Default Minimum Display Value</b>	Sets the default minimum scale value for this measurement when it is initially selected in some data displays.
<b>Default Maximum Display Value</b>	Sets the default maximum scale value for this measurement when it is initially selected in some data displays.
<b>Error Bar Type</b>	Select the type of error bar to display. Choose Default Error to set the length of the error bars equal to the default absolute error of the data. Choose Percent of Value to set the length of the error bars equal to the specified percentage of each data point value. Choose Fixed Range to set the length of the error bars to equal specific val-

Error Type	Description
	ues. Choose Measurement to use a User Entered Data Set or other Measurement to define error values for individual data points.
<b>Reset</b>	Resets the error settings for this measurement to defaults.


## How do I lock and unlock the axes of a graph?

PASCO Capstone has several options for locking the axes of a graph from Panning and Scaling.

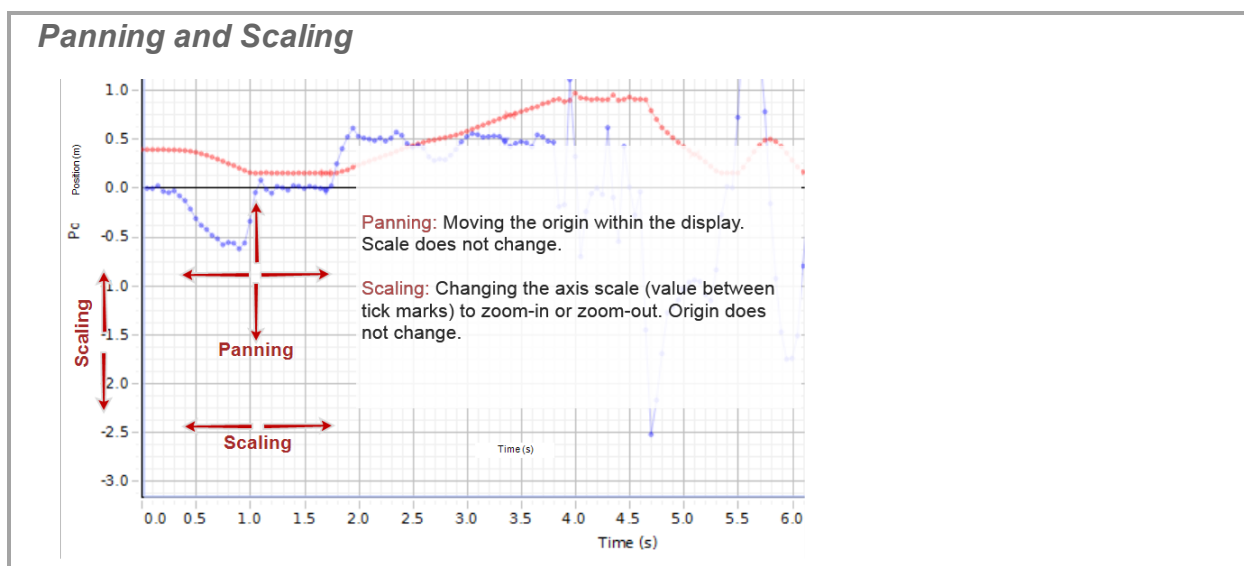


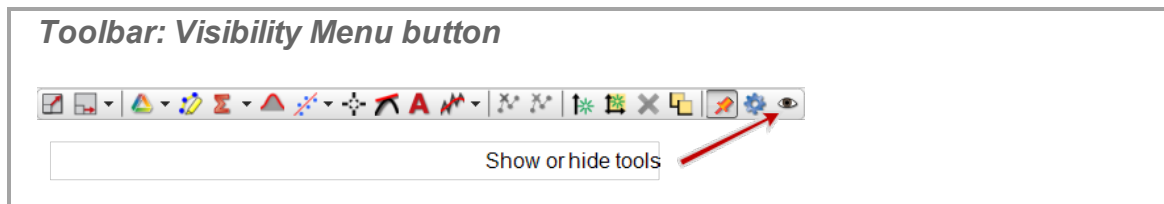
## Lock panning or scaling from both user *and* data changes


1. Deselect **Scaling Method**  in the Display Toolbar.

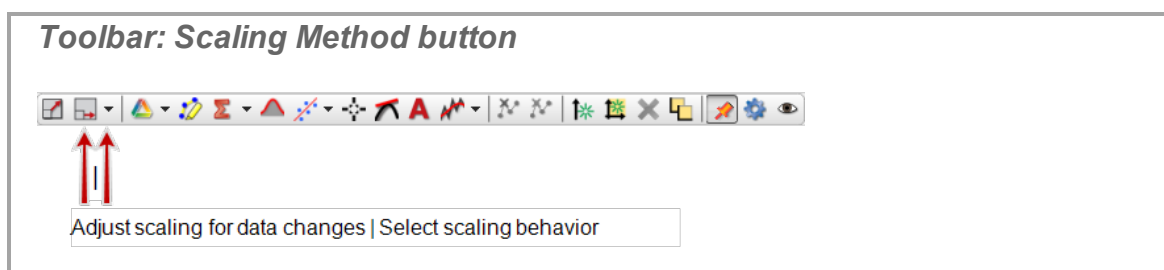

- ⌚ Capstone will not scale to the data.
2. Right click on an axis and select one or both:
  - **Lock Axis Panning**
  - **Lock Axis Scale**


- ➡ Users cannot manually pan or scale.
- 3. Optional: repeat for additional axes.
- 4. Click **Show or hide tools**  in the Display Toolbar.



- 5. Deselect **Scale to Fit**  in the list.
- ➡ Users cannot scale to fit.
- 6. Click anywhere in Capstone to close **Show or hide tools**.


### Lock scaling from data changes only


Deselect **Scaling Method**  in the Display Toolbar.

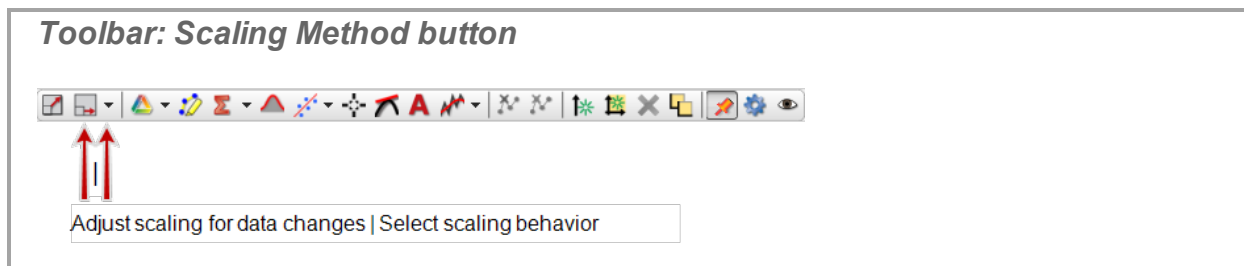


➡ Capstone will not scale to the data.


### Lock panning or scaling from user changes only

- 1. Right click on an axis and select one or both:
  - **Lock Axis Panning**
  - **Lock Axis Scale**
- ➡ Users cannot manually pan or scale.
- 2. Optional: repeat for additional axes.
- 3. Click **Show or hide tools**  in the Display Toolbar.

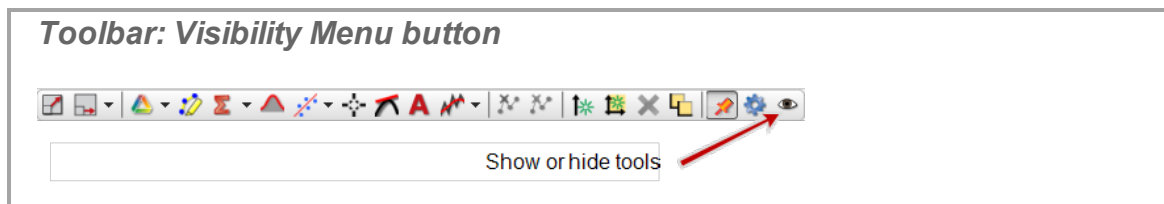

4. Deselect **Scale to Fit**  in the list.
  - ➔ Users cannot scale to fit.
5. Click anywhere in Capstone to close **Show or hide tools**.

## How do I change layer order of a display?

1. Click on the display to make it active. ( See "[How do I select active data?](#)": 111)
2. Right-click on the border of the display.
3. Select: **Send To Back** or **Bring To Front**.
  - ➔ Send To Back = places display at the bottom of the stack.
  - ➔ Bring To Front = places display at the top of the stack.



## How do I... Tables

"How do I reorder columns in a table?": 124

"How do I use a table to build data in Keep Mode?": 124


### How do I reorder columns in a table?

1. Click on an empty part of the column header.
2. Drag until the blue column marker indicates the desired location for the column.
3. Release the mouse button.


### How do I use a table to build data in Keep Mode?



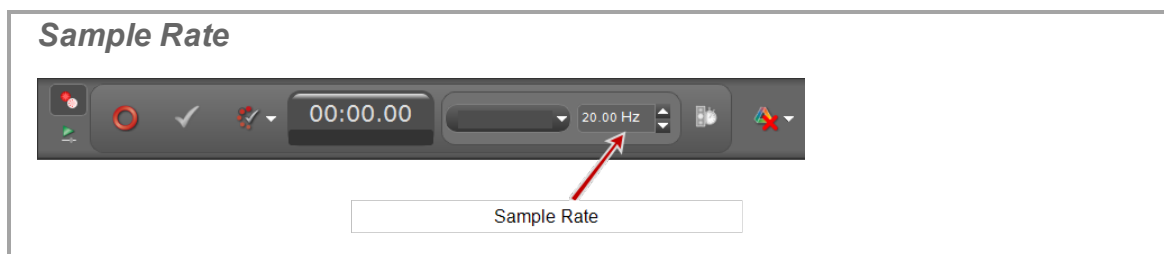

#### Procedure


1. Create a table for the measurement to capture. (See "Adding a display ": 56)
2. Optional: Create User-Entered Data into one column of the table.
3. Click in the first cell to record data.
4. Select **Keep Mode** in the Controls Palette.
5. Adjust the sample rates of each of the measurements so that the sensors are sufficiently responsive for desired data.

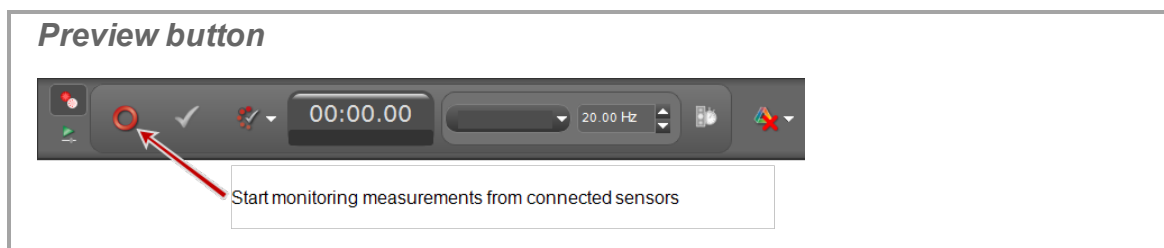




6. Click **Preview** in the Controls Palette.



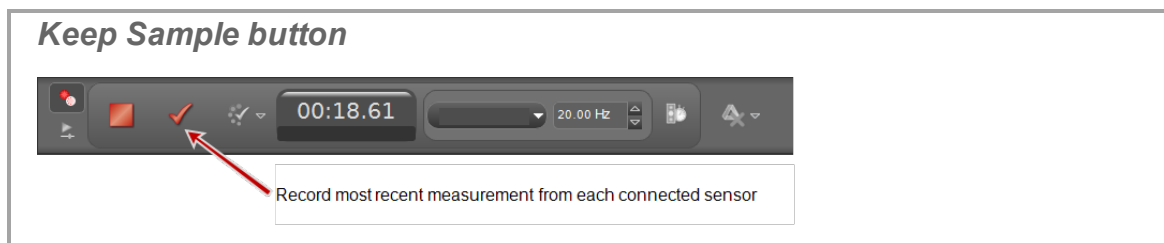

- ➔ Capstone begins monitoring data and displays previews.
  - ➔ **Preview**  toggles to **Stop** .
  - ➔ **Keep Sample**  button is active.
7. Click **Keep Sample**  to record the data currently monitored.
- ➔ One instance of all data is recorded.
  - ➔ The cursor advances to the next table cell.



8. Repeatedly click **Keep Sample**  to record data.
- ➔ One instance of all data is recorded each time **Keep Sample**  is clicked.
  - ➔ The cursor advances to the next table cell on each **Keep Sample**.
9. Optional: Replace values in a table. (See ["Replacing a current value in a table": 125.](#))

**Note:** You can replace data in tables while in preview mode for the run (before clicking **Stop** .

 See ["How do I edit a measurement?": 87](#) to edit data after you click **Stop** .

10. Click **Stop**  in the Controls Palette to stop observing the data.
- ➔ Collected data appear in all displays of the current page.
  - ➔ The data run is listed in the legend for each display.
  - ➔ Recorded data appear on the **Sensor Data Summary** tab () in **Data Summary**, and is selectable from any data display with a default name **Run #%1**, where **%1** is an automatically generated run number.


### Replacing a current value in a table



1. While previewing data, click in the cell to change.
- ➔ The current monitored value is displayed with the existing value.


*Replace data in table cell*



	2.926	
	3.013	
3.062		3.087
<b>1</b>	3.384	<b>2</b>
	3.627	
	3.731	
	3.883	

1 Existing value  
2 New value

Click **Keep Sample**  to replace the existing value with the new value.

2. Click **Keep Sample**  to replace the existing value with the new value.
  - ➔ The cursor advances to the next table cell.
3. Perform one of the following options:
  - Continue to replace values.
  - Click in empty cell to record additional data.
  - Click **Stop**  in the Controls Palette to stop observing the data.

**Note:** You can replace data in tables while in preview mode for the run (before clicking **Stop** .

 See "How do I edit a measurement?": 87 to edit data after you click **Stop** .

## How do I... Text Boxes

"How do I enter subscripts, superscripts, Greek letters, and symbols?": 127

"How do I enter an equation into a text display?": 127

### How do I enter subscripts, superscripts, Greek letters, and symbols?


#### New: enter subscripts, superscripts or symbols

1. Right-click in any area that accepts text such as: text box, text entry, name, or calculator fields.
2. Select one of the following:
  - **Insert Superscript**
  - **Insert Subscript**
  - **Insert Upper Case Greek Letter**
  - **Insert Lower Case Greek Letter**
  - **Insert Special Character**
3. Select the character to insert from the drop-down list.

#### Existing: subscript or superscript selected text (text and text entry boxes only)


1. Highlight existing text.
2. Right-click and select one of the following:
  - **Superscript Selected Text**
  - **Subscript Selected Text**

### How do I enter an equation into a text display?


1. Create the equation using an application capable of exporting [MathML](#) as plain text.
2. Paste into a Text Box or Text Entry Box in PASCO Capstone. ( See "Adding a display ": 56)

## Using Microsoft Word\* (2007/2010)

1. Switch to MathML mode (if needed):
  - a. Click **Insert > Equation > Insert new equation**.
  - b. Select the arrow at the lower right of the tools part of the ribbon.

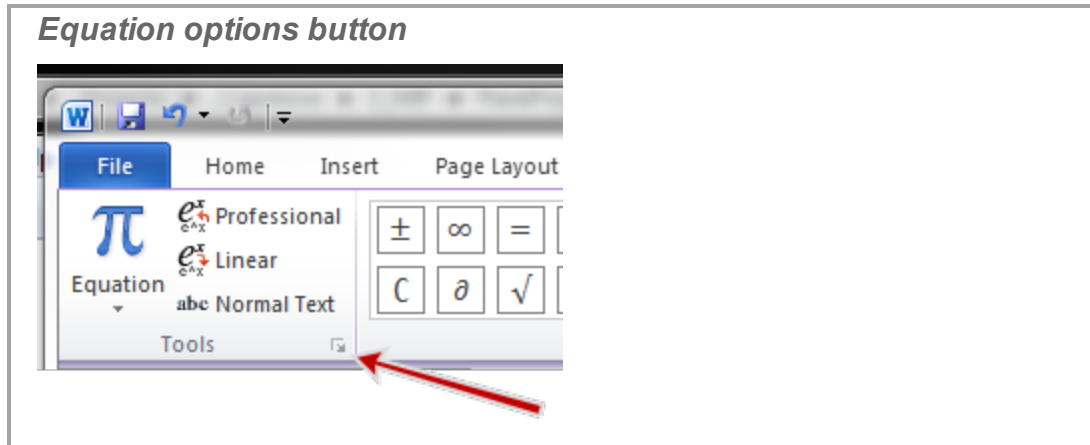

- c. Select **Copy MathML to the clipboard as plain text**

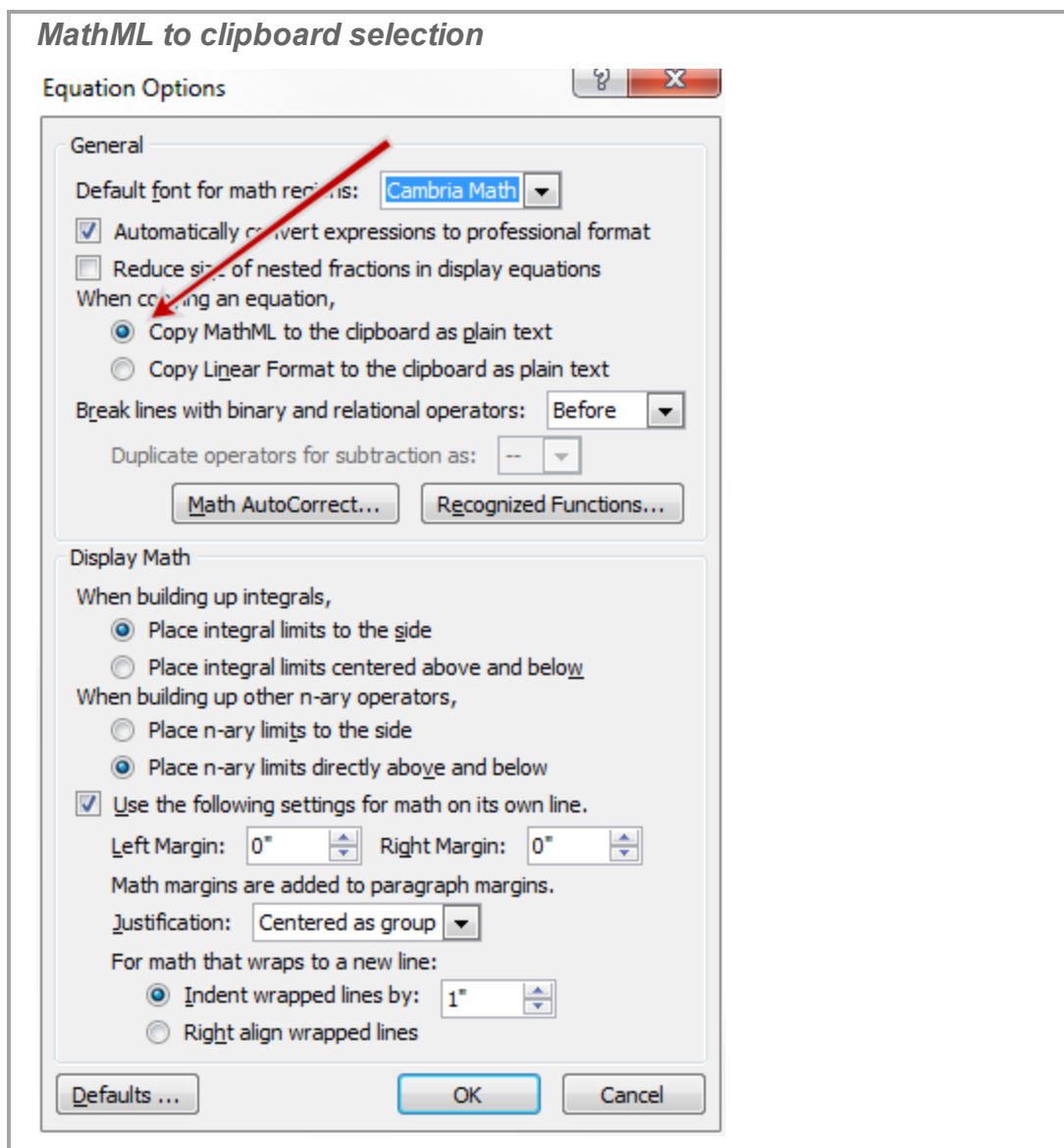

- d. Click **OK**.

2. Type the equation in the programs equation editor.
3. Highlight the equation.
4. Press **CTRL+C** to copy.
5. Switch to a Text Box or Text Entry Box in PASCO Capstone.
6. Place the cursor and press **CTRL+V** to paste.

## How do I... Scope

"How do I compare scope traces from the same sensor when monitoring?": 130

"How do I select the scope trigger?": 131

"How do I pre-trigger?": 134

"How do I freeze and record the scope display?": 135

### How do I compare scope traces from the same sensor when monitoring?


1. Select one of the traces that you would like to compare with later traces by clicking on the symbol corresponding to that plot in the legend.

**Legend: selecting the active plot**


Symbol and color in legend matches plot symbol and color. Click legend

- 1 to make active. (The active plot is shaded in legend and darkest in display.)
- 2 Click in the upper left corner of the legend to clear all active runs.

2. Click **Snapshot**  in the toolbar to capture the visible trace.

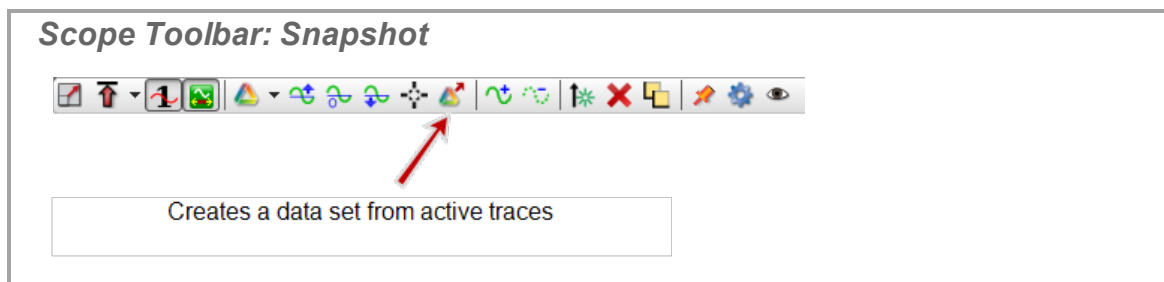

3. Click **Run Selection** .
  - ➡ Multiple runs can display simultaneous when the button is depressed.
4. Click the pull-down arrow (▼) next to **Run Selection** in the toolbar.

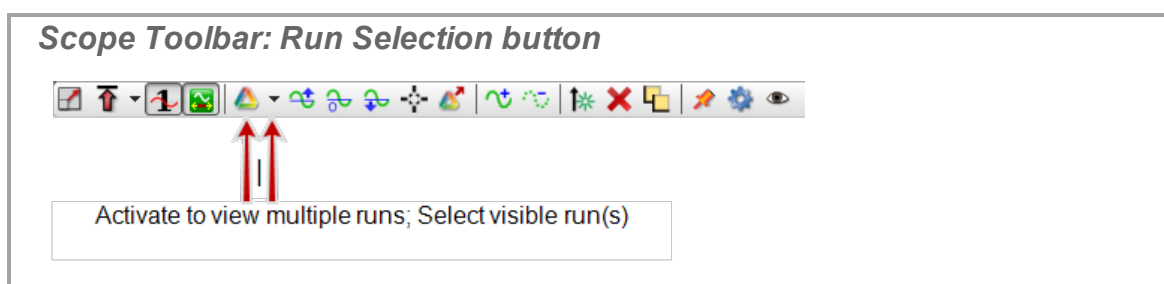

5. Select the measurement trace snapshots that you would like to display.

See "How do I record data in PASCO Capstone? ": 73

## How do I select the scope trigger?


**Note:** If the trigger button in the scope toolbar is depressed, the display of data starts when the trigger condition is satisfied. If the trigger button is not depressed, the display of data starts immediately.

The trigger in the scope applies to the rightmost measurement on the left of the scope plot area.

### I. Select Measurement for Trigger:




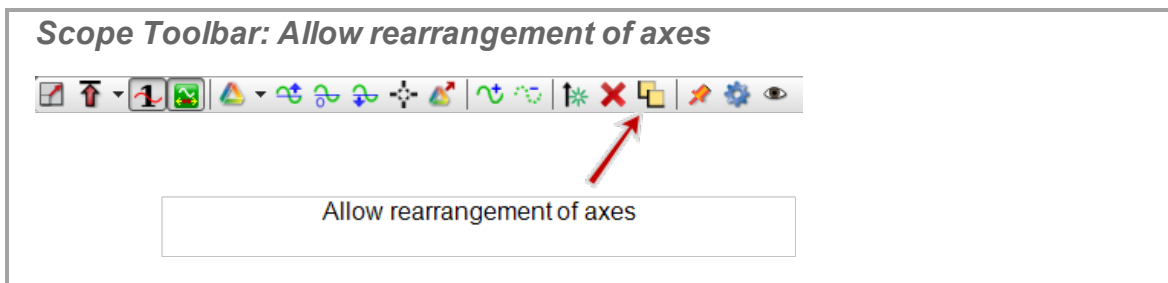
1. Click on the measurement name on the immediate left of the scope.

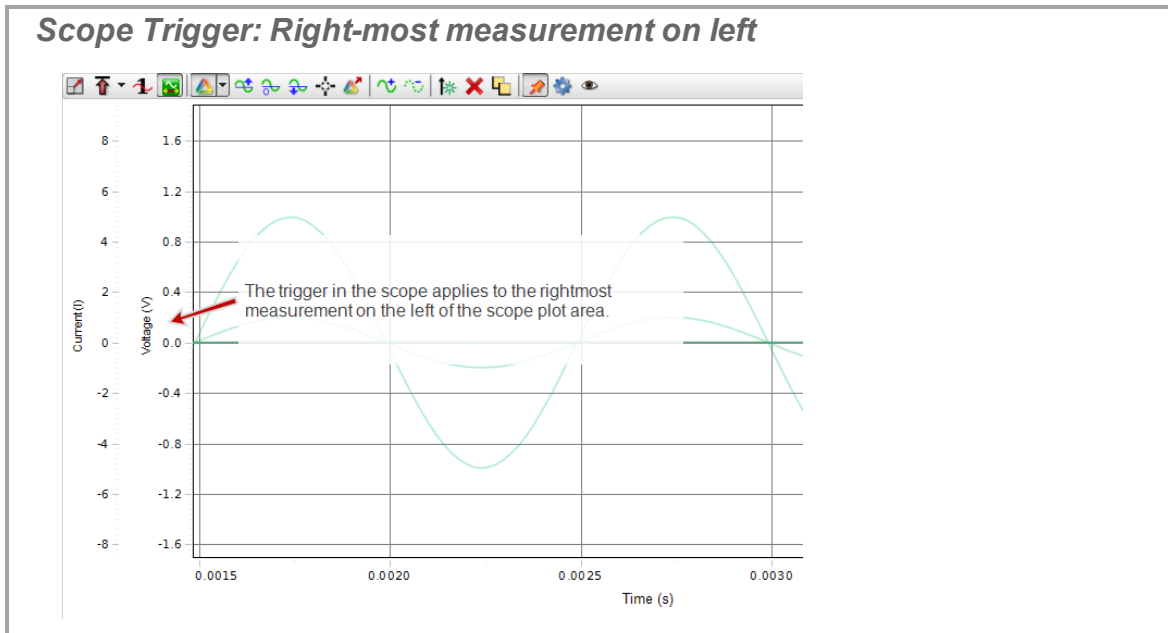


2. Select the measurement for the trigger from the drop-down list.

–Or–

1. Click **Allow rearrangement of axes**  in the toolbar.





2. Click and drag the measurement to use as the trigger to the immediate left of the scope.

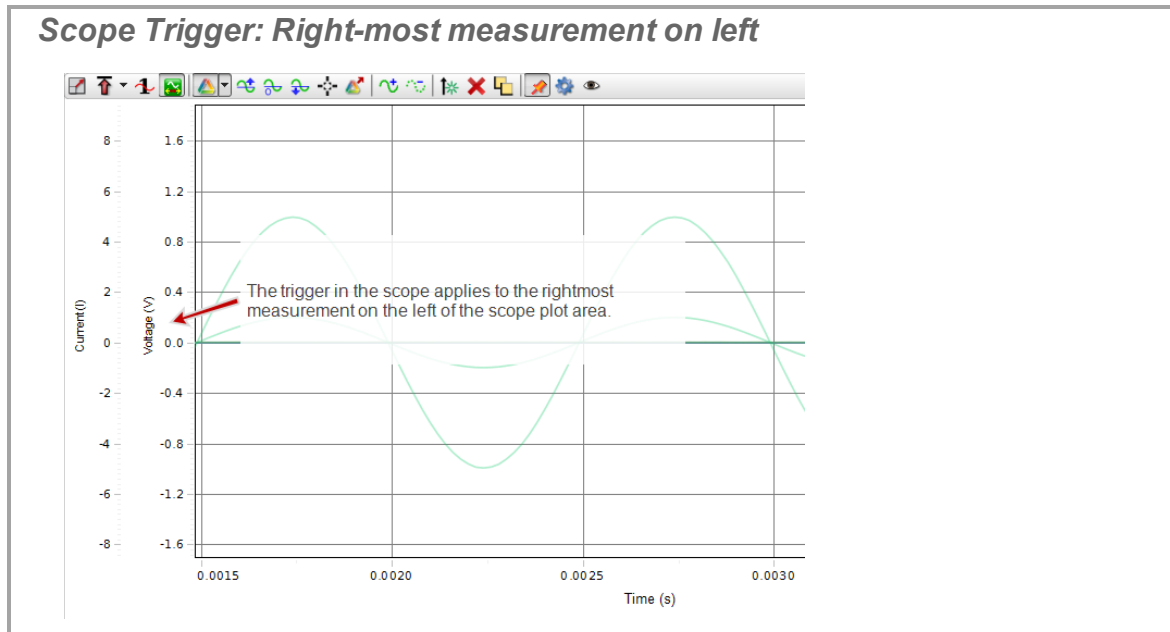

## II. Turn on Trigger:


1. Click **Trigger**  in the toolbar.

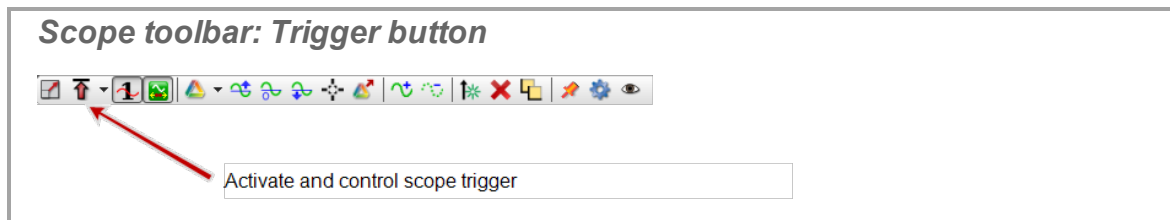

2. Click the pull down menu of the trigger .
  3. Select:
    - **Positive Edge**
    - **Negative Edge**
- ➡ A trigger indicator is displayed on the plot.

### Trigger Indicator


- Indicators point up for Positive Edge and down for Negative Edge triggers.
- Trigger indicator color matches run color.

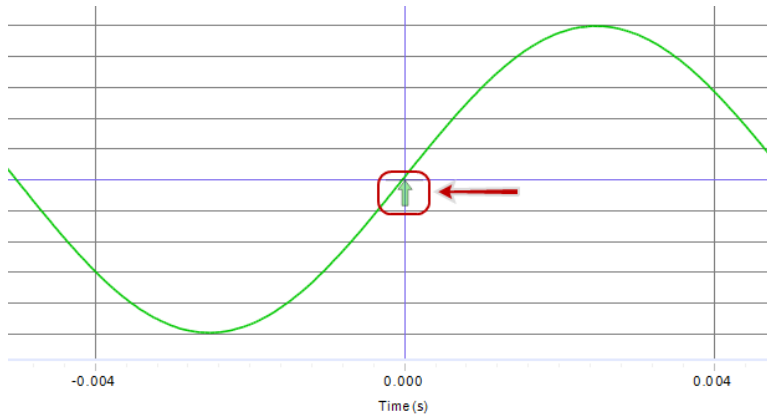


## How do I pre-trigger?

**Note:** An 850 Universal Interface is required to use the pre-trigger functionality.


- Click **Trigger**  in the toolbar.

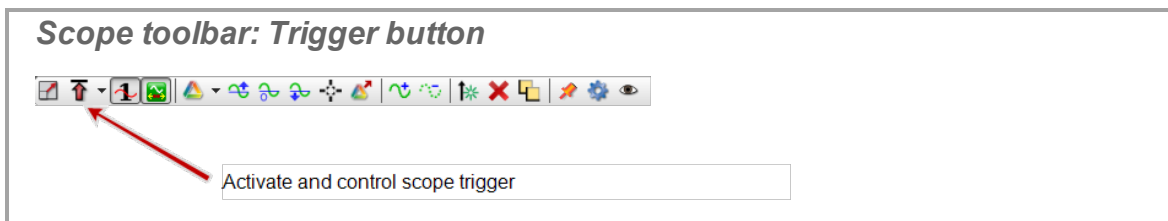

- Click the pull down menu of the trigger .
- Select:
  - Positive Edge**
  - Negative Edge**
- ➡ A trigger indicator is displayed on the plot.

### Trigger Indicator

- Indicators point up for Positive Edge and down for Negative Edge triggers.
- Trigger indicator color matches run color.

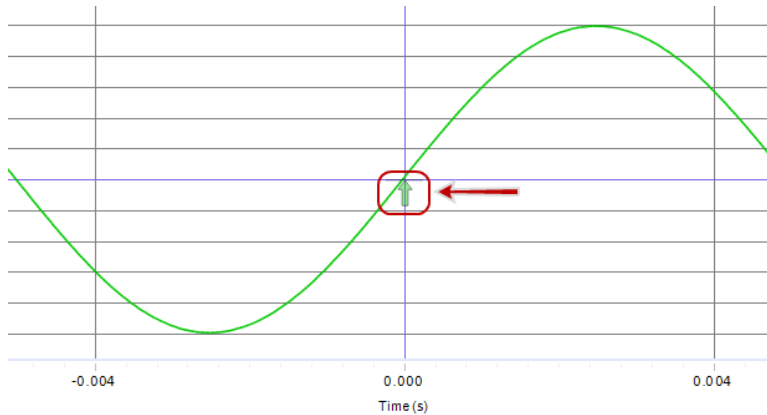


- Pan the scope display to the right, so that the left edge of the viewing area is at the time before the trigger. (See "How do I pan and scale?": 110)

## How do I freeze and record the scope display?

(See "How do I record data in PASCO Capstone? ": 73 for information on Fast Monitor Mode.)


- Select **Fast Monitor Mode** in the Controls Palette.
- Adjust the sample rates of each of the measurements so that the sensors are sufficiently responsive for desired data.

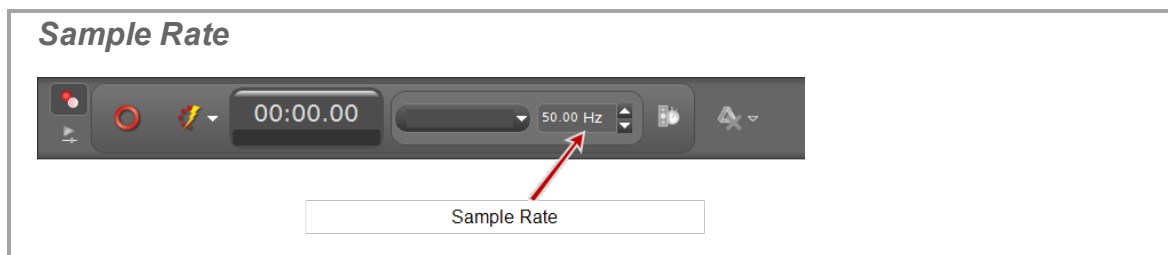


- Select one or more scope or FFT displays. (See "Adding a display ": 56)


**Note:** To record a *single event*, use the scope display and click the single-shot button **1**. (See "What does the Single Shot button do?": 178)


- Click **Monitor** in the Controls Palette.

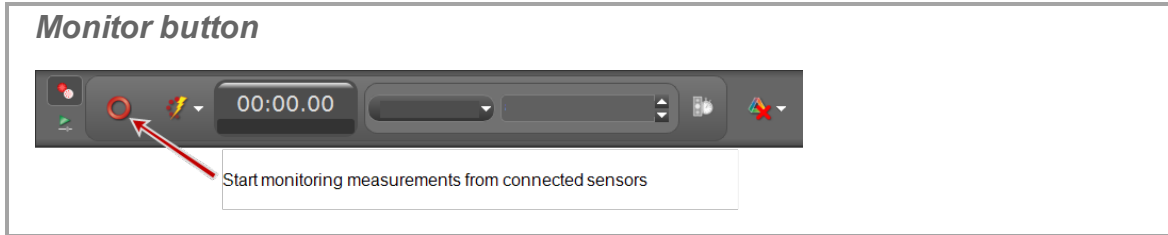

- Click **Stop**  in the Controls Palette to stop observing the data.

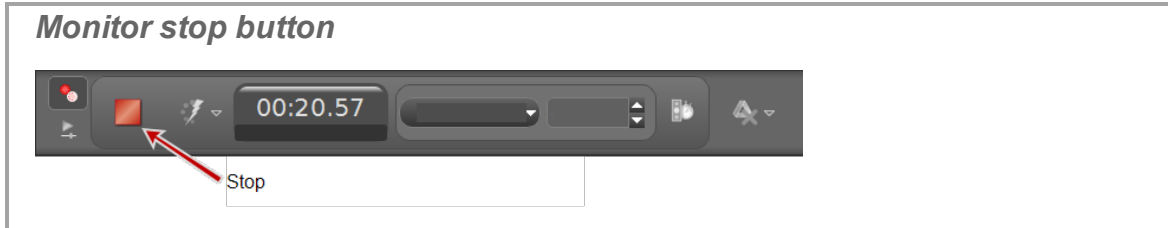

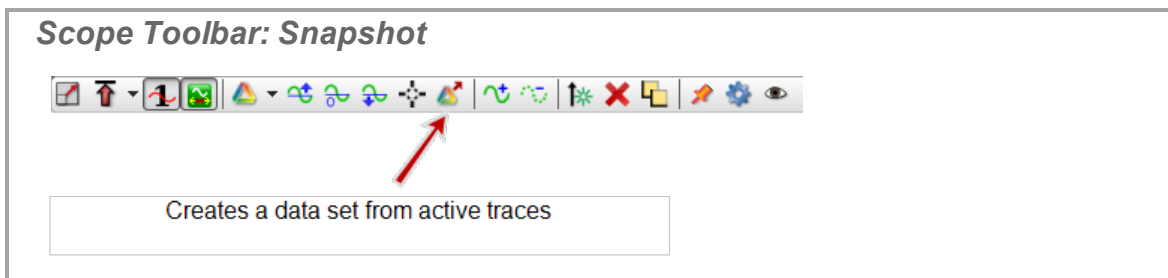

- Capstone freezes the display.
- Scale the data to fit the data to record within the displays. ([See "How do I pan and scale?": 110](#))
- Click **Snapshot**  in the scope toolbar.



- Capstone records the data.

**Notes:**

Only the data that is visible in the plot area is recorded as a run when **Snapshot**  is clicked.

Recorded data appear on the sensor tab in the Data Summary tab () and is selectable from any data display with a default name **Monitor Run--ScopeSnapshot%1** where **%1** is an automatically generated run number.


## How do I... Video

- "How do I add a video source?": 137
- "How do I analyze a video?": 138
- "How do I manually synchronize video with data?": 140
- "How do I adjust the frame rate of a recorded video?": 141
- "How do I adjust the resolution of a video?": 141
- "How do I make files with videos smaller?": 141
- "How do I troubleshoot video problems?": 142

### How do I add a video source?

PASCO Capstone supports most driverless (UVC 1.1) USB video sources, which includes almost all newer web cameras, USB microscopes, and telescopes, and many older USB video sources that require drivers.

#### Procedure

1. Connect a driverless (UVC 1.1) USB video source, such as a web camera, microscope, or telescope.
2. Drag and drop a **Movie Display**  from the **Displays Palette** to a workbook page. (See "Adding a display ": 56)
3. Select one of the four choices presented in the display:

Option	Definition
<b>Open Movie File</b>	Use for Prerecorded QuickTime compatible videos.
<b>Capture Image*</b>	Use to capture a still image. (Do not use for video analysis.)
<b>Record Explanatory Movie*</b>	Use for videos that will not be synchronized to data. (The video records without data recording.)
<b>Record Movie with Synced Data*</b>	Use for a movie that you wish to keep synchronized to sensor data that will be recorded simultaneously. (The video records with data recording.)

\*A live video is displayed in Movie Display.

**Note:** While recording, sound output from the camera is muted by default in order to prevent feedback loops between the microphone and computer speakers. You may

unmute the recorder in the Movie Display properties. There are separate volume/muting controls available for playback and recording.

## How do I analyze a video?

You can manually track the path of objects in any recorded or compatible pre-recorded movie.

### Tips for good results

- PASCO Capstone analyzes in two dimensions. Position the camera plane parallel to the plane of motion. (That is, point the camera perpendicular to the plane of motion.)
- Place the camera approximately at the center of the path of the object of interest.
- Use intense lighting (such as sunlight) to avoid motion blurring that might obscure the image.

### Procedure

1. Add a camera. (See "How do I add a video source?": 137)
2. Add a Movie Display (See "Adding a display ": 56)
3. Select a type of Movie Display:

Option	Definition
<b>Open Movie File</b>	Use for Prerecorded QuickTime compatible videos.
<b>Capture Image*</b>	Use to capture a still image. (Do not use for video analysis.)
<b>Record Explanatory Movie*</b>	Use for videos that will not be synchronized to data. (The video records without data recording.)
<b>Record Movie with Synced Data*</b>	Use for a movie that you wish to keep synchronized to sensor data that will be recorded simultaneously. (The video records with data recording.)
*A live video is displayed in Movie Display.	

➡ If using **Open Movie File**, skip to step 5.

4. Prepare recording.
  - a. Place a physical ruler (or object of known length) in the movie scene to determine scale.
  - b. Ensure the plane of action is parallel to the plane of the lens of the camera.

5. Record movie.

Using **Record Explanatory Movie**:


- a. Start Recording: click **Record**  in the Display toolbar.
- ➔ The video records without data recording.



- b. Stop Recording: click **Record**  in the Display toolbar.


Using **Record Movie with Synced Data**:

- a. Start Recording: click **Record**  in the Controls Palette.
- ➔ The video records with data recording.

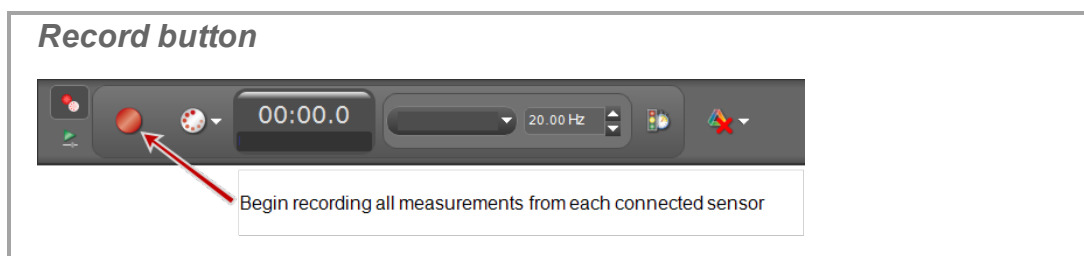


- b. Stop Recording: click **Stop**  in the Controls Palette.

6. Click **Enter video analysis mode**  in the Display toolbar.

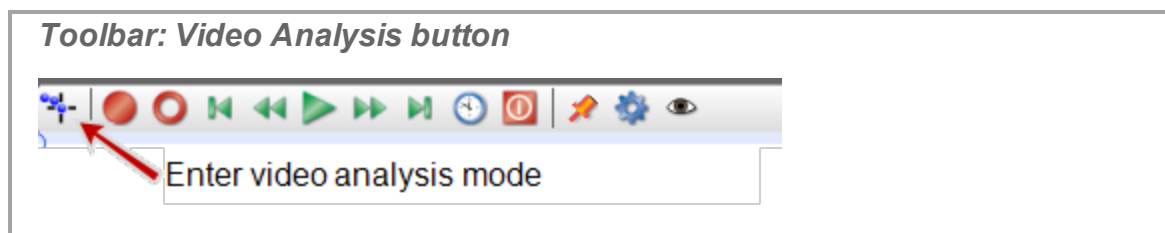



7. Select the data run to use.






8. **Optional: Change the units or length of the scale.**




- a. While in **Video Analysis** mode, click toolbar properties .
- b. Click **Enter video analysis mode**.



- c. Modify as needed: **Real World Length** (Sets Real World Length In Specified Units).
  - d. Modify as needed: **Data Units** (Sets Length Units).
9. Adjust the scale tool to match the scale of the ruler (or object of known length) in the scene.
  10. Click on the object that you would like to track.
    - ➡ Capstone places a + at the point of click and the video will advance to the next frame.
  11. Repeat until the object has been completely tracked.
    - ➡ The video analysis data appear on the **Video Data Summary** tab () in the **Data Summary**, and are selectable from any data display.
  12. Open a Display (such as a Table, Graph, or Histogram). ( See "Adding a display": 56)
  13. Click **<Select Measurement>** in the display and select a video analysis measurement:
    - x-Position
    - y-Position
    - x-Velocity
    - y-Velocity
    - x-Acceleration
    - y-Acceleration

## How do I manually synchronize video with data?

1. Add a Video Display. ( See " How do I add a video source?": 137)
2. Open and existing video recording or record new video.
3. Add a table or graphic display with the data to use. ( See "Adding a display ": 56)
  - ➡ The page contains two displays: video of experiment and numerical data.
4. Click **Playback Mode**  in the Controls Palette.
5. Click the **Play**  in the Controls Palette.
6. Playback the recorded data until an identifiable event occurs. (For example, the point of collision.)
7. Click **Manually synchronize movie with data**  in the Video toolbar.

8. Click **View/adjust current movie playback position**  in the Video toolbar.  
  - ➔ A slider appears under .
9. Adjust the video slider until the video frame corresponds to the time of the identifiable event from step 6.
10. Click **Manually synchronize movie with data**  in the Video toolbar.

## How do I adjust the frame rate of a recorded video?

Frequently high-speed or low-speed video files contain false frame rate metadata. Camera manufacturers deliberately do this so that the video plays back at a easily-viewable frame rate. For PASCO Capstone to synchronize the video correctly, the video rate in the file must be accurate.


### Procedure

Change frame rate metadata of the video file, using a free video editor:

1. Download and install [Avidemux](#).
2. Right click on your video file and select **Open with...**
3. Select **Avidemux**.
4. In Avidemux, select **Video**.
5. Select **Frame Rate**.
6. Check **Use Custom Frame Rate**  and set the desired frame rate.
7. Save the file.
8. For adjusting the phase between the video and the data, see [Tech Note #493](#).

## How do I adjust the resolution of a video?



The settings for all movies are 320×240×30 fps in order to keep file sizes low. A USB 2.0 camera is unlikely to drop frames at this rate. Higher resolutions or frame rates increases the probability of dropped frames.

If the camera supports higher resolutions, you may increase the resolution. ( See "[How do I adjust the frame rate of a recorded video?](#)": 141)


## How do I make files with videos smaller?

PASCO Capstone saves files as uncompressed. Select a compression codec to reduce the file size of recorded videos. The available codecs vary by computer, and some codecs may result in a movie that is not playable within PASCO Capstone.

## Procedure

1. Add a movie to the workbook. ( See " How do I add a video source?": 137)
2. Click on properties  in the video toolbar.
3. Click **Movie Recording**.
4. Click the **Video Compression** drop-down.
5. Select a video compression algorithm.

**Note:** The compression algorithms available depend on the operating system and the software installed by the video camera. Different compression algorithms affect different properties of the video file including frame rate and frame size.

6. Click **OK**.
7. Test: play the video in Capstone.
  -  If the video no longer plays in Capstone, repeat steps and select a different algorithm.

## How do I troubleshoot video problems?

### Dropped Frames and Automatic Synchronization


The video sensor attempts to synchronize the movie with the data. This might not work if too many operating system processes are running simultaneously. (Windows and MAC are not a real time operating systems, and modern installations might have up to hundreds of processes running. The operating system can decide to let other process to have priority, resulting in dropped frames or synchronization issues.)

To improve performance of the auto-synchronization algorithm:

- Increase the sensor sample rate within Capstone
- Terminate unused processes in the operating system
- Temporarily disable virus scanners in operating system
- Turn off the internet connection in operating system

### Capture Sound Issues

Some cameras are not automatically able to assign a matching microphone. If you are missing sound:

1. Click on properties  in the video toolbar.
2. Click the **View/adjust current movie playback position** drop-down.

Choose a microphone. Some computers may not automatically choose the correct microphone. In these cases, you may select a camera and a microphone, and then choose "Save Defaults" to associate the camera and microphone.

Click **OK**.

3. Test video recording and playback.

If no audio, repeat steps to select a different Microphone source.

## Windows Drivers

Open **Device Manager**: **My Computer** > **Properties** > **Manage** > **Device Manager** > **Imaging Devices**

If the camera does not appear in the **Device Manager**, try another USB port or disconnecting and reconnecting the camera to the same USB port.

If the camera appears in the **Device Manager**, click on it and press the delete button, unplug and replug the device into the USB port. Install the latest device driver provided by the manufacturer if requested.

- Devices that support the UVC (USB Video Class 1.1) standard do not need to have any special drivers installed to function in XP-SP3 or Windows Vista/7/8. Such devices appear automatically in **Device Manager** shortly after being connected.
- Older devices will need a driver installed.
- Older 16 Bit Devices (VFW devices) are not plug and play, and will appear to still exist after they have been physically disconnected. To uninstall such devices, right click on the device in Device Manager and select "uninstall driver"

## Testing in Windows

### Windows Vista/Windows 7 Testing

Windows Vista/7 does not include a built-in video viewer. A simple utility that allows you to test a webcam in Windows Vista/7 can be found here:

<http://www.bustatech.com/webcamviewer-a-simple-viewer-for-webcam/>.

You can also use "Skype" or any other video-chat capable program to test a webcam.

### Windows XP Testing

---

**Note:** SP3 (Service Pack 3) is required.

---

1. Close PASCO Capstone
2. Open Windows Explorer.
3. Look under **Scanners and Cameras**.
4. Click on the camera in use.

- ➔ A window opens displaying the camera output.
5. Take a snapshot to test the camera.

## 4 Frequently Asked Questions (FAQ)

### PASCO Capstone

"What is PASCO Capstone?": 147

"Where can I find help documentation and curriculum for PASCO Capstone?": 147

"What are the new features in the latest PASCO Capstone?": 148

"Can PASCO Capstone open DataStudio, SPARKlab, and GLX files?": 151

"Can Capstone files be opened on computers with different operating systems?": 151

"How does Capstone compare to other PASCO applications? ": 152

### Installation and Updates

"What licenses are required?": 160

"Is an MSI available for PASCO Capstone?": 160

"Do I need to update the interfaces?": 161

"Do I need to update the sensor datasheets?": 161

### Setup

"What do I need to have a complete measurement system? ": 162

"Can sensors sample at different rates?": 167

### Data files

"Where is my data saved?": 165

"What is the file export format?": 165

"What is the format for text import?": 166

### Data and Measurements

"Which data is recorded?": 167

"Can sensors sample at different rates?": 167

"What is Continuous Mode?": 168

"What is Fast Monitor Mode?": 169

"Is there a way to delete data points?": 170

"Which data is active?": 172

"How are the curve fit and smoothing calculations determined?": 172

## Calculations and Calibration

"What is the format for calculations?": 174

"What are the equations for calibration?": 174

"Can multiple measurements be calibrated simultaneously?": 175

## Displays

"What is panning and scaling?": 176

"Why are the display toolbars not active?": 177

## Scope

"What is the scope?": 178

"What does the Single Shot button do?": 178

"What is pre-triggering in the scope?": 179

## Video and Text

"What cameras and video formats are supported?": 180

"Why does the text color remain black after picking a color?": 180

## FAQ... PASCO Capstone

"What is PASCO Capstone?": 147

"Where can I find help documentation and curriculum for PASCO Capstone?": 147

"What are the new features in the latest PASCO Capstone?": 148

"Can PASCO Capstone open DataStudio, SPARKlab, and GLX files?": 151

"Can Capstone files be opened on computers with different operating systems?": 151

"How does Capstone compare to other PASCO applications? ": 152

### What is PASCO Capstone?

PASCO Capstone is a feature-rich data acquisition, visualization, and analysis program designed by PASCO scientific for advanced science education. In addition, Capstone contains the tools needed for the creation and execution of structured scientific inquiries, by allowing simple drag-and-drop integration of rich text, graphics, video, and data displays on multiple page displays called workbooks.

Capstone is completely compatible with all PASCO USB Interfaces, including the 850 Universal Interface, Xplorer GLX, SPARKlink, SPARKlink Air, SPARK SLS, USB Link and older interfaces such as the ScienceWorkshop 750 or 500.

 See "Tour the Interface": 1


### Where can I find help documentation and curriculum for PASCO Capstone?

#### Where can I find Capstone documentation?

PASCO Capstone is designed to be as intuitive as possible. In addition the following Help is available:



- **Tool Tips:** hover over a button to view the associated tool tip.

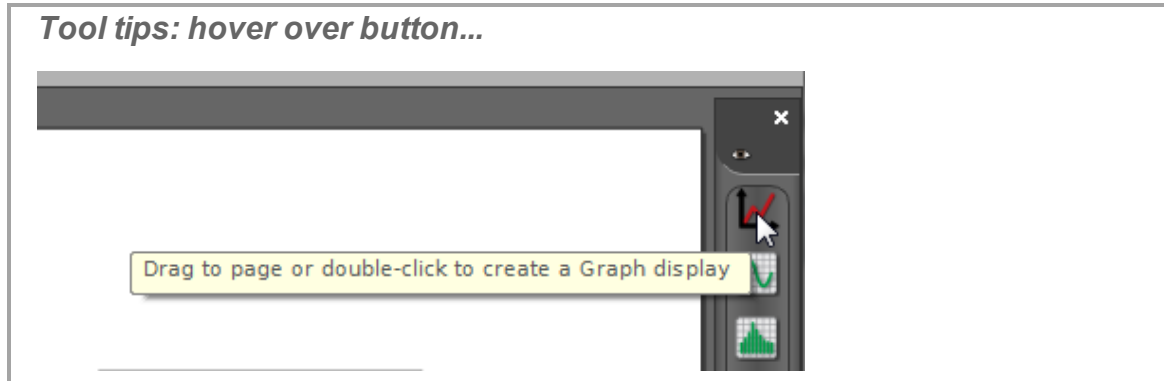


- **Context Help:** helpful information at the bottom of dialog windows. (See ["Tour Context Help": 14](#))
- **Online Help:** searchable instructions on working with PASCO Capstone, frequently asked questions, and links to instructional videos.  
Click **Help > PASCO Capstone Help** (or **F1**)
- **Teacher Support:** for specific questions that are not answered by these resources. (See ["Teacher and Technical Support": ii](#))

### What curriculum is available for Capstone?

A number of Capstone experiments designed for the 850 Universal Interface can be downloaded. (See ["Downloading Capstone experiments": 61](#))

In addition, PASCO Capstone can open and convert many DataStudio 1.9.x and GLX files. (See ["How do I open DataStudio experiments and data?": 68](#))

### What videos are available for Capstone?

A number of Capstone experiments are on the PASCO  channel. (See <http://www.youtube.com/user/pascoscientific>)

### What are the new features in the latest PASCO Capstone?

The following features are new to version 1.1:

- Localization (Input Method Editor support)
- Support for MacOS 10.6 +
- Histogram Display

- Different modes: side-by-side, overlay, stacked
- Statistics
- Coordinates tool
- Video Analysis
  - Manual tracking of objects
  - Video analysis can be done on a movie recorded as “sensor” data or an opened movie file. Can make measurements on a “Captured” image.
- Expansion Port Support
  - Control the visibility of the UI 850 expansion port by clicking on the icon in upper right of Hardware Setup.
  - Click on an analog port and turn it into a Voltage or Peak Amplitude input. These ‘measurements’ will show up in the Data Summary under 850 Expansion Port. For the digital ports, the user can configure either an input or an output.
  - An input produces a State measurement and can also be used as an input for a Custom Timer. An output can be manipulated using a new Digital Output tool that appears.
- Journal
  - Two new buttons on main toolbar: one to take the Snapshot and the other to show/hide the journal.
  - Snapshot options include taking a snap of entire page with palettes, the white content area, or the currently selected display.
  - Within the journal, one can edit the title or annotation, rearrange snapshots, view full-screen, export to HTML, print, delete, etc.
  - The journal can also be re-sized or floated.
- Exclusions/Deletions
  - Can be done either in the Table or Graph.
  - Exclusions remove data from statistics, curve fits, and auto-scaling behavior but are still visible as and X on a graph or as stricken-through numbers in the Table display.
  - Deleting data eliminates the data visually and removes it from any analysis.
  - Excluded or deleted data may be recovered at any time before the Capstone workbook is closed.
- GLX File Import
  - One can either use File Open, or the GLX File Management dialog you get when you click the folder icon next to a picture of the Xplorer GLX in the Hardware Setup window to import data from the Xplorer GLX into Capstone.

- Import/Export Data
  - An Import data... and Export data... options appear in the File menu, allowing the import and export of tab-delimited text. Successfully imported text will appear in the user-entered data tab of the Data Summary window.
- User-defined error bars
  - One can now set error values for individual data points.
  - One can set different positive and negative error values by specifying different measurements for positive and negative (defaults to assuming you want the same for both).
  - One could even use a calculation on a measurement to specify error values.
- Customization options
  - Workbook page shape options, including new default best fit.
  - Can set default font size for workbook displays.
  - Application preferences, including putting your preferences in an external, shared file.
- Improved Print quality
- Linear fits weighted based on error bar values.
- Moving digital probes and analog sensors (via Hardware Setup)
- New Add Similar Measurement menu on graph measurement picker. The ability to add multiple measurements to the same axis has been available for drag/drop but this was not very visible to the user. This new menu item makes that more visible. For example, add eight Force measurements to the same axis. Conserves space and assures that all measurements get scaled with the same axis. The legend becomes the means for identifying what is what. Tries to match the generic measurement type and unit of measurements.
- Graph smoothing slider now has a pop-up with on/off option.
- Graph Curve Fit button has on/off.
- Improved and additional formatting options for the graph coordinates tool.
- Added Peak Amplitude Sensor and Frequency Sensor to the 850 Signal Generator outputs. This feature is useful for LRC sweep of Peak Amplitude versus Frequency.
- Added a Peak Amplitude Sensor to analog inputs for 850, which can be used simultaneously with a Voltage sensor on the same channel.
- Added a new “special” calculation, diff, which, for a single measurement, returns a set of the current value –  $n^{\text{th}}$  previous value.
- Added support for the AirLink2 and SPARKLink Air.

- Displays can 'preview' data in when you have Start condition and Keep sampling.
- Help (English only, initially oriented toward Tasks and Goals)

## Can PASCO Capstone open DataStudio, SPARKlab, and GLX files?

### Can PASCO Capstone open DataStudio files?

**Yes.**

PASCO Capstone can open and convert most DataStudio 1.9 files to Capstone format.

 See ["How do I open DataStudio experiments and data?"](#): 68

### Can PASCO Capstone open SPARKlabs?

**No.**

PASCO Capstone cannot open SPARKlabs.

### Can PASCO Capstone open saved GLX files?

**Yes.**

PASCO Capstone can open and convert most Xplorer GLX files to PASCO Capstone format.

 See ["How do I open DataStudio experiments and data?"](#): 68

### Can PASCO Capstone transfer files from the Xplorer GLX?

**Yes.**

PASCO Capstone can transfer most Xplorer GLX files for use in Capstone.

 See ["How do I import Xplorer GLX data?"](#): 71

## Can Capstone files be opened on computers with different operating systems?

**Yes.**

PASCO Capstone files (.cap) that are created on either a Mac or PC can be opened on another computer running another Capstone-compatible operating system (Mac OS, Windows) with the same version or later of Capstone.

## How does Capstone compare to other PASCO applications?

Specification	PASCO Capstone	DataStudio 1.8.9r10	SPARKvue 2.0
Availability	<a href="http://www.pasco.com">www.pasco.com</a>	<a href="http://www.pasco.com">www.pasco.com</a>	Computer version: <a href="http://www.pasco.com">www.pasco.com</a> Pad version: Apple App Store / Google Play Store
PASCO Part Numbers	UI-5400 UI-5401	CI-6870 CI-6859C CI-6871	Computer version: PS-2400 PS-2401 Pad version: N/A
Development Status	Active	Inactive	Active
Supported Languages	English español de América latina (Latin American Spanish) 简体字 (Simplified Chinese) le français (French) العربية (Arabic)	العربية (Arabic) 繁體字 (Traditional Chinese) 简体字 (Simplified Chinese) Dansk (Danish) English Suomi (Finnish) le français (French) Deutsch (German) italiano (Italian) 日本語 (Japanese) Norsk (Norwegian) português brasileiro (Brazilian Portuguese) Русский (Russian) español cas-	English Dansk (Danish) Deutsch (German) español de América latina (Latin American Spanish) español de México (Mexican Spanish) español castellano (Spanish) français canadien (Canadian French) le français (French) italiano (Italian) Nederlands (Dutch) Norsk (Norwegian) português brasileiro (Brazilian Por-

Specification	PASCO Capstone	DataStudio 1.8.9r10	SPARKvue 2.0
			tugese) português de Portugal (Portuguese) Suomi (Finnish) Svenska (Swedish) Türk (Turkish) Český (Czech) tellano (Spanish) Русский (Russian) Svenska (Swedish) Казак (Kazakh) Türk (Turkish) العربية (Arabic) 繁體字 (Traditional Chinese) 简体字 (Simplified Chinese) J日本語 (Japanese) 한국의 (Korean) ไทย (Thai)
SUPPORTED OPERATING SYSTEMS	Windows XP 32-bit Windows 7 32/64-bit Windows 8 32/64-bit MacOS 10.6-10.8	Windows XP 32-bit Windows 7 32/64-bit (requires special installation) MacOS 9.x, 10.7	Windows XP 32-bit Windows 7 32/64-bit Windows 8 32/64-bit MacOS 10.5-10.8 Apple iOS 5.1 Apple iOS 6.0 Android
INTENDED AUDIENCE	Post-secondary / Advanced High-School Science Educators	Post-secondary / Advanced High-School Science Educators	Elementary/Middle School/ Introductory Secondary Science

Specification	PASCO Capstone	DataStudio 1.8.9r10	SPARKvue 2.0
			Educators
<b>Supported Hardware</b>			
PASPORT			
PS-2100(A) USB Link	YES	YES	YES
PS-2001 PowerLink	YES	YES	YES
PS-2009 SPARKlink	YES	YES	YES
PS-2000 Xplorer	ACQUIRE ONLY	YES	YES
PS-2002 Xplorer GLX External Sensors	YES	YES	ACQUIRE ONLY
PS-2002 Xplorer GLX Sound/Sound Level	YES	YES	NO
PS-2008(A) SPARK SLS	ACQUIRE ONLY	YES	ACQUIRE ONLY
<b>WIRELESS</b>			
PS-2005 AirLink SI	NO	YES	NO
PS-2010 AirLink2	YES	YES	NO
PS-2011 SPARKLink Air	YES	NO	NO
Universal (PASPORT/ScienceWorkshop)			
UI-5000 850 interface	YES	NO	NO
ScienceWorkshop			
ScienceWorkshop 300	NO	YES	NO
CI-6760 ScienceWorkshop 500	via CI-6759A only	Serial / CI-6759A	NO
CI-6560 ScienceWorkshop 700	NO	YES	NO
CI-7500 ScienceWorkshop 750 –SCSI	via CI-6759A only	SCSI / Serial / CI-6759A	NO
CI-7599 ScienceWorkshop 750 – USB	YES	YES	NO
<b>GENERIC DEVICES</b>			
USB video cameras	YES	NO	YES
Microphones	850 or SW750 + sound sensor	Waveport + Built-in Microphone / ScienceWorkshop 750 + sound sensor	Computer version: YES Pad version: As part of video sensor only
<b>THIRD-PARTY HARDWARE</b>			
Ocean Optics Red Tide/Amadeus/ USB 2000 or 4000 spectrometers	NO	NO	Computer version: YES Pad version: NO

Specification	PASCO Capstone	DataStudio 1.8.9r10	SPARKvue 2.0
Ohaus Scout Pro balances	NO	YES	NO
<b>Software Features</b>			
<b>ACQUISITION</b>			
Timer Wizards	YES	NO	NO
Independent sampling rates	YES	YES	NO
Per-page sample rates	YES	NO	NO
Directly Import Data From SPARK SLS	NO	NO	NO
Directly Import Data From Xplorer GLX	YES	YES	NO
<b>EXPORT</b>			
Copy & Paste	From Table	From Summary	NO
Menu Option	All Measurements	One Measurement	All Measurements
<b>IMPORT</b>			
Copy & Paste	To Table	To Summary	NO
Menu Option	YES	YES	NO
<b>CALIBRATION</b>			
Calibration	ScienceWorkshop & select PASPORT Sensors	ScienceWorkshop & select PASPORT Sensors	Select PASPORT Sensors
Calibration types	1 pt slope, 1 pt offset, 2 pt linear	1 pt slope, 1 pt offset, 2 pt linear	1 pt slope, 1 pt offset, 2 pt linear
Ability to return to factory defaults.	YES	NO	NO
<b>Analysis</b>			
Calculations	YES	YES	YES
Immediate evaluation of constant expressions	YES	YES	YES
Reorder calculations	YES	NO	NO
Right-click selection of measurements	YES	NO	NO
Right-click selection of models	YES	NO	NO
Right-click selection of functions	YES	NO	NO
Right-click selection of symbols	YES	NO	NO
Right-click selection of sub-/superscripts	YES	NO	NO
<b>Curve Fits</b>			
Standard Fits	YES	YES	YES



Specification	PASCO Capstone	DataStudio 1.8.9r10	SPARKvue 2.0
Error-weighted Linear Fit	YES	NO	NO
Editable Initial Guesses	YES	YES	NO
Lockable Parameters	YES	NO	NO
User-Defined Functions	YES	YES	NO
Data Manipulation			
Delete Data	YES	YES	NO
Exclude Data	YES	NO	NO
DISPLAYS			
Virtual Oscilloscope Display	YES	YES	NO
Fast Fourier Transform	YES	YES	NO
Automatic video synching	YES	NO	NO
Re-title displays.	YES	YES	NO
X-Y GRAPH			
QuickCalc	YES	NO	NO
Smooth in place	YES	NO	NO
Zooming with the scroll wheel	YES	NO	NO
Snap to origin when panning.	YES	YES	NO
Click/Touch & drag panning	YES	NO	YES
Select measurements from graph labels	YES	YES	NO
Multiple y-axis measurements	YES	YES	YES
Manage y axes	YES	NO	NO
Prediction Tool	NO	YES	YES
Copy a portion of a plot.	NO	YES	NO
Scrolling chart option	YES	YES (UNSMOOTH)	NO
Delete selected data points.	YES	YES	NO
Exclude selected data points.	YES	NO	NO
Change unit directly in the display	YES	NO	NO
Delta Tool	YES (Coordinate Tool Option)	YES	YES
Translucent data values	YES	NO	NO
Preview Data in Keep Mode	YES	YES	YES
TABLE			
Index column	YES	NO	YES
Support for Text Entry	YES	NO	YES
Out-of-order data entry	YES	NO	YES
Non-numeric data entry	YES	NO	YES
Automatic data type switching	YES	NO	NO

Specification	PASCO Capstone	DataStudio 1.8.9r10	SPARKvue 2.0
Cell insertion/deletion	YES	NO	NO
Row insertion/deletion	NO	YES	NO
Drag and Drop to Column Reordering	YES	NO	NO
Cut/Copy/Paste Values	YES	NO	NO
Rename measurement directly	YES	NO	NO
Change unit directly	YES	NO	(MULTIPLE STEPS)
Out-of-order Sampling	YES	NO	NO
Selection Synchronized with other Displays	YES	YES	NO
Insert/delete columns	YES	NO	(MULTIPLE STEPS)
Row Banding	NO	NO	YES
<b>SCOPE</b>			
Un-triggered Sampling	YES	YES	NO
Trigger	YES	YES	NO
Pre-trigger	850 only	NO	NO
Continuous Sampling	YES	NO	NO
Sample Rate Control	YES	NO	NO
Export Trace	YES	YES	NO
Smart Cursor	YES	YES	NO
Multiple Measurements	YES	YES	NO
Multiple Axes	YES	NO	NO
Trace Offset Controls	YES	YES	NO
Compare Traces from the same sensor	YES	NO (Requires graph.)	NO
<b>Histogram</b>			
Count & bin measurement values	YES	YES	NO
Side-by-Side Runs/Sets	YES	YES	NO
Overlay Runs/Sets	YES	YES	NO
Stack Runs/Sets	YES	YES	NO
Autoscaling	YES	YES	NO
Smart Cursor	YES	NO (Does not snap to value.)	NO
Statistics	YES	NO	NO
<b>Bar Graph</b>			
Display Values	NO	NO	YES
Acquire Values Automatically	NO	NO	YES

Specification	PASCO Capstone	DataStudio 1.8.9r10	SPARKvue 2.0
Rename Bins	NO	NO	YES
FFT			
Automatic resolution adjustment	YES	NO	NO
WORKBOOK			
Universal Undo	YES	NO	NO
Data Replay	YES	(Movie Only)	NO
Automatic hiding of tool bars option	YES	NO	NO
Automatic rescaling of workbook	YES	NO	NO
Reordering workbook pages	YES	NO	NO
Live microscope/camera support	YES	NO	YES
Automatic display fitting	YES	NO	YES
Re-sizable displays	YES	YES	NO
Auto-scaling Workbooks	YES	NO	NO
Sensor-Independent Measurement ID	NO	NO	YES
Drag and Drop Image Files	YES	YES	NO
TEXT			
Rich Text Tools	YES	NO	NO
Subscript/Superscript Tools	YES	NO	YES
Symbol insertion	YES	NO	Computer version: NO Pad version: YES
Copy and paste rendered calculations from Microsoft Word.	YES	NO	NO
Drag and drop rendered calculations from calculator	YES	NO	NO
Bulleted list support	YES	NO	NO
<b>Video</b>			
Play pre-recorded video	YES	YES	NO
Capture still images	YES	NO	YES
Acquire and automatically-sync	YES	NO	NO
Automatic playback option	YES	NO	NO
Manually track objects	YES	NO	NO
Automatically generate position, velocity and acceleration data from calibrated track	YES	NO	NO

Specification	PASCO Capstone	DataStudio 1.8.9r10	SPARKvue 2.0
points.			
Measure lengths	YES	NO	NO
Measure angles	YES	NO	NO
Measure radii	YES	NO	NO
JOURNAL			
Snapshots	YES	NO	YES
Reorder Pages	YES	NO	YES
Annotate snapshots	YES	NO	YES
Export as HTML	YES	NO	YES
Print Journal	YES	NO	YES
<b>Function Generation</b>			
850 Universal Interface Support	YES	NO	NO
PI-8127 Support	YES	NO	NO
Peak Amplitude Sensor	850 only	NO	NO
Frequency Sensor	850 only	NO	NO
ScienceWorkshop 750 support	YES	YES	NO
Sweep Frequency	850 only	NO	NO
Voltage Offset	850 only	NO	NO
Generator output based on a measurement calculation	YES (850 or 750)	On/Off only via outputswitch(x) function	NO
Peak Amplitude Virtual Sensor	YES	NO	NO
Frequency Sensor	YES	NO	NO
<b>PROGRAM PREFERENCES</b>			
Adaptive Workbook Page Shape	YES	NO	NO
Custom default display font size.	YES	NO	NO
Application preferences	YES	YES	NO
HELP	YES	YES	YES

## FAQ... Installation and Updates

"What licenses are required?": 160

"Is an MSI available for PASCO Capstone?": 160

"Do I need to update the interfaces?": 161

"Do I need to update the sensor datasheets?": 161

### What licenses are required?

#### What licenses are available for PASCO Capstone?

- Capstone Single User License (UI-5401)
- Capstone Site License (UI-5400)

#### Does the license expire?

**No.**

The license does not expire and entitles you to minor updates, such as the update from version 1.0 to 1.1.

#### Are upgrades free?

**Yes, minor upgrades are free.**

Minor upgrades and bug fixes for the licensed version of PASCO Capstone do not require additional purchase. They may be obtained by downloading the latest trial version from the PASCO scientific website. If the previous version of PASCO Capstone was licensed, the minor upgrade will accept the license key stored in your computer.

If you do not have a valid license key, the trial will expire in 60 days. A new license key may be purchased and entered from the **About PASCO Capstone** menu:

1. Click **Help > About PASCO Capstone**
2. Click **Buy Key** or **Enter Key**

### Is an MSI available for PASCO Capstone?

An independent msi is not available for PASCO Capstone. The Capstone msi is wrapped by InstallShield\*; parameters may be passed to the msi with command line switches. ([🔗 See "How do I automate a network installation?": 66](#)) for information on passing parameters to the msi.

## Do I need to update the interfaces?

PASCO Capstone will prompt to update the firmware if a new version exists. Follow the on-screen instructions.

However, you may encounter a problem with the interface requiring a manual firmware re-installation. ([See "How do I force an update to an interface?": 66](#))

## Do I need to update the sensor datasheets?

**No.**

PASCO Capstone obtains default sensor information from a file that is installed and updated with Capstone.

(DataStudio and SPARKvue do obtain information from the sensors; however, PASCO Capstone ignores any information from the sensor other than the ID.)

## FAQ... Setup


"What do I need to have a complete measurement system? ": 162

"Can sensors sample at different rates?": 167

### What do I need to have a complete measurement system?

#### Required Components

A complete measurement system consists of the following components:

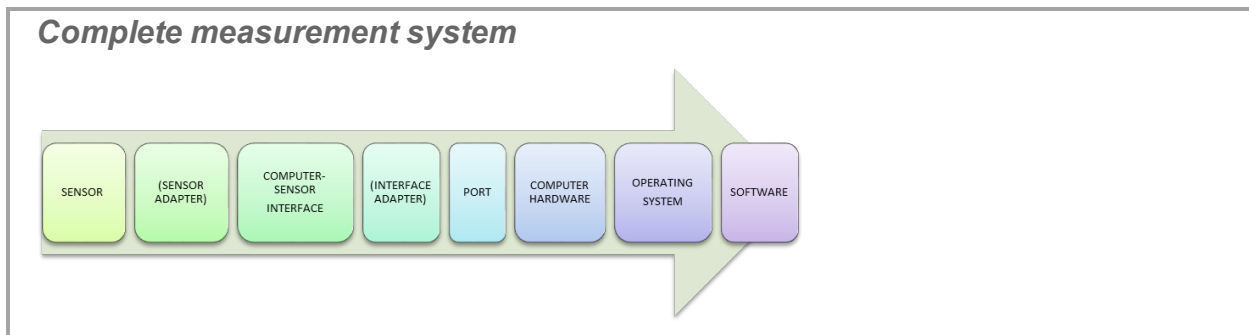


The equipment required depends on the type of sensor used.

#### PASPORT Sensor Connections

PASPORT sensors have part numbers that are prefixed with PS-. These sensors feature auto identification, digital communication, and the ability to measure multiple quantities simultaneously (on some sensors).


Sensors connect as follows:

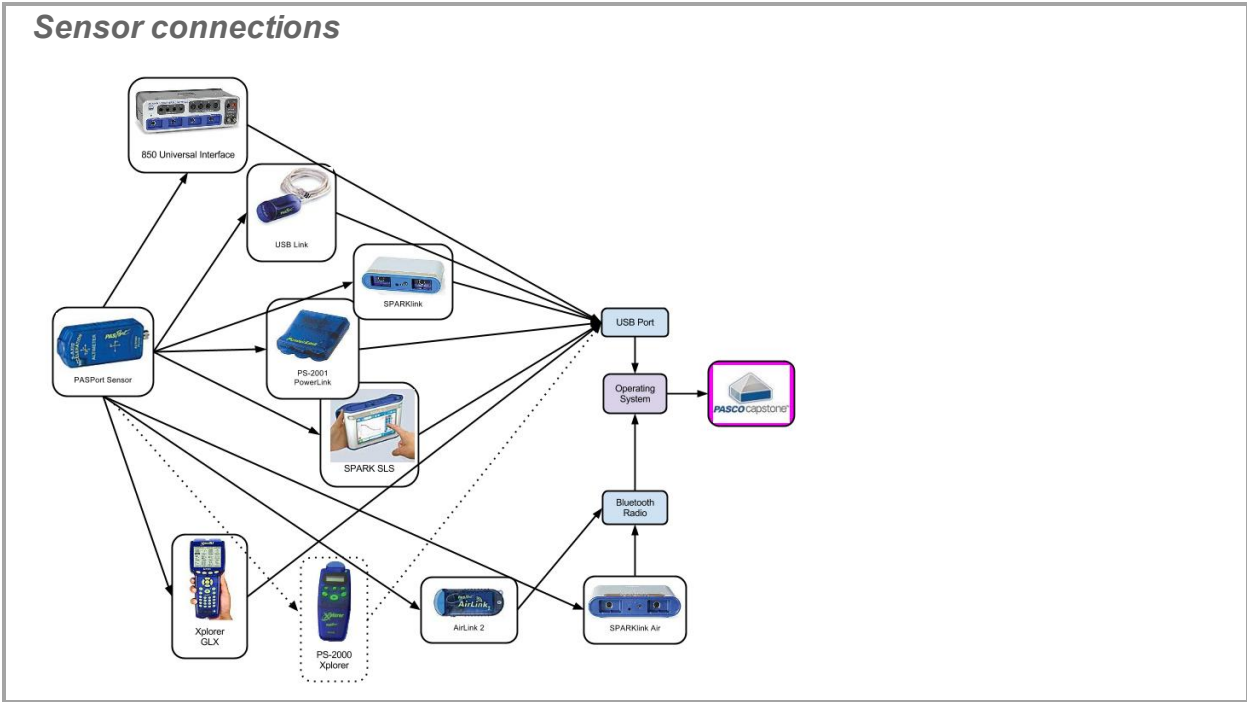


## ScienceWorkshop Sensor Connections

ScienceWorkshop sensors have part numbers that are prefixed with CI-. There are two classes of ScienceWorkshop sensors: analog and digital. The analog sensors may be identified by their round 8-pin DIN connectors. The digital sensors are identified by their 3.5 mm Tip-Ring-Sleeve (headphone) connectors.

These sensors connect as follows:

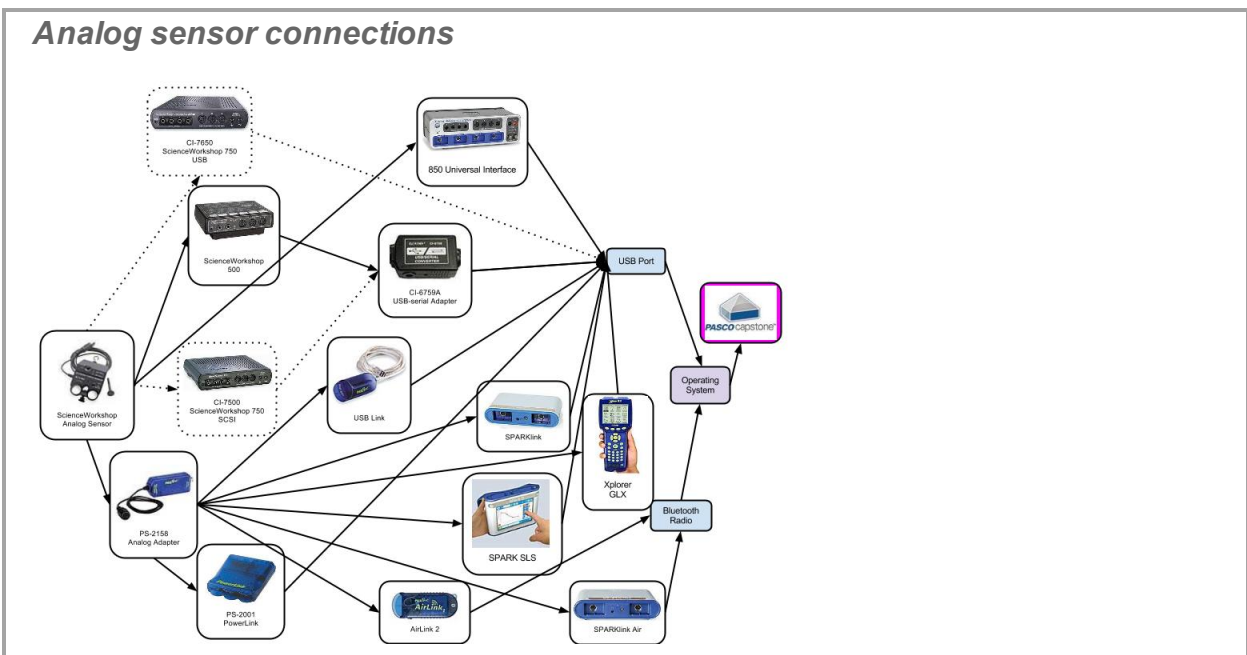



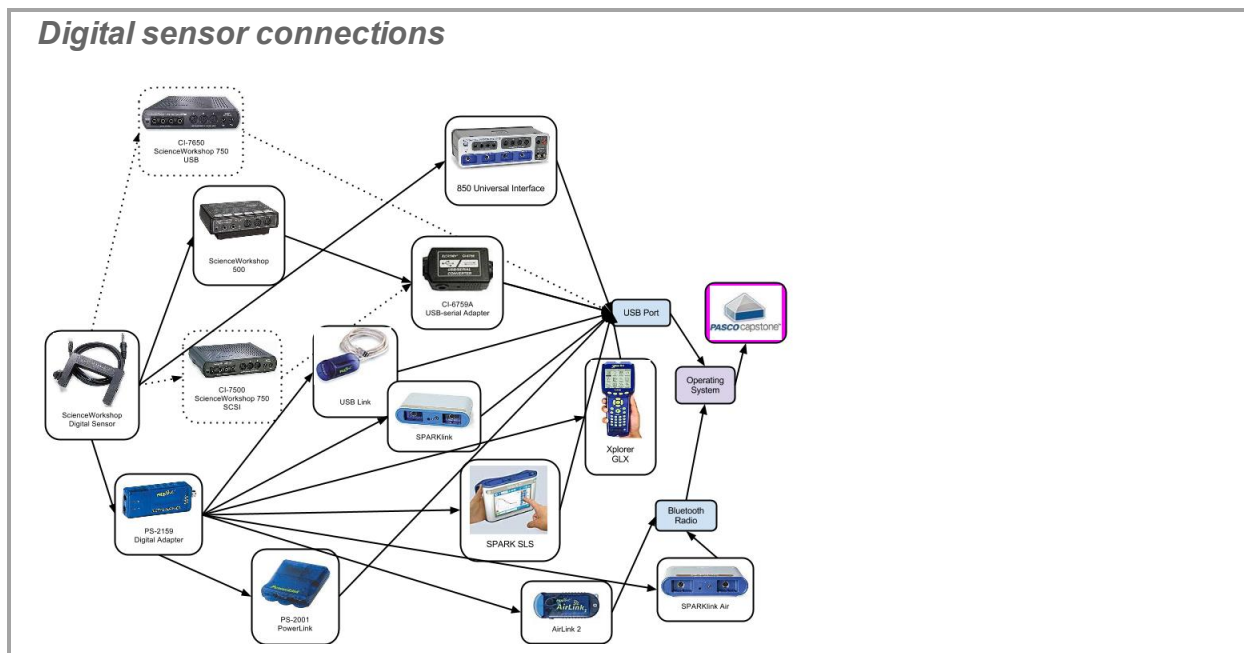

## Can sensors sample at different rates?

**Yes.**

PASPORT sensors may have sampling rates different from the sampling rates of other PASPORT or ScienceWorkshop sensors.

**Note:** All ScienceWorkshop sensors sample at the same rate. PASPORT sensors can sample at different rates.

## FAQ... Data files

"Where is my data saved?": 165

"What is the file export format?": 165

"What is the format for text import?": 166

### Where is my data saved?

Recorded data is contained within the PASCO Capstone file (identified by the **.cap** extension).

**Note:** Data is not permanently saved unless the Capstone file is saved or the data exported. If PASCO Capstone closes, without saving or exporting the data, all data will be lost.

### What is the file export format?

PASCO Capstone exports data in tab or comma delimited files with the following information:

```
Run [ID]/Set [ID]
[Measurement] ([Unit])
Value
Value
⋮
```

**Note:** The precision of the exported data matches the numerical properties of the measurement. (See ["How do I change the numerical format of data?": 81](#))

#### Example

Run #1	Run #1	Run #1	Run #2	Run #2	Run #2
Time (s)	Force (N)	Position (m)	Time (s)	Force (N)	Position (m)
0.000	0.61	0.16	0.000	0.63	0.17
0.050	0.42	0.26	0.050	0.44	0.26

## What is the format for text import?

Imported data are listed in the **User-Entered Data Summary** tab () of **Data Summary**.

The recommended format for file import is tab-delimited text with the measurement names and units on the first line:

$$MN_1(u_1) \rightarrow MN_2(u_2) \dots \rightarrow MN(u_n) \text{ [CR]}$$

$$M_1V_1 \rightarrow M_2V_1 \rightarrow \dots M_nV_1 \text{ [CR]}$$

$$M_1V_2 \rightarrow M_2V_2 \rightarrow \dots M_nV_2 \text{ [CR]}$$

$$\vdots$$

$$M_1V_m \rightarrow M_2V_m \rightarrow \dots M_nV_m \text{ [CR]}$$

**Where:**

$MN_n$   $\equiv$  measurement name of the  $n^{\text{th}}$  measurement

$(u_n)$   $\equiv$  unit of measure of the  $n^{\text{th}}$  measurement—leave off if no units

$M_nV_m$   $\equiv$   $m^{\text{th}}$  value of the  $n^{\text{th}}$  measurement

$\rightarrow$   $\equiv$  tab

**[CR]**  $\equiv$  carriage return

**Note:** PASCO Capstone may not recognize data that does not exactly fit the recommended file import format.

**Example text file:**

<b>animal</b>	<b>family</b>	<b>speed (km/h)</b>
Peregrine Falcon	Falco peregrines	389
Golden Eagle	Aquila chrysaetos	320
Gryfalcon	Falco rusicolus	208

 See "How do I import data from Microsoft Excel\*? ": 69

## FAQ... Data and Measurements

"Which data is recorded?": 167

"Can sensors sample at different rates?": 167

"What is Continuous Mode?": 168

"What is Fast Monitor Mode?": 169


"Is there a way to delete data points?": 170


"Which data is active?": 172

"How are the curve fit and smoothing calculations determined?": 172

### Which data is recorded?

PASCO Capstone records data for each available measurement (*All data is recorded for immediate or future use in displays, calculations, and analysis.*) and associated time data at the defined sample rate(s).

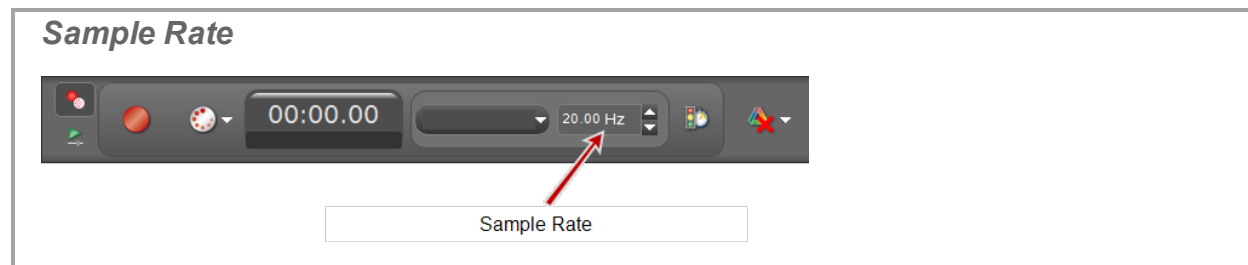

Data appear on the sensor tab in the Data Summary tab () and is selectable from any data display with a default name **Run #%1**, where %1 is an automatically generated run number.

 See "How do I record data in PASCO Capstone? ": 73

### Can sensors sample at different rates?

Yes.


PASPORT sensors may have sampling rates different from the sampling rates of other PASPORT or ScienceWorkshop sensors.


**Note:** All ScienceWorkshop sensors sample at the same rate. PASPORT sensors can sample at different rates.

## What is Continuous Mode?

In this mode (See ["How do I record data in PASCO Capstone? "](#): 73):

- Data continuously records as soon as **Record**  is pressed in the Controls Palette.

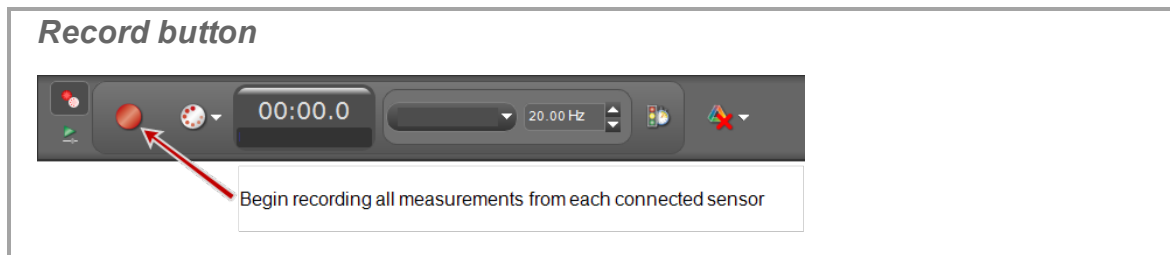

- Data displays at the specified sampling rate(s) either until the hardware is no longer able to process the data (at which time data collection will stop with an error message) or until the **Stop**  is pressed.

### Maximum Frame Rate

The maximum **frame rate**<sup>1</sup> is directly related to the processing speed of the host and inversely related to the number of samples per frame:  $S_F$ .

**Where:**

$$S_F = R_s \times T$$



$S_F$   $\equiv$  Samples per frame

$R_s$   $\equiv$  Sampling rate (displayed in the Controls Palette)


$T$  = time specified on the right of the plot area window

### Automatic Sample Rate ( $R_s$ )

Select the auto-sample-rate-adjustment button  to enable automatic sampling rate:


- The sampling rate will increase when zooming into the time axis or when increasing the number of samples per frame by using the increase-samples-per-trace button .
- The sampling rate will decrease when zooming out of the time axis or when decreasing the number of samples per frame by clicking the decrease-samples-per-trace button .

### Manual Sample Rate ( $R_s$ )

Deselect the auto-sample-rate-adjustment button  to disable automatic sampling rate:

<sup>1</sup>The rate at which these frames of data are displayed.

- The sampling rate is set in the Controls Palette.

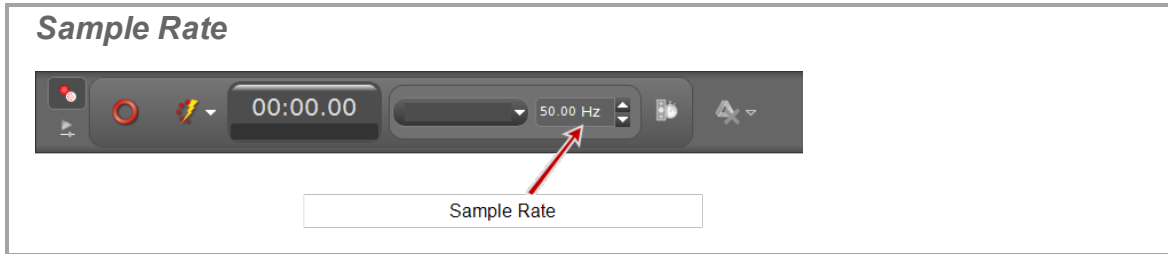

## What is Fast Monitor Mode?

In this mode (See "How do I record data in PASCO Capstone? ": 73):

- The scope display behaves like a digital storage oscilloscope.
- Data is manually recorded using **Snapshot** (See "How do I record data in PASCO Capstone? ": 73)

### Notes:

Only the data that is visible in the plot area is recorded as a run when **Snapshot** is clicked.

Recorded data appear on the sensor tab in the Data Summary tab, and is selectable from any data display with a default name **Monitor Run--ScopeSnapshot%1** where **%1** is an automatically generated run number.

- The scope will not drop samples within a frame due to an excessive sample rate.
- The trigger, by definition, occurs at a scope time of 0.00.
- The scope displays complete frames of data. (The maximum **frame rate**<sup>1</sup>, depends on the processing speed of the host computer.)

### Maximum Frame Rate

The maximum **frame rate**<sup>2</sup> is directly related to the processing speed of the host and inversely related to the number of samples per frame:  $S_F$ .

**Where:**

$$S_F = R_s \times T$$

$S_F \equiv$  Samples per frame

$R_s \equiv$  Sampling rate (displayed in the Controls Palette)



$T =$  time specified on the right of the plot area window

### Automatic Sample Rate ( $R_s$ )


<sup>1</sup>The rate at which these frames of data are displayed.

<sup>2</sup>The rate at which these frames of data are displayed.


Select the auto-sample-rate-adjustment button  to enable automatic sampling rate:

- The sampling rate will increase when zooming into the time axis or when increasing the number of samples per frame by using the increase-samples-per-trace button .
- The sampling rate will decrease when zooming out of the time axis or when decreasing the number of samples per frame by clicking the decrease-samples-per-trace button .

### Manual Sample Rate ( $R_s$ )

Deselect the auto-sample-rate-adjustment button  to disable automatic sampling rate:

- The sampling rate is set in the Controls Palette.

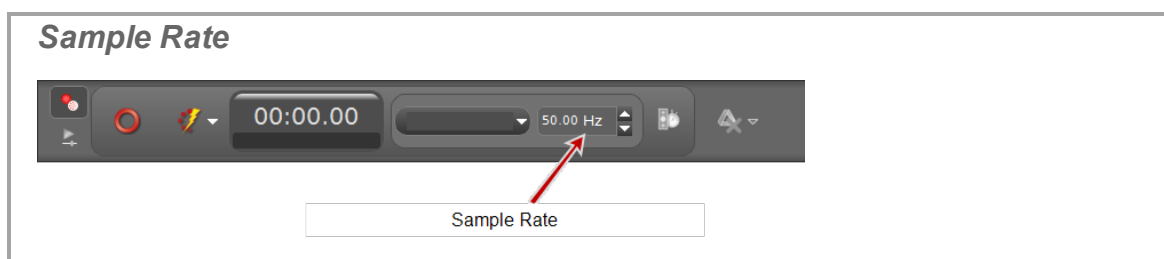


### Is there a way to delete data points?

Yes.

In PASCO Capstone version 1.1 or later, data points can be deleted or excluded.

- Neither excluded nor deleted points are included in curve fits or statistical calculations.
- If data points are excluded, they remain in the displays and are marked:

- with an X in the plot

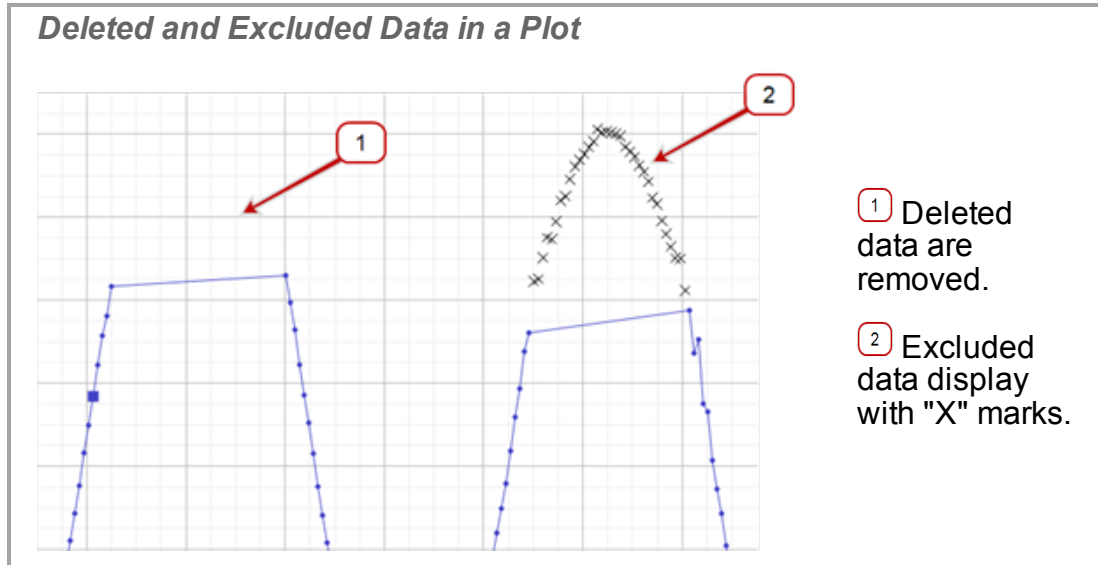

- with a strike-through in the table

**Deleted and Excluded Data in a Table**

0.900	0.30
0.950	<del>0.26</del>
1.000	<del>0.25</del>
1.050	<del>0.19</del>
1.100	<del>0.18</del>
1.150	<del>0.18</del>
1.200	<del>0.18</del>
1.250	<del>0.21</del>
1.300	<del>0.28</del>
1.350	0.35
1.400	0.42
2.200	0.42


1 Deleted data are removed. Note the gap in time data.

2 Excluded data display with strikethrough.

- If data points are deleted, they disappear from all of the displays. Either of these functions may be reversed with the undo button.

[See "How do I exclude, delete, and restore selected data?": 84](#)



**Note:** Deleted data are removed from the data set and the data cannot be restored at a later time. You can immediately undo a delete action by clicking **Undo** .

**Note:** Excluded data are masked in the data set and are not used by Capstone for data operations including display and calculations. Excluded data can be restored at any time.

## Which data is active?

By default, the last selected measurement of the last data run is the active data for statistics and other actions.

 See "How do I select active data?": 111

## How are the curve fit and smoothing calculations determined?

### How are the errors in the parameters of the linear curve fits calculated?

The linear model error estimations were derived assuming that there is no error in the reported measurements and tends to underestimate the actual error in the parameters. The errors are calculated as follows:

$$\Delta m = \sigma \sqrt{\frac{N}{\Delta}}$$

$$\Delta b = \sigma \sqrt{\sum \left( \frac{x_i^2}{\Delta} \right)}$$

**Where:**

$m \equiv$  slope

$b \equiv$  intercept

$N \equiv$  number of values  $(x_i, y_i)$

...and...

$$\sigma \equiv \frac{\sqrt{\sum (y_i - b - mx_i)^2}}{N - 2}$$


$$\Delta \equiv N \times \sum (x_i^2) - \sum (x_i)^2$$

### What is the algorithm used by the graph smoothing tool?

The graph smoothing tool uses a center-weighted [Savitzky–Golay smoothing](#) algorithm of degree 4. The the value displayed in the Data Smoothing Slider **n** is an odd value

representing the total number of values in the original data set that are used in determining each smoothed value.

**Data Smoothing Slider**



1 Slider

2 **n** (used in the [Savitzky–Golay smoothing](#) algorithm)

Coefficient weights are used, such that the values closest to the value being smoothed usually have a greater weight, and thus a greater influence on the determining the resulting smoothed value. These coefficients are symmetric unless there are fewer than  $\frac{n-1}{2}$  unsmoothed values to both sides of the smoothed value.

## FAQ... Calculations and Calibration

"What is the format for calculations?": 174

"What are the equations for calibration?": 174

"Can multiple measurements be calibrated simultaneously?": 175

### What is the format for calculations?

Calculations have the format:

**(variable name) = (valid functional expression)**

The variable name may contain symbols, common subscripts and superscripts, and spaces.

If you enter a functional expression without **(variable name) =**, Capstone adds a generic name and "=". For example: **Calc1=** for variables and **Constant1=** for constants.

Calculations that involve only constants are immediately evaluated in the user-entry area. Calculations involving measurements are rendered in MathML. ([📎 See "How do I see the result of a calculation?": 103](#))

#### Examples

**Constant1=2\*3**

**Calc1=[Acceleration (m/s<sup>2</sup>) ] ^2**

**r^2=sin ( [Angle (rad) ] )**

**ϕ=integral ( [Force (N) ] , [Time (s) ] )**

**Resultant Acceleration=sqrt ( ( [Acceleration- x (m/s<sup>2</sup>) ] ^2+[Acceleration- y (m/s<sup>2</sup>) ] ^2+[Acceleration- z (m/s<sup>2</sup>) ] ^2 ) )**

### What are the equations for calibration?

The actual value,  $y(x)$  is calculated using measured value  $x$  and the calibration points  $(x_i, y_i)$ , where  $x_i$  is the  $i^{\text{th}}$  measured value and  $y_i$  is the corresponding actual (standard) value:

### 1 Point Offset

$$y(x) = (y_1 - x_1)x$$

**Examples:** used internally by the force sensors, accelerometers

### 1 Point Slope

$$y(x) = \left(\frac{y_1}{x_1}\right)x$$

**Examples:** dissolved oxygen, type-K temperature, conductivity, drop counter

### 2 Point Linear

$$y(x) = \left(\frac{y_2 - y_1}{x_2 - x_1}\right)x + \frac{x_2 y_1 - y_2 x_1}{x_2 - x_1}$$

**Examples:** pH and other ion selective electrodes

## Can multiple measurements be calibrated simultaneously?

**Yes.**

You may calibrate multiple measurements simultaneously if they are of the same type and exposed simultaneously to the same standards when performing the calibration.

**Note:** Simultaneous calibration is not recommended for pH, ISE, ORP, Conductivity, stainless-steel temperature sensors without insulating sleeves, or Dissolved Oxygen sensors because these sensors will not read correctly if the probes are placed in the same solution simultaneously.


 See "How do I calibrate a sensor?": 94

## FAQ... Displays

"What is panning and scaling?": 176

"Why are the display toolbars not active?": 177

### What is panning and scaling?

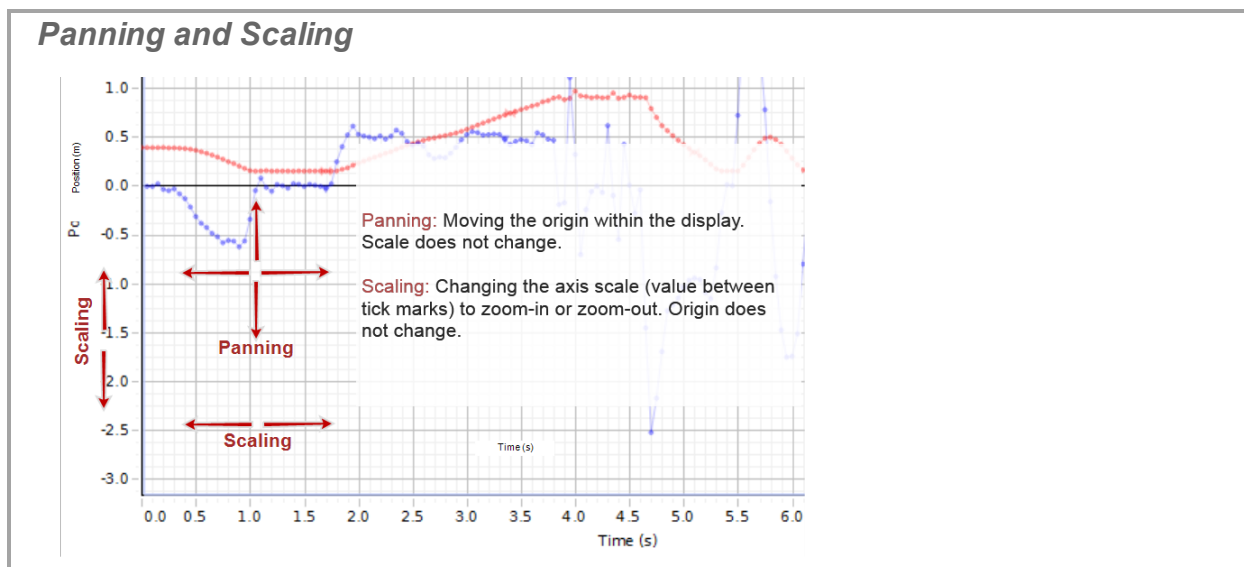

PASCO Capstone supports several panning and scaling methods:

- user (manual) **Scaling<sup>1</sup>** and **Panning<sup>2</sup>**
- scaling based on maintaining the visibility of new data
- scaling and panning the data to maximize the visibility of existing data

See "How do I pan and scale?": 110.

**What is Panning** *Moving the origin within the display. Scale does not change.*

Panning is the process of changing the value of the constant P in the Scaling and Panning Equation.

<sup>1</sup>Changing the axis scale (value between tick marks) to zoom-in or zoom-out. Origin does not change.

<sup>2</sup>Moving the origin within the display. Scale does not change.

**What is Scaling** *Changing the axis scale (value between tick marks) to zoom-in or zoom-out. Origin does not change.*

Scaling is the process of changing the value of S in the Scaling and Panning Equation. The process of increasing the value of the scaling factor S, is said to be zooming in and decreasing the value of the scaling factor is said to be zooming out.

**What is Zooming-in/out around a point** *Changing the scale and origin around a single point.*

You can simultaneously scale and pan such that a point other than the origin (for example:  $(m_{i0}, m_2)$ ) remains stationary in the plot area when scaling. This process is the equivalent of scaling by S and panning by  $m_{i0} \times S$  for each coordinate.

### Scaling and Panning Equation

$$x_i = (m_i - m_{i0}) \times S + P$$

**Where:**

$x_i$   $\equiv$  virtual value of the  $i^{\text{th}}$  coordinate within the plot area

$m_i$   $\equiv$  the  $i^{\text{th}}$  measurement value

$m_{i0}$   $\equiv$  the coordinate value of the scaling stationary point

S  $\equiv$  scaling factor

P  $\equiv$  panning factor

### Why are the display toolbars not active?

Toolbars are only active when a single display is active, and will only apply to the active display.

For displays with multiple plots such as for Graph or FFT displays, the toolbar actions apply to the active plot within the display.

 See ["How do I select active data?": 111](#)

## FAQ... Scope

"What is the scope?": 178

"What does the Single Shot button do?": 178

"What is pre-triggering in the scope?": 179

### What is the scope?



The scope is designed to view high-frequency oscillating data or single high-speed events that the graph display may not be able to capture.

**Maximum Sampling Rate:** The maximum supported sampling rate is a function of the computer-sensor interface that is connected to the computer (for example: 850 Universal Interface → 10 MHz, ScienceWorkshop 750 → 250 kHz).

Features	See
Continuous Mode and Fast Monitor Mode	"What is Continuous Mode?": 168 "What is Fast Monitor Mode?": 169
Single shot	"What does the Single Shot button do?": 178
Multiple triggers	"How do I select the scope trigger?": 131
Pre-trigger	"How do I pre-trigger?": 134
Image freeze	"How do I freeze and record the scope display?": 135


### What does the Single Shot button do?


The Single Shot button is useful for fast, non-repetitive events, such as measuring the time between an impulsive sound and its echo to determine the speed of sound.

When the Single Shot button  is depressed on the scope toolbar and **Monitor**  is pressed on the Controls Palette, the scope displays one frame of scope data and then stops monitoring.



**Monitor button**

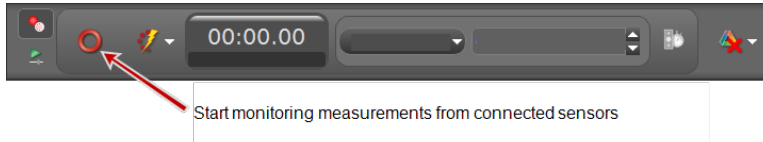

**Note:** If the trigger button  in the scope toolbar is depressed, the display of data starts when the trigger condition is satisfied. If the trigger button is not depressed, the display of data starts immediately.

**Note:** Fast Monitor Mode is required for the Single Shot function.

**Note:** Single Shot is limited by the hardware buffer size of the interface (~ 2000 samples for the 750 interface and ~ 16000 sample for the 850 interface).

## What is pre-triggering in the scope?

The 0.00 second measurement on the scope is referenced to the time of the scope triggering event. Recording and displaying data that happened before a triggering event is called pre-triggering.

**Note:** An 850 Universal Interface is required to use the pre-trigger functionality.

 See "How do I pre-trigger?": [134](#)



## FAQ... Video and Text

"What cameras and video formats are supported?": 180

"Why does the text color remain black after picking a color?": 180

### What cameras and video formats are supported?

#### Supported cameras

- PASCO Capstone supports the UVC (USB Video Class 1.1) standard. Cameras that support UVC do not require additional drivers.
- UVC 1.1 cameras are recommended; however, some older cameras may work with Capstone with the drivers developed by the manufacturer.

 See ["How do I troubleshoot video problems?"](#): 142

#### Supported video formats

- PASCO Capstone will record and playback video recorded in Capstone without any supporting software.
- Capstone requires the [latest version of the Apple QuickTime\\* player](#) to playback movies that were not recorded in Capstone. Once QuickTime is installed, Capstone can playback all movies with the **mov** file extensions and most movies with the **mp4** or **avi** extensions.

### Why does the text color remain black after picking a color?

PASCO Capstone uses the Hue-Saturation-Value (HSV) model to determine the color. The default value of the color value (brightness) defaults to **0** and the color displays black.

To display the selected color: move the brightness slider for the desired brightness.